DECLARACIÓN DE CONFORMIDAD CE  
CE DECLARATION OF CONFORMITY  
DÉCLARATION DE CONFORMITÉ CE

El fabricante | The manufacturer | Le fournisseur:

AUXILAB, S.L.

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Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita vortex mixer 681/10. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

**TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.**

Please bear in mind the following:

- ◆ This manual is inseparable from the Nahita oscillating stirrer, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the equipment avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the equipment to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the oscillating stirrer.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to [asistencia@auxilab.es](mailto:asistencia@auxilab.es) or by telephone: +34 807 117 040 (0.30 Euros/min).
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the equipment warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the warranty.
- ◆ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Please do not forget filling the warranty certificate and send it before 15 days after the date of purchase.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

**ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

#### INSTRUCTIONS ON ENVIRONMENT PROTECTION

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

- ◆ When you are not using the mixer for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ◆ Never try to repair the stirrer by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the mixer (see Maintenance).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

## 5. MAINTENANCE AND CLEANING

To get the best results and a higher duration of this equipment it is essential to follow the processes of use.

**Note:** All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ The brushless motor is maintenance-free.
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- ◆ The mixer is supplied with a Schuko standard wire. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.
- ◆ Never try to repair it by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the mixer (burns, hurts...) or damages in nearby equipments.
- ◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.

### Cleaning

- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the mixer and produce an early ageing of the equipment
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

## INDEX OF LANGUAGES

Spanish .....	2-7
English .....	8-13
French.....	14-20

## INDEX OF CONTENTS

1. USES OF THE INSTRUMENT .....	9
2. DESCRIPTION .....	9
3. TECHNICAL SPECIFICATIONS .....	10
4. INSTALLATION / SETTING UP .....	10
5. MAINTENANCE AND CLEANING .....	12
ANNEX I: CE CERTIFICATE .....	20

## 1. USES OF THE INSTRUMENT


The vortex mixer for multiple tubes allows a fast and efficient mixing of solutions thanks to its vibratory movement. It is an ideal equipment for intermittent stirring of tubes of different sizes as well as small Erlenmeyers, beakers, flasks etc.

The three suction feet assures a good stability during operation.

## 2. DESCRIPTION

- 1.1. Stirring platform
- 1.2. Power on/off
- 1.3. Base with suction feet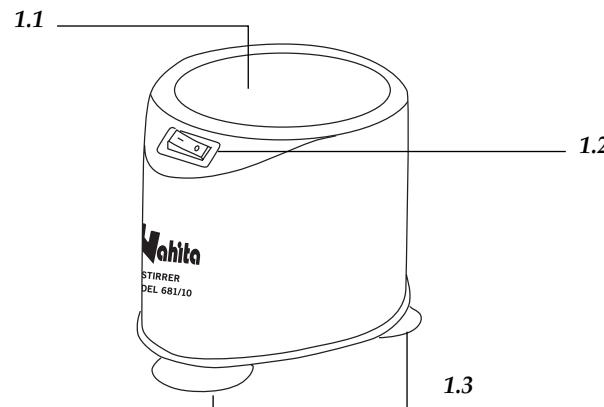

### 3. TECHNICAL SPECIFICATIONS

<b>Code</b>	50681010
<b>Movement</b>	Vibratory
<b>Stirring amplitude</b>	6.2 mm
<b>Stirring speed</b>	2600 rpm
<b>Useful platform diameter</b>	100 mm
<b>Consumption</b>	25 W
<b>Power</b>	220 V / 50 Hz
<b>Dimensions</b>	150x130x140 mm
<b>Weight</b>	2.3 Kg

### 4. INSTALLATION / SETTING UP


#### Preliminary inspection

Unwrap the vortex mixer, take off the involving plastic and take off the polyspan protection in which it comes fitted. Take off all protections and without connecting the mixer to the power supply, make sure that it does not present any damage because of the shipment. In case it presents any damage, tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.

Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.

Please check that all the accessories are enclosed with the equipment:

- 3 suction feet
- User's manual
- Warranty certificate

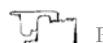We will only accept any equipment return within 15 days after delivery and provided it comes in its original wrapping.

#### *Installation*

Before using the vortex mixer, it is convenient for you to familiarize with its components and basic essentials, as well as with its control functions.

**TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.**

- ◆ Put the vortex mixer on top of a horizontal, plane and stable table making a free space at least at 30 cm per side. Do not put the equipment near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc.

- ◆ During operation, flammable liquids must be out of the working area.

- ◆ Screw the suction feet supplied into the corresponding holes at the base of the equipment. Put the mixer on the suction feet so as it stays perfectly stable.

Please insert the Schuko wire that feeds the AC electric current in the base of current 220 V, 50 Hz ± 10% provided with earth wire.

**Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to persons suffered by the improper use of the electric connection. The tension should be 220 V, 50 Hz ± 10%.**

#### *Setting up*

- ◆ Turn the vortex mixer on by pressing power on/off (1.2).
- ◆ Put the tube with the solution to be mixed on the stirring platform (1.1). **Note: tubes must be perfectly closed to avoid liquid spillages on the equipment.**
- ◆ Once solution is mixed turn the equipment off by pressing power on/off (1.2).

If you are not using the vortex mixer for a long period of time please make sure it is disconnected from the net and protected from dust (this way you will avoid accidents and will extend its working-life).

#### *Security*

- ◆ The vortex mixer for multiple tubes must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ◆ Put the mixer on top of a horizontal, plane and stable table making a free space at least at 30 cm per side.
- ◆ Do not put the mixer near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc. Avoid vibrations, dust and dry environments.
- ◆ During operation, flammable liquids must be out of the working area.