ANEXE I: CERTIFICAT CE

AUXILAB S.L.

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ANNEX I: CE CERTIFICATE

AUXILAB S.L.

CE DECLARATION OF CONFORMITY
OVERHEAD STIRRER NAHITA 686 of AUXILAB, S.L. for the Directive of Machines (89/392/CEE modified) and the regulations adopted for their transposition

NAME OF THE MANUFACTURER / IMPORTER:

AUXILAB, S.L.

ADDRESS:

**Polígono Morea Norte, 8
 31191 Beriáin (Navarra)**

WE STATE THAT:

**OVERHEAD STIRRER MODEL 686
 CODES 50686001 and 50686002**

Are designed and manufactured according to:

- ♦ Directive 89/392/CEE, including the modifications and the national regulations that transpose them.
 - ♦ Directive 73/23/CEE modified over the electric security.
 - ♦ Directive 89/336/CEE modified over the electromagnetic compatibility.
- And that the following harmonized rules have been applied (or part of them):
- UNE 292-1/-2/-2/A1, UNE-EN 1050, UNE-EN 614-1, UNE-EN 953, UNE-EN 294, UNE-EN 418, UNE-EN 894-1, UNE-EN 894-2, UNE-EN 954-1, UNE-EN 60555-2-3, UNE-EN 60204-1, UNE 50082-1, UNE 55014-1, UNE 61010-1/A2, UNE-EN61010-2-051.

BERIAIN 6th October 2005

Signed by: ALFONSO AINCIBURU SANZ
 DIRECTOR/MANAGER

Polígono Morea Norte, 8 31191 Beriain (Navarra) - Spain. Tel. 948 310 513 Fax 948 312 071
Internet: www.auxilab.es · Email: correo@auxilab.es

Cleaning

- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the overhead stirrers and produce an early ageing of the equipment.
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

ATTENTION!!
IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY
WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

6. TROUBLESHOOTING

PROBLEM	CAUSE	SOLUTION
The overhead stirrer doesn't work	Insure: - Insert well the wire. If you pressing the security switch in position " I ". - If you pressing  .	In the event of breakdown please contact your distributor.
The stirring isn't homogeneous	Insure: - The paddle stirrer is well place. - The paddle stirrer is centre and - The paddle stirrer isn't best or worn out.	- Read the instructions: assembly the paddle stirrer - Centre the paddle stirrer - Change the paddle stirrer

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Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita Overhead Stirrers 686. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user. The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of *Maintenance*, where we expose the security notes.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

Please bear in mind the following:

- ♦ This manual is inseparable from the Overhead stirrers 686, so it should be available for all the users of this equipment.
- ♦ You should carefully handle the overhead stirrers 686 avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ♦ Never dismantle the different pieces of the overhead stirrers 686 to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the overhead stirrers 686.
- ♦ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ♦ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to asistencia@auxilab.es or by telephone: +34 807 117 040 (0,30 Euros/min).
- ♦ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ♦ Overhaul is not covered by the overhead stirrers 686 warranty.
- ♦ Operations made by non-qualified staff will automatically produce a loss of the overhead stirrers 686 warranty.
- ♦ Neither fuses (1A) nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ♦ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ♦ Please do not forget filling the warranty certificate and send it before 15 days after the date of purchase.

ATTENTION!!

IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

INDEX OF LANGUAGES

Spanish	2-9
English	10-17
French	18-24

Security

- ♦ The overhead stirrers must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ♦ You should put the overhead stirrers in a horizontal plane stable table, having a safety area of at least 30 cm per side.
- ♦ Do not place the overhead stirrers near any warm supply (burners, blowlamps, etc), nor expose it directly to the sun. Avoid vibrations, dust and dry environments.
- ♦ During its functioning dangerous materials such as flammable or pathological substances must be out of the safety area.
- ♦ When you are not using the overhead stirrers for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ♦ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ♦ Never try to repair the overhead stirrers by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ♦ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the overhead stirrers (see Maintenance).
- ♦ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.


5. MAINTENANCE AND CLEANING

To get the best results and a higher duration of this equipment it is essential to follow the processes of use.

Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

- ♦ Please follow the processes of use of this manual.
- ♦ This manual should be available for all users of this equipment.
- ♦ Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- ♦ The overhead stirrers is supplied with a Schuko standard wire. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.
- ♦ Never try to repair the overhead stirrers by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the overhead stirrers (burns, hurts...) or damages in nearby equipments.
- ♦ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.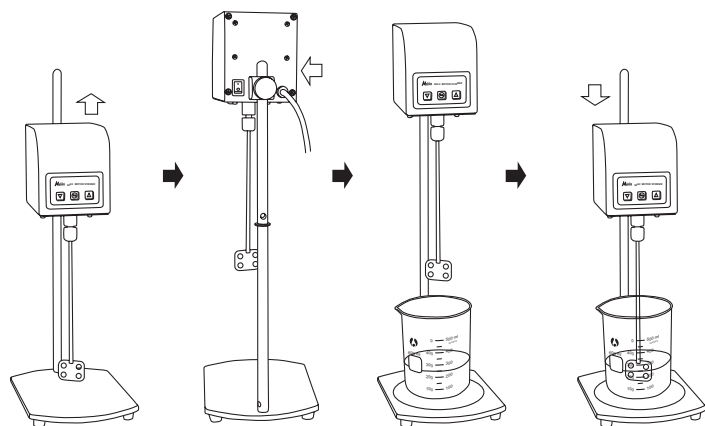

♦ Please insert the wire that feeds the AC electric current in the base of current 220V 50Hz $\pm 10\%$ provided with earth wire and to the other end to the overhead stirrer connector.

Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to persons suffered by the improper use of the electric connection. The tension should be 220V 50Hz $\pm 10\%$.

♦ If you are not using the overhead stirrers for a long period of time please make sure it is disconnected from the net and protected from dust (this way you will avoid accidents and will extend its working-life).

♦ Pressing the security switch 1.7 into position "I".

♦ Pressing  (1.8); the stirrer begins to work. the speed which it begins to turn the paddle stirrer is rpm average.

♦ Set the speed pressing   (1.9).

Note: In the model 686/2 sound on bip by pulse.

The range speed 200-1000 rpm in model 686/1 y 200-2000 r.p.m in model 686/2. If viscosity increases due to the process, the speed reduces gradually and viceversa. Please note that fluctuations in mains voltage within the permitted tolerances may also cause minor fluctuations in speed.

♦ When you want to stop, pressing  (1.8). If you want again to start pressing  (1.8); the rpm speed is the previous.

♦ If we the stirrer from the security switch or first we turn off from  (1.8) and soon pressing the security switch the speed which it begins to turn the paddle stirrer is rpm average.

♦ When you finish the work:

- Pressing the security switch 1.7 to the position "O". Unplug the stirrer of the network.
- Lift the stirrer until above of the plate stand
- Remove the paddle stirrer, We relaxed the screw we held the superior part and we turned with the other hand (Clockwise). Clearing and drying the paddle stirrer.


INDEX OF CONTENTS

1. USES OF THE INSTRUMENT	11
2. DESCRIPTION	11
3. TECHNICAL SPECIFICATIONS	12
4. INSTALLATION / SETTING UP	12
5. MAINTENANCE AND CLEANING	15
6. TROUBLESHOOTING	16
ANNEX I: CE CERTIFICATE	17

1. USES OF THE INSTRUMENT

The overhead stirrers Nahita is suitable for stirring and mixing fluids with low viscosity. It is designed for use in laboratories.

2. DESCRIPTION

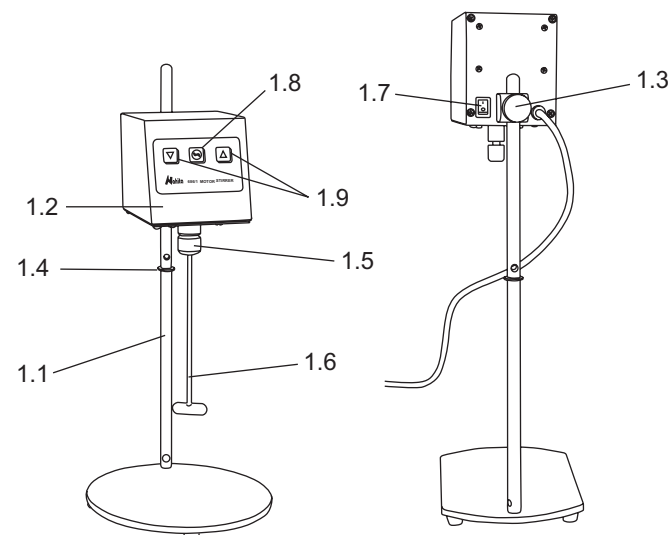- 1.1. Plate stand
- 1.2. Stirrer board
- 1.3. Thread or screw
- 1.4. Limit
- 1.5. Screw


- 1.6. Paddle stirrer
- 1.7. Security switch
- 1.8. Knob 
- 1.9. Increase/decrease Knob of speed


3. TECHNICAL SPECIFICATIONS

Code	50686001	50686002
Stirring quantity max. H ₂ O	5 L	10L
Speed range	200-1000 rpm	200-2000 rpm
Speed control	Digital	Digital
Extension arm diameter	4 mm	5 mm
Out power	20W	29 W
Input power	100 W	100W
Dimensions	16x16x37 cm	16x24x50 cm
Weight	2,75 Kg	5,50 Kg

686/1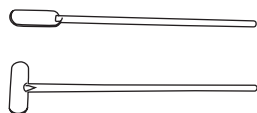

686/2

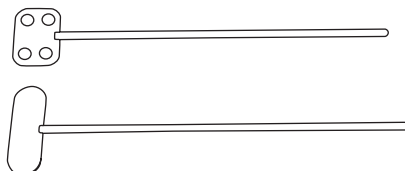Paddle stirrer	Shaft length	Stirrer diameter
Paddle stirrer (686/1)	160 mm	10 mm
Paddle stirrer (686/1)	160 mm	30 mm
Paddle stirrer (686/2)	240 mm	60 mm
Paddle stirrer with 4 holes (686/2)	240 mm	40 mm

4. INSTALLATION / SETTING UP

Preliminary inspection

- ◆ Unwrap the overhead stirrer, take off the involving plastic and take off the polispap protection in which it comes fitted. Take off all the protective items and, **without connecting the equipment to the net**, make sure that it does not present any damage because of the shipment. In case the overhead stirrer presents any damage tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.
- ◆ Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.
- ◆ Please check that all the accessories are enclosed with the equipment:
 - Paddle stirrer
 - Plate stand
 - Schuko cable
 - Anti skind base

We will not accept any equipment in return period unless it comes in its original wrapping.

Installation

- ◆ Before using this instrument, it is convenient for you to familiarize with its components and basic essentials.

PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT.




- ◆ Please put the overhead stirrer on top of a horizontal, plane and stable table making a free space at least at 30 cm per side. Do not put the overhead stirrers near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc.
- ◆ Avoid inflammable or toxic substances in the working area.
- ◆ The overhead stirrer is supplied with a Schuko standard wire.
- ◆ The stirrer 1.2 is fixing in the stand plate 1.1. It has a thread to fix the stirrer 1.3 The height to select will depend to the container use The rod-stand has limit 1.4, it is optimal height to the work. Depending the thickness of the base-container. In order to adapt the height of the rod you have to consider the characteristics to the container: thickness, curvature, etc. the paddle stirrer never must touch the bottom of the container.

Assembly of the paddle stirrer

- ◆ Relax screw 1.5 We help the superior part and turned with the other hand (clockwise)
- ◆ Insert the paddle stirrer 1.6 in the screw of the stirrer, to do this, press the extension arm until it catches in the hole please do not use force to do this. Relax the screw. Apretar the screw we help the superior part and turned with the other hand (opposite clockwise)
- ◆ Check that the extension arm is firmly seated.

Note: When you place the paddle stirrer, check that the extensión arm is firmly seated and centered, distribute the loads of stirring the uniform.

- ◆ Take up the stirrer until above of rod-stand 1.1.
- ◆ Put the container with the product that goes to stirrer.

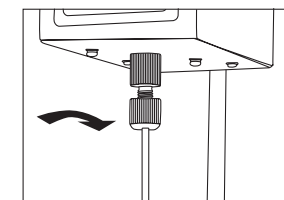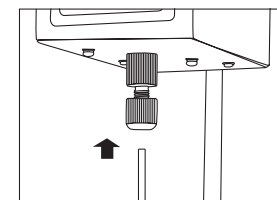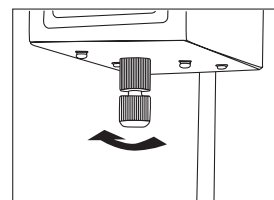- ◆ Take down the stirrer until height interest The rod-stand has limit 1.4, it is optimal height to the work. Depending the thickness of the base-container. In order to adapt the height of the rod you have to consider the characteristics to the container: thickness, curvature, etc. the paddle stirrer never must touch the bottom of the container.