DECLARACIÓN DE CONFORMIDAD CE
CE DECLARATION OF CONFORMITY
DÉCLARATION DE CONFORMITÉ CE

El fabricante | The manufacturer | Le fournisseur:

AUXILAB, S.L.

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Thank you for choosing this equipment. We sincerely wish that you enjoy your balance Nahita model 515X. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT

Please bear in mind the following:

- ◆ This manual is inseparable from the balance Nahita model 515X, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the balance avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the balance to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle it.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to asistencia@auxilab.es or by telephone: +34 807 117 040 (0.35 Euros/min).
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the balance warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the balance warranty
- ◆ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Please do not forget filling the warranty certificate and send it before 15 days after the date of purchase.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

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<i>Problem</i>	<i>Cause</i>	<i>Solution</i>
Value of weight which appears in the display is wrong	1 Balance is not calibrated 2 Balance has not been tared before weighing	1 Calibrate the balance 2 Tare the balance before each weigh
Balance is not calibrated	1 Calibration weight is not right for the balance 2 Zero is not adjusted	1 Use right weight 2 See Tare

INSTRUCTIONS ON ENVIRONMENT PROTECTION

Nahita digital balances serie 515X work with batteries; deposit batteries when exhausted in special containers or collection points. Batteries contain heavy metals (zinc, lithium, mercury, etc) if these metals are released into nature, they can contaminate water, land, etc..

Do not throw out this balance in the ordinary garbage; take it to a collection point for recycling electrical and electronic equipments. Does not contain dangerous or toxic elements to humans.

Materials are recyclable. Recycling is an important contribution to environmental protection.

Please contact the administration of your community for be advised on the collection points.

ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

INDEX OF LANGUAGES

1. Spanish	2-10
2. English.....	11-19
3. French.....	20-28

INDEX OF CONTENTS

1. USES OF THE INSTRUMENT	12
2. DESCRIPTION	12
3. TECHNICAL SPECIFICATIONS	13
4. INSTALLATION / SETTING UP.....	14
5. BALANCE FUNCTIONS	15
6. MAINTENANCE AND CLEANING	17
7. TROUBLESHOOTING.....	18
ANNEX I: CE CERTIFICATE	29


1. USES OF THE INSTRUMENT

Nahita electronical balances are a basic instrument in any laboratory to carry out weigh process. Developed with the latest technology, great reliability and durability , its solid, clean and ergonomic design is very suitable for industrial use.

2. DESCRIPTION

1. Plate
2. Digital display
3. ON / OFF
4. UNIT: selecting weigh units
5. COUNT: count pieces
6. TARE: tare button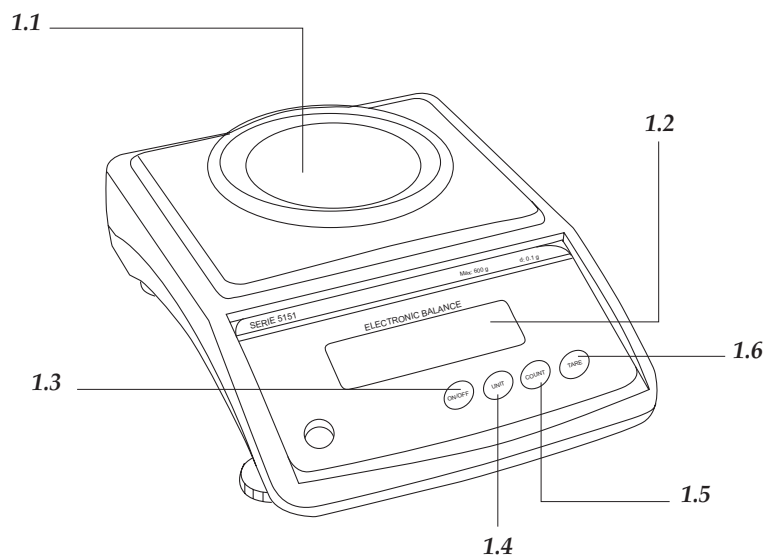

3. TECHNICAL SPECIFICATIONS

We emphasize the following features of electronic balance Nahita serie 515X:

- ◆ Fast start-up and stabilization (gauge “o” on display)
- ◆ Easy use, clear reading LCD screen.
- ◆ Features tare, parts counting and calibration
- ◆ Selecting units of weigh: grams (g), grains (GN), carats (ct), pennyweights (dwt), pounds (lb), ounces (oz) and American ounce (ozt)
- ◆ Bubble level and threaded legs for perfect balance of the equipment.
- ◆ Keyboard waterproof resistant to falling liquid.
- ◆ Connected to the network by adapter (optional) or battery operation (not included).

Code	55151260	55152230
Model	5151	5152
Capacity	600 g	300 g
Sensibility	0.1 g	0.1 g
Tara Range	0-600 g	0-300 g
Plate diameter	115mm	
Power	Batteries (6x 1.5 V AA) or Adapter 220V (optional)	

- ◆ To clean the equipment we recommend using a fluff-free cloth, moistened with soapy water. Do not use abrasive products.

ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

NOTE: According to present legislation for “Not automatic weight instruments” (balances included), by Order December 22th, 1994 (B.O.E. March 1st 1995) these balance can not be used for:

- Carrying out business transactions
- Calculating rates, tariffs, taxes, remunerations, compensations and other kind of similar canons.
- Experts judicial reports.
- Pharmaceutical preparations, ordered medicines and analysis carried out in medical and pharmaceutical laboratories.
- etermination of price or total amount in direct sales and preparation of pre-packed.

7. TROUBLESHOOTING

<i>Problem</i>	<i>Cause</i>	<i>Solution</i>
Display does not show digits	1 Adapter is not connected 2 Exhausted batteries	1 Connect the adapter to the power supply 2 Change batteries
Display does not show values	1 Low battery	1 Change batteries
Value of display is unstable	1 Adapter broken	1 Check operation with batteries, if it works, change the adapter. If the problem goes go, contact with alter sales service

main supply to prevent possible accidents.

- ◆ For cleaning, verification or replace any component (ex: replace fuse) it is obligated to turn off the equipment and disconnect it to the power supply.
- ◆ Do not try to repair by yourself; you will lose warranty and may cause damage in the overall functioning of the equipment, personal damages (burns, injuries...) and damages to the wiring.
- ◆ Avoid getting water on the table controls, even if it is adequately insulated. If for any reason you suspect that water or other liquid have come inside, quickly switch off the equipment (check maintenance).

6. MAINTENANCE AND CLEANING

For a proper functioning of the balance is necessary to follow certain recommendations.

Note: All rules for use have no value unless there is a continuous maintenance work

- ◆ Follow the instructions and warnings indicated in this manual.
- ◆ Keep this manual for consulting.
- ◆ Avoid abrupt movements, knocks, direct sunlight or air currents. Handle the balance carefully, like a precision instrument.
- ◆ Optionally, an adaptor is available. Adapter must be connected to a power supply ground wired to disconnect it in an emergency.
- ◆ Always switch off the adapter by pulling its base, never pulling the wire.
- ◆ Leave a free space around the balance for proper ventilation. Never cram the balance (ex: into a shelf).
- ◆ Do not use sharp objects such as pens, etc, to press buttons of front panel, use your fingers.
- ◆ Do not place on weigh plate an object heavier than range of balance, sensor can be damaged.
- ◆ Do not immerse or throw liquid on the balance.
- ◆ If it works with batteries, remove them when the balance is not going to be used for a long time.
- ◆ If for any reason liquid comes in contact with electrical parts, turn off and switch off the balance immediately. Send the equipment to the after sales service as soon as possible for its review and adjustment.
- ◆ Always use original spare parts and components. Other devices can be similar, but they can damage the equipment.

Cleaning

- ◆ To clean metal parts, stainless steel, aluminium, paints, etc never use scourers or products that may scratch the surface, because they deteriorate the balance.

4. INSTALLATION / STARTING-UP

Preliminary inspection

Unpack the balance, remove the plastic and the porexpan protection. Remove all protections and don't connect the balance to the power supply. Check there is no transport damage, if there is, please contact as soon as possible with the carrier or your distributor so he can make the appropriate complaint.

Keep the packaging, because if it is necessary returns this must be made in its original packaging with all supplied accessories.

Check you have received all the accessories:

- weights calibration (serie 5152)
- windbreaks (serie 5152)
- supply adapter
- owner's manual
- warranty

We only accept refunds of equipments within the first fifteen days after receiving and whenever they come complete in their original packaging.

Installation

Before using the equipment, we recommend you, review components as well as the functions of its controls.

READ THIS MANUAL CAREFULLY BEFORE OPERATING TO GET THE MAXIMUM PERFORMANCE AND LONGER DURATION

1. Put the balance on an horizontal table, flat and stable, making a free space of at least 30 cm by each side. Do not put the equipment near of heat sources (burners, torches, etc), and do not expose directly to sunlight.
 2. Level the balance thanks to the threaded legs and center level bubble correctly.
 3. Nahita electronic balances serie 515X can work with batteries or switch to the power supply with an adapter CA.
 4. Place the AC adapter.
- Insert the switch of the adaptor in the holes that you will find in the back of the balance and connect to the power supply with grounded wire.

Neither the manufacturer nor the distributor assume any responsibility for damages in the equip, installations or injuries to persons due to the failure of the proper procedure for electrical connection. Voltage must be 220V 50Hz+10%

5. Place the batteries
- ◆ Put the balance upside down, remove the cap battery compartment and loosen the

screws.

- ◆ Use six alkaline batteries 1.5V type AA, be careful with the polarity (indicated inside).
- ◆ Replace the battery cover fixing the screws.

Getting started

- ◆ Press button ON/OFF (1.3) for several seconds to switch the balance.
- ◆ When display indicates “0.0”, the balance is ready to weigh.
- ◆ Press button UNIT “(1.4) to choose weigh unit you desire.
- ◆ Press button ON/OFF (1.3) to turn off the balance.

Abbreviation	Name	Conversion
Kg	Kilogram	1000 g
g	Gram	1 g
mg	Milligramme	0.001 g
GN	Grain	0.064799 g
ct	Carat	0.1999694 g
dwt	Pennyweights	1.55517 g
lb	Libras	453.592 g
oz	Ounce	28.3495231 g
ozt	American ounce	31.1034 g

5. BALANCE FEATURES

Tare feature

Tare is used when you want to weigh substances but without the container, for example when you want to weigh liquids or solid reagents that after they are going to dissolve.

- ◆ Put the empty container on the plate and press button Tare (1.6), display will indicate “0.0”. when you the symbol of stabilized “0”, balance is tared and ready to weigh the substance.

NOTE: Balance presents a range of tare equal to the maximum capacity of the equipment; take into account that maximum because if we use the maximum capacity tare will not be able to weigh anything else. (Example: if we carry out a tare of 450 g in a balance of capacity maximum 500 g, we will be able to weigh only 50 g more).

Calibration feature

Before using the balance and when significant changes are observed in wrong weigh or weighs we should calibrate the balance. Factors such as variation of the gravity depending on latitude, changing workplace or a rude and inappropriate used, indicate that the cali-

bration must be carry out.

Important! Not move the balance during calibration

Turn on the balance pressing ON/OFF (1.3) and be sure there is nothing on the plate (1.1).

Press button Tare (1.6) and when display indicates “0.0g” press button UNIT (1.4) for three seconds. Display will blink indicating CAL-0 and then 500.0g (Serie 5152:200.00 g). Place calibration weight on the balance.

When stabilized symbol (“o”) appears in the display and the value indicated is right, calibration is ended (if the value is not right, repeat the calibration process).

Code	Calibration weights
55151260	500 g (not included)
55152230	200 g (included)

Before calibration the balance can be used normally

Counting part feature (“Mode COUNTING”)

In mode “counting” the balance calculate number of pieces in the plate. To do this proceed as follows:

Press ON/OFF (1.3) to turn on the balance and wait until value “0.0g” stabilizes in the display.

Place reference samples in the balance and press COUNT (1.5)

Screen displays default value “10” that indicates number of pieces assigned to the sample.

If you want to assigned others reference values press successively button COUNT. Display will show consecutively the values 10, 20, 50 and 100.

Remove the sample (value “O” in display) and place the samples for counting.

Display will show the corresponding value .

Safety

Balance should be used by qualified personal. This personal should know the equipment and how to use it thanks to the manual.

Place the balance on an horizontal, flat and stable table, creating a free space of at least 30 cm by each side.

Do not place the balance near heat sources (burners, blowtorch...), do not expose the equipment directly to sunlight. Avoid vibrations, dust and dry environments.

When equipment is not going to be used for a long time, be sure it is disconnected from