DECLARACIÓN DE CONFORMIDAD CE
CE DECLARATION OF CONFORMITY
DÉCLARATION DE CONFORMITÉ CE

El fabricante | The manufacturer | Le fournisseur:

AUXILAB, S.L.

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This manual should be available for all users of these equipments. To get the best results and a higher duration of this equipment it is advisable to read carefully this manual and follow the processes of use.

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
ANNEX I: WIRING DIAGRAM

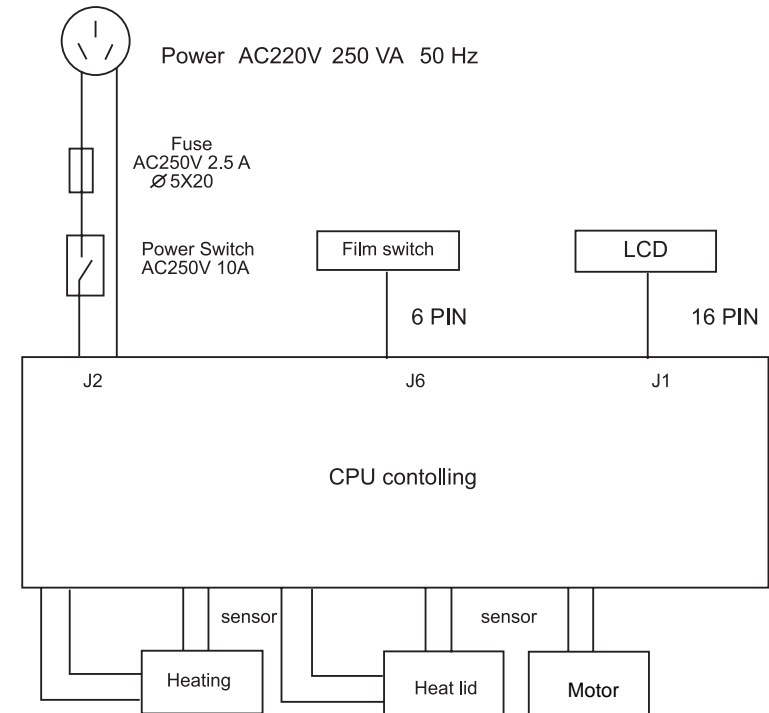(This diagram is just for you reference. Manufacturer reserves the right to modify or improve it).

◆ Never try to repair the equipment by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the thermoshaker (burns, hurts...) or damages in nearby equipments.

◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.

Cleaning

◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the equipment and produce an early ageing of the equipment.

◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

INSTRUCTIONS ON ENVIRONMENT PROTECTION

Do not dispose of this product in the usual household garbage at the end of it's life cycle; hand it over at a collection point for the recycling of electrical and electronic appliances. The symbol on the product, the instructions for use or the packing will inform about the methods for disposal.

The materials are recyclable as mentioned in it's marking. By recycling, material recycling or other forms or re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

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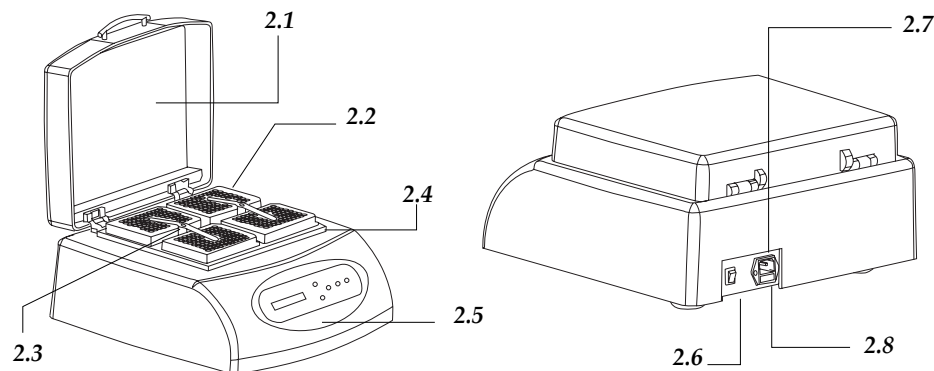
Fixing of the microtest plates (only 603/104 model).

Turns on the cover, the two spring patch according to anticlockwise revolves to the highest position, moves the spring patch the level location. Put the microtest plates in the heating plate. Rotates spring patch 90°, put them on the top of microtest plates. Rotating clockwise direction. Tightens the knob around, contracts the spring patch to suppress the microtest plates. The microtest plates do not skid up to.

Security

- ◆ The thermoshaker must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ◆ You should put the equipment in a horizontal plane stable table, having a safety area of at least 30 cm per side.
- ◆ Do not place the thermoshaker near any warm supply (burners, blowlamps, etc), nor expose it directly to the sun. Avoid vibrations, dust and dry environments.
- ◆ During its functioning dangerous materials such as flammable or pathological substances must be out of the safety area.
- ◆ When you are not using the equipment for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ◆ Never try to repair the thermoshaker by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the equipment (see Maintenance).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

5. MAINTENANCE AND CLEANING

To get the best results and a higher duration of this equipment it is essential to follow the processes of use.

Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.

The thermoshaker is supplied with a Schuko standard wire. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.

The temperature of "P:" and "AdjTemp" is current temperature 20.5 °C. When the temperature reaches 40 °C, the temperature of "P" is current temperature of block while the temperature of "AdjTemp" is still 40 °C with cursor. The "Adj" and "*" is flickering.

P: 40.1	Adj*
AdjTemp=	40.0C

◆ After 20 min, the actual temperature of thermometer is 38.8 °C. Press SET to make the cursor rest on the digital you want to change. Press increasing or decreasing button to adjust and press START / STOP to confirm.

P: 40.1	Adj*
AdjTemp=	38.8C

The system will save the data and increase to 100 °C automatically. The cursor will disappear as showed in chart

P: 40.1	Adj*
AdjTemp=	40.1C

◆ When the system temperature reaches 100 °C, the temperature of P is actual temperature and the temperature of "AdjTemp" is 100 °C with the cursor. The "Adj" and "*" is flickering.

P: 100.1	Adj*
AdjTemp=	100.0C

◆ After 20 min, the actual temperature of thermometer is 98.0 °C. Press SET to make the cursor rest on the digital you want to change. Press increasing or decreasing to adjust.

P: 100.1	Adj*
AdjTemp=	98.0C

◆ After the finish of adjustment, press increasing button and SET at the same time and keep off at once. After 8s the cursor display and the system accepts the adjustment. The screen returns to the display as the chart.

P: 98.5	1000	00:00
S: 100.0	1000	00:00

After the adjustment, the temperature displayed is the same with the actual temperature.

P: 100.0	1000	00:00
S: 100.0	1000	00:00

Note: During the adjustment, press the increasing and decreasing buttons contemporary to cancel the adjustment. The system keeps the last adjustment, so don't press increasing button and SET at the same casually.

Exchange of the metal block (only 603/100 model)

Open the transparent lid and pull out the four screws which fix the block to the heating board with the screwdriver, close the lid, and put out the block from the main engine. Take another model of block, steadily lays aside on the main engine. Put in the instalment holes the screws, and fix with the spanner the metal block on the instrument.

The following blocks are available:

Code	Capacity
90603021	96 wells of 0.2 mL
90603022	54 x 0.5 mL
90603023	35 x 1.5 mL
90603024	35 x 2.0 mL

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◆ The operation runs, long time press START / STOP, and the operation stops. The temperature will be decreasing, the speed displays 0, the time displays 00:00. Presse again, the operation runs again.

P: 55.5	0	00:00
S: 60.0	1000	20:00

3. Instantaneous mixing functions

◆ Press SET and keep off at once, after the "S:" there is a cursor "_", press SET to remove the cursor, and remove the cursor under the time column, press increasing or decreasing button, the time function can be changed 00:00.

P: 55.5	0	00:00
S: 60.0	1000	20:00

◆ Press START / STOP and keep off at once, the instrument will be mixing 6s and then stop.

P: 59.5	0	00:00
S: 60.0	1000	20:00

Temperature error adjustment (only 603/100 model)

The temperature of the Instrument has been adjusted before it is sold out. But if there is deviation between the actual temperature and the displayed temperature due to some reasons, you can do as follows to correct the error.

Notes: The Instrument uses double temperatures adjustment to ensure its veracity. This means it is linearly adjusted on 40 °C and 100 °C two points. The temperature veracity will be within ± 0.5 °C after the double temperature adjustment.

Both circumstances and the block temperature should be lower than 35 °C.

Adjustment methods as follows:

◆ After the start up of the Instrument, it enters waiting interface. Make sure the temperature in display is below 35 °C. If the temperature is higher than 35 °C, you should wait until the temperature is below of 35 °C.

◆ Inject olefin oil into one of the cone-shaped wells, and then put a thermometer into this well (Make sure the precision of the thermometer should be within 0.1 °C and the temperature ball should be absolutely immersed into the cone-shaped well). Heat insulation material is needed on the block to separate it from the circumstance.

Notes: Please read the actual value after 20minutes' constant temperature to ensure the adjustment precision.

◆ Press increasing button and SET button contemporary to enter the temperature adjustment interface as in the drawing. The temperature begins to rise to 40.0 °C automatically and "*" is flickering in the right corner.

P: 20.5C	Adj*
AdjTemp=	20.5 C

The equipment is supplied with a Schuko standard wire

Please insert the wire that feeds the AC electric current in the base of current 220 V, 50 Hz ± 10% provided with earth wire and to the other end to the thermoshaker connector.

Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to persons suffered by the improper use of the electric connection. The tension should be 220 V, 50 Hz ± 10%.

If you are not using the equipment for a long period of time please make sure it is disconnected from the net and protected from dust (this way you will avoid accidents and will extend its working-life).

Setting up

1. Temperature, speed and time set

◆ The LCD will display the picture as the chart when the instrument powers on and the instrument goes into the initial state with the sound of "du...";

System-Testing		
--------------------------------	--	--

◆ About 6s later, "P:" line will display: 25.5 is the heating block's current temperature; 0 is the current speed; 00:00 is the current run time. "S:" line will display: 60.0 is the former set temperature; 1000 is the former set speed; 20:00 is the former set time.

P: 25.5	0	00:00
S: 60.0	1000	20:00

◆ Press the button of SET and keep off at once, after the "S:" there is a cursor "_", press SET to remove the cursor, and press increasing or decreasing button to change the digital on the cursor into the one required.

◆ If will shut off temperature function, press SET and the the cursor "_" will be removed the temperature column, and press OFF temperature column will be displayed "OFF". It's the same as shut off the speed and time function.

P: 25.5	0	00:00
S: OFF	1000	20:00

Notes: 1) Time function set "OFF", the time is without end.

2) Temperature, speed and time can't be set simultaneously "OFF"

3) The instrument starts to heat and according to the previous set temperature after the power on.

◆ After pressing "start" and keeping off at once, the operation runs as the set.

2. Temperature, speed and time set

◆ Press START / STOP and keeping off at once, the operation runs as the set.

P: 25.5	0	00:00
S: 60.0	1000	20:00

Note: When the operation is finished, the buzzer alarms.

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4. INSTALLATION SETTING UP

Preliminary inspection

Unwrap the equipment, take off the involving plastic and take off the polispán protection in which it comes fitted. Take off all the protective items and, without connecting the equipment to the net, make sure that it does not present any damage because of the shipment. In case the thermoshaker presents any damage tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.

Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.

Please check that all the accessories are enclosed with the equipment:

- Net wire
- Instruction manual
- Warranty certificate

We will only accept any equipment return within 15 days after delivery and provided it comes in its original wrapping.

Installation

Before using this instrument, it is convenient for you to familiarize with its components and basic essentials.

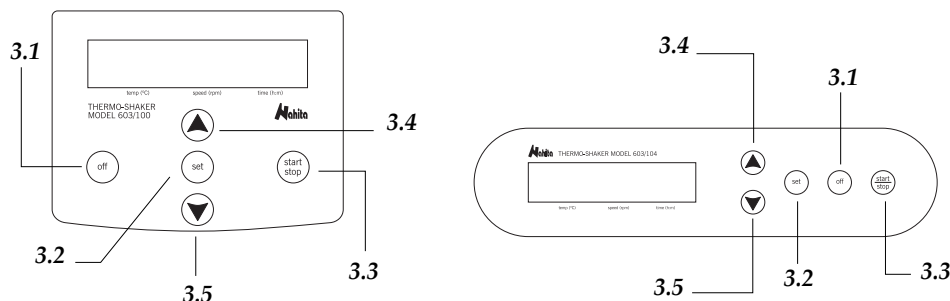
PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT

◆ Please put the thermoshaker on top of a horizontal, plane and stable table making a free space at least at 30 cm per side. Do not put the thermoshaker near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc.

◆ Avoid inflammable or toxic substances in the working area.

Operation panel and display diagram:


Button instruction:

- 3.1 OFF:** Shut or open temperature and speed and time function
- 3.2 SET:** Set button: used together with increasing and decreasing button to set the operation temperature, speed and time required
- 3.3 START / STOP:** Run / stop button, the long time presses this key to stop when the instrument runs.
- 3.4 ▲ :** Increasing button of the temperature or speed or time
- 3.5 ▼ :** Decreasing button of the temperature or speed or time

Display diagram

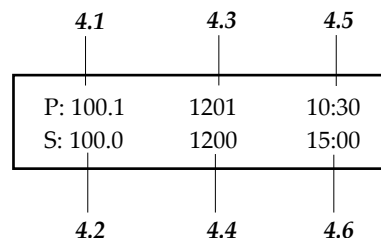- 4.1 Display of the current temperature
- 4.2 Display of the set temperature
- 4.3 Display of the current speed
- 4.4 Display of the set speed
- 4.5 Display of the run time
- 4.6 Display of the set time

3. TECHNICAL SPECIFICATIONS

Normal working condition:

The using environmental temperature: 5 C - 30 C
 The relative humidity: ≤ 70%
 The using power: AC220V 22V 250VA 50Hz 1Hz

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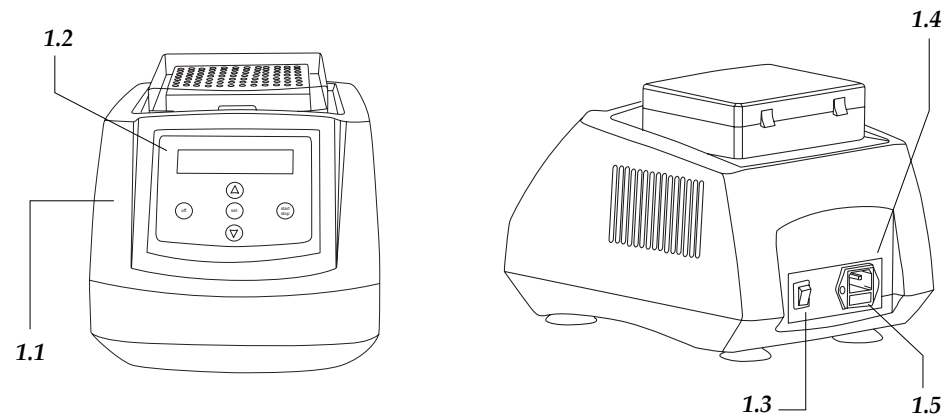
2. DESCRIPTION


Modelo 603/100

- 1.1 Block
- 1.2 Operation panel
- 1.3 Power switch
- 1.4 Power socket
- 1.5 Fuse socket

Modelo 603/104

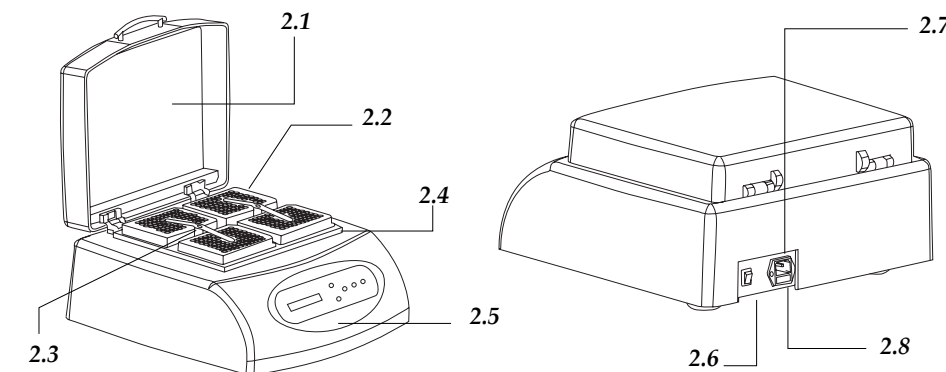- 2.1 Hot lid
- 2.2 Microtest plates
- 2.3 Pressure flake
- 2.4 Heating plate
- 2.5 Operation panel
- 2.6 Power switch
- 2.7 Power socket
- 2.8 Fuse socket

 **ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

INDEX OF LANGUAGES

Spanish 2-12
 English 13-23
 CE certificate 22

INDEX OF CONTENTS

1. USES OF THE INSTRUMENT 14
 2. DESCRIPTION 15
 3. TECHNICAL SPECIFICATIONS 16
 4. INSTALLATION / SETTING UP..... 17
 5. MAINTENANCE AND CLEANING 21
 ANNEX I: WIRING DIAGRAM 23

1. USES OF THE INSTRUMENT

The Thermo-shaker is an adequate instrument for intensive mixing of samples in the regulated temperature conditions.

Mixing and heating modes can be used both simultaneously and independently; the device can work as shaker and as a thermostat. The main body of the Mixing Block can be used with different kinds of blocks. It is applicable for DNA analysis, DNA library creation, PCR amplification, pre-denaturation in electrophoresis, etc.

Features of this product as follows:

1. Microprocessor controlled time, speed and temperature.
2. Simultaneous display of set and actual time, temperature and speed.
3. Beep-signal / Stop after program completion.
4. External power supply AC 220V, 2.5 A.

Model 603/100: The main body of the Mixing Block can be used with different kinds of blocks. The equipment is supplied without blocks. As it is equipped with various optional mixing blocks, the instrument can adapt to different tubes & wells to cope with experimental needs. It is easy to replace the metal blocks and is very simple to clean and sterilize. Customized blocks are available to suit your specific demands. Furthermore, temperature can be calibrated to meet user's needs.

Model 603/104: it is designed for shaking from 1 to 4 standard 96 well microplates.

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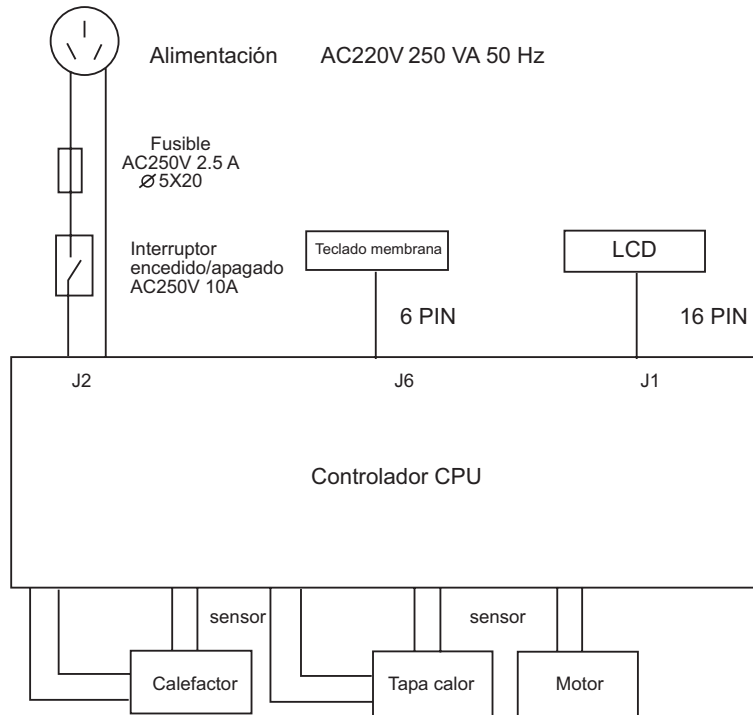
Thank you for choosing this equipment. We sincerely wish that you enjoy your thermoshaker Nahita. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

Please bear in mind the following:

- ◆ This manual is inseparable from the thermoshaker Nahita, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the equipment avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the thermoshaker to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the equipment.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to asistencia@auxilab.es or by telephone: +34 807 117 040 (0.30 Euros/min).
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the thermoshaker warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the equipment warranty.
- ◆ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Please do not forget filling the warranty certificate and send it before 15 days after the date of purchase.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.