**AUXILAB S.L.**

**CE DECLARATION OF CONFORMITY**  
**NAHITA CENTRIFUGES of AUXILAB, S.L. for the Directive of Machines (89/392/CEE modified) and the regulations adopted for their transposition**

NAME OF THE MANUFACTURER/IMPORTER:  
**AUXILAB, S.L.**

ADDRESS: **Polígono Morea Norte, 8**  
**31191 Beriáin (Navarra)**

WE STATE THAT:

**NAHITA CENTRIFUGES SERIES 2690**  
**CODES 52690000 AND 52690050**

Are designed and manufactured according to:

- ◆ Directive 89/392/CEE, including the modifications and the national regulations that transpose them.
  - ◆ Directive 89/336/CEE modified over the electromagnetic compatibility.
  - ◆ Directive 73/23/CEE modified over the electric security.
- And that the following harmonized rules have been applied (or part of them):

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BERIAIN 7th July 2004

Signed by: ALFONSO AINCIBURU SANZ  
 DIRECTOR/MANAGER

Polígono Morea Norte, 8 31191 Beriain (Navarra) - Spain. Tel. 948 310 513 Fax 948 312 071  
Internet: www.auxilab.es · Email: correo@auxilab.es

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*This manual should be available for all users of these equipments. To get the best results and a higher duration of this equipment it is advisable to read carefully this manual and follow the processes of use.*

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BERIAIN 1st April 2004

Signed by: ALFONSO AINCIBURU SANZ  
DIRECTOR/MANAGER

Polígono Morea Norte, 8 31191 Beriain (Navarra) - Spain. Tel. 948 310 513 Fax 948 312 071

Internet: www.auxilab.es · Email: correo@auxilab.es

## ANNEX I: CE CERTIFICATES

AUXILAB S.L.

## CE DECLARATION OF CONFORMITY

**NAHITA CENTRIFUGES of AUXILAB, S.L. for the Directive of Machines (89/392/CEE modified) and the regulations adopted for their transposition**

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- ◆ Directive 89/392/CEE, including the modifications and the national regulations that transpose them.
- ◆ Directive 89/336/CEE modified over the electromagnetic compatibility.
- ◆ Directive 73/23/CEE modified over the electric security.

And that the following harmonized rules have been applied (or part of them):

UNE-EN 1050, UNE-EN 292-1, UNE-EN 292-2, UNE-EN 292-2/11, UNE-EN 953, UNE-EN 894-1, UNE-EN 894-2, UNE-EN 894-2, UNE-EN 981, UNE-EN 60204-1.

BERIAIN 12th January 2004

Signed by: ALFONSO AINCIBURU SANZ  
DIRECTOR/MANAGER

Polígono Morea Norte, 8 31191 Beriain (Navarra) - Spain. Tel. 948 310 513 Fax 948 312 071  
Internet: www.auxilab.es · Email: correo@auxilab.es

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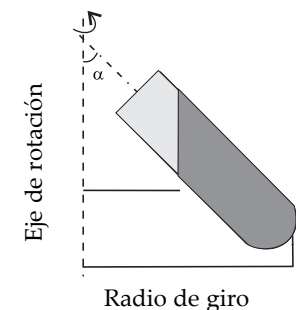
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- ◆ The analogic centrifuges is supplied with a Schuko standard wire and it should be connected to a current wire provided with an earth wire, it should be handy to be disconnected in case of emergency.
- ◆ Never try to repair the analogic centrifuges by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the analogic centrifuges (burns, hurts...) or damages in nearby equipments.
- ◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.
- ◆ Use tubes that are resistant to the R.C.F, check that at the bottom of the adapter there are some shock absorbers that are used as anti-shock.
- ◆ Balance the centrifuge placing tubes that are equal in weight, shape and size in symmetrical positions. If it is necessary, we will use tubes filled with water when one of the holes is empty. In a centrifuge it is very important the concept of weight, then it is advisable to have a balance to adjust the weights of the different tubes.
- ◆ This is absolutely necessary since if the rotor is not well balanced it can produce shakings, if these shakings are small it is difficult to notice them, then little by little the centrifuge wears away and then the breaking of the tube holders increases and it produces a worse sedimentation. It is advisable to change all the tube holders at the same time.
- ◆ Do not stop the centrifuge by the application of any external force, in this way you will avoid the possible mixture of the separate parts.
- ◆ Fuse change: According model remove the cover which is in blackboard.

- ◆ In case of damage or the change of the brushes, you have to get in touch with your supplier for the repair by the Nahita Technical Service.
- ◆ Change the carbon brush of the motor has been used for accumulative time of 1000 hours it should be replace.
- ◆ When revolving speed exceeds 4500 r.p.m, use nylon centrifugal tubes instead of glass tubes to avoid the possible break-up of glass.


#### Cleaning

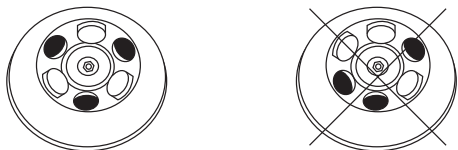- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the analogic centrifuges and produce an early ageing of the equipment.
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.
- ◆ With high g-values, there is a possibility that tube breakage will occur. Should this happen, the centrifuge, rotor, buckets, adapters and the rotor chamber must be thoroughly cleaned and all broken particles removed immediately.

**ATTENTION!!**  
**IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY**  
**WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

### Power

- ◆ Turn on the switch (1) and the light is lit to indicate that the power is connected.(Only on models 2650, 2698)

Adjust the speed and time with knob respective: speed knob (2) and time knob (3).

### Starting the centrifuge

- ◆ When adjust the time to start the centrifuge. If you want to continuous run turn the knob in opposite clockwise to the stop limit.(according to the model)

### Stop

- ◆ Once it past time the rotor stop. If you want to interrupt a centrifuge run, turn the knob timer (3) in clockwise.

### Security

- ◆ The centrifuge must not be operated by unqualified persons not familiar with the correct use and intended purpose of the machine.
- ◆ You should put the Schuko standard wire in a horizontal plane stable table, having a safety area of at least 30 cm per side.
- ◆ Do not place the analogic centrifuges near any warm supply (burners, blowlamps, etc), nor expose it directly to the sun. Avoid vibrations, dust and dry environments.
- ◆ During its functioning dangerous materials such as flammable or pathological substances must be out of the safety area.
- ◆ When you are not using the analogic centrifuges for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ◆ **It is essential to have the equipment switched off and unplugged from the net** before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ◆ Never try to repair the analogic centrifuges by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

## 6. MAINTENANCE AND CLEANING

*To get the best results and a higher duration of this equipment it is essential to follow the processes of use.*

Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.

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## 5. INSTALLATION / SETTING UP

### Preliminary inspection

- ◆ Unwrap the analogical centrifuges, take off the involving plastic and take off the polistan protection in which it comes fitted. Take off all the protective items and, without connecting the analogical centrifuges to the net, make sure that it does not present any damage because of the shipment. In case the analogical centrifuges present any damage tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.
- ◆ Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.
- ◆ Please check that all the accessories are enclosed with the equipment:
  - Tube holders
  - Schuko standard wire
  - Instruction manual
  - Warranty

*We will not accept any equipment in return period unless it comes in its original wrapping.*

### Installation

- ◆ Before using this instrument, it is convenient for you to familiarize with its components and basic essentials.

**PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT.**

- ◆ Please put the analogical centrifuges on top of a horizontal, plane and stable table making a free space at least at 30 cm per side. Do not put the analogical centrifuges near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc.
- ◆ Avoid inflammable or toxic substances in the working area.
- ◆ The analogical centrifuges is supplied with a Schuko standard wire.

Please insert the wire that feeds the AC electric current in the base of current 220V 50Hz  $\pm 10\%$  provided with earth wire and to the other end to the analogical centrifuges connector.

*Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to persons suffered by the improper use of the electric connection. The tension should be 220V 50Hz  $\pm 10\%$ .*

If you are not using the analogical centrifuge for a long period of time please make sure it is disconnected from the net and protected from dust (this way you will avoid accidents and will extend its working-life).

### Setting up

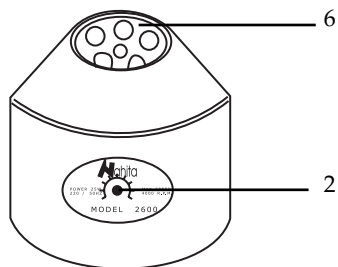
#### Cover's opening

(According to the model) press the button which is in right-lateral and open the cover. Put the test sample into centrifuge tubes carefully.


**IMPORTANT:** Make sure that the samples placed in the symmetrical tubes are of equal weight to avoid serious shaking due to big difference in weight, close the cover.

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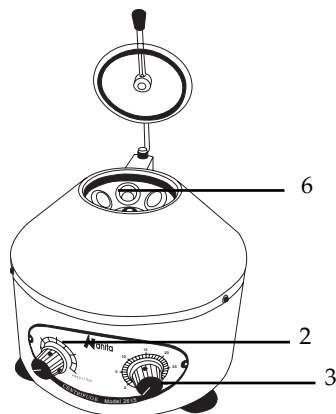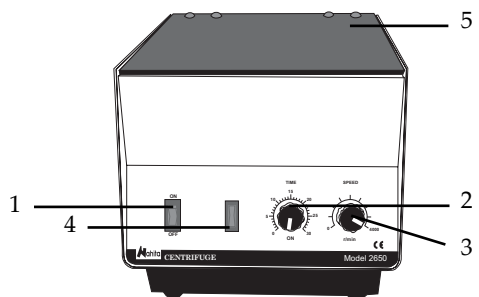
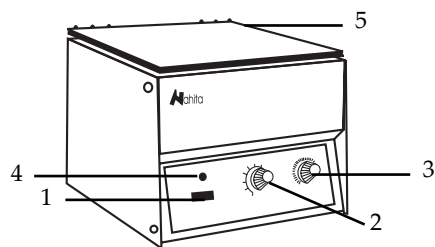
**1. USES OF THE INSTRUMENT**

Compact centrifuges, they sturdy in design and can be used for a variety of applications. Models with mechanical regulation have head for tubes of different volumes and with system of elastic suspension.

There are some examples of different activities in which this centrifuge is used as a technique of separation:

- Separation of serum or plasma.
- Cell concentration.
- Separation of substances.
- Biochemical laboratories.

**2. BASIC THEORY**

The centrifuge is an instrument that provides a technique of separation that is based on the movement of the particles. In this way the particles are moved towards the end of the rotating axis depending on their different masses and shapes.

The *Relative Centrifuge Force* (R.C.F.) is the required force so that the separation can be carried out. The units of this force are expressed in number of times higher than the gravity (g).

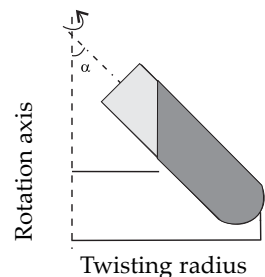
The R.C.F. is calculated by means of the following formula:

$$R.C.F. = 1,118 * 10^{-5} * r * n^2$$

Being  $1,118 * 10^{-5}$  a constant.

r = horizontal distance of the radius (in centimetres) from the centre of rotation up to the bottom of the tube.

n = speed of rotation expressed in revolutions per minutes (r.p.m.).

Thank you for choosing this equipment. We sincerely wish that you enjoy your analogical Nahita centrifuges. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of *Maintenance*, where we expose the security notes.

**TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.**

Please bear in mind the following:

- ◆ This manual is inseparable from the analogical centrifuges Nahita, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the analogical centrifuges avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the analogical centrifuges to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the analogical centrifuges.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to [asistencia@auxilab.es](mailto:asistencia@auxilab.es) or by telephone: +34 807 117 040 (0,30 Euros/min).
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the analogical centrifuges warranty.

Operations made by non-qualified staff will automatically produce a loss of the analogical centrifuges warranty.

Neither fuses (1A) nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.

- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Please do not forget filling the warranty certificate and send it before 15 days after the date of purchase.

Manufacturer reserves the right to modify or improve the manual or equipment.

**ATTENTION!!  
IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY  
WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

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