**AUXILAB S.L.**

DECLARATION OF CONFORMITY
Nahita centrifuge Digimax of Auxilab, S.L for the Directive
of Machines (89/392/CEE modified) and the regulations
adopted for their transposition

NAME OF THE MANUFACTURER / IMPORTER:
AUXILAB, S.L.

ADDRESS: **Polígono Morea Norte, 8**
31191 Beriáin (Navarra)

WE STATE THAT:
CENTRIFUGE DIGIMAX MODEL 2705/1
CODE 52705010

Are designed and manufactured according to:

- " ♦ Directive 89/392/CEE, including the modifications and national regulations that transpose them.
- " ♦ Directive 89/336/CEE modified over the electromagnetic compatibility.
- " ♦ Directive 73/23/CEE modified over the electric security.

And that the following harmonized rules have been applied (or a part of them):
 UNE 292-1/-2/-2/A1, UNE-EN 1050, UNE-EN 614-1, UNE-EN 1037, UNE-EN 1088, UNE-EN 547, UNE-EN 953, UNE-EN 294, UNE-EN 418, UNE-EN 894-1, UNE-EN 894-2, UNE-EN 954-1, UNE-EN 981, UNE-EN 60204-1, UNE 61010-1/A2, UNE-EN 61010-2-051.

BERIAIN, 15th March 2007

Signed by: **ALFONSO AINCIBURU SANZ**
DIRECTOR/MANAGER

Polígono Morea Norte, 8 31191 Beriain (Navarra) - Spain. Tel. 948 310 513 Fax 948 312 071
Internet: www.auxilab.es · Email: correo@auxilab.es

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This manual should be available for all users of these equipments. To get the best results and a higher duration of this equipment it is advisable to read carefully this manual and follow the processes of use.

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◆ Never force the stoppage of the centrifuge, it is a risk of accident and debilitation and it may produce as well a mixing of the separated parts.

◆ Change of fuse: the fuse compartment is located at the base of the socket.

◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.

◆ The centrifuge is supplied with a Schuko standard wire. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.

◆ Never try to repair the centrifuge by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the centrifuge (burns, hurts...) or damages in nearby equipments.

◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.

Cleaning

◆ When working at high speed there is a risk of rupture of the tubes. If this happens, the stainless steel rotor, the adapters and the tube holders must be cleaned and disinfected with a disinfectant agent and the broken parts must be replaced; otherwise different components of the centrifuge will suffer important damages and due to the lack of cleaning mistakes in subsequent analysis will take place.

◆ Pay special attention on aluminium alloy components; periodic treatment with antioxidant agents will prolong its life-span.

◆ For the cleaning of the different parts of the equipment, we recommend the use of neutral products with a pH value not higher than 6.

◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the centrifuge and produce an early ageing of the equipment.

◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

INSTRUCTIONS ON ENVIRONMENT PROTECTION

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location

Note: To stop the centrifuge before the selected time has passed, press STOP button (2.6). Never try to open the centrifuge with the rotor in movement, the lid is blocked.

Turn the equipment off by pressing the general power switch (1.4) and disconnect it from the net.

Security

- ◆ The centrifuge must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ◆ Nahita centrifuges are not suitable for treatment of flammable liquids, explosive samples or samples capable of forming explosive mixtures.
- ◆ When working with toxic or pathologic substances close hermetically the tubes and strictly observe the safety regulations corresponding to such substances.
- ◆ When you are not using the centrifuge for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ◆ Never try to repair the centrifuge by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the centrifuge (see Maintenance).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

6. MAINTENANCE AND CLEANING

To get the best results and a higher duration of this equipment it is essential to follow the processes of use.

Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ Use the appropriate tubes resistant to the R.C.F. used.
- ◆ Balance the centrifuge by using tubes with the same weight, shape and size and put them in geometrically symmetrical positions. If it is necessary, use tubes filled with water. In a centrifuge it is very important the concept of weight so it is advisable to have a balance to compensate the weight of the different tubes. This is absolutely necessary because an upset of the rotor may produce vibrations that debilitate the centrifuge and headstocks increasing the rate of broken tubes and producing a worse sedimentation.

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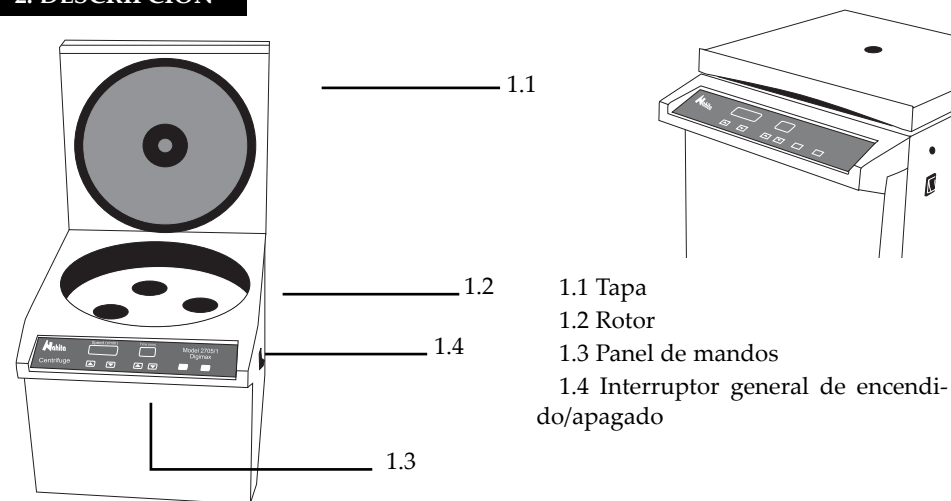
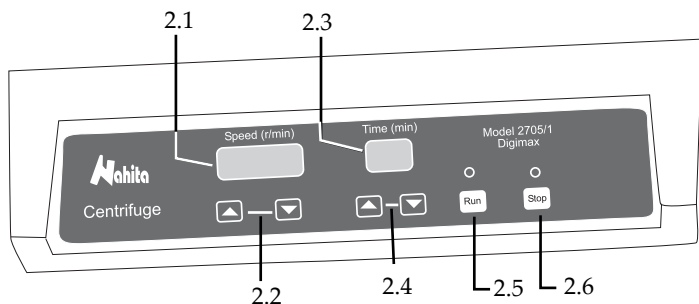
Note: When the equipment is disconnected from the net, the lid is automatically blocked until connecting the centrifuge again.

To open the lid manually without connecting the centrifuge, proceed as follows:

Remove the rubber caps located at both sides of the centrifuge (1.7) (SEE DRAWING). Insert into the holes a thin wire and press the electromagnetic device at the same time. The lid will open.

- ◆ Put the adapters to be used and make sure they are well placed. Check that the adapters swing.
- ◆ Put the tubes into the tube holders and make sure they are well placed.

NOTE: Balance the centrifuge by using tubes with the same weight, shape and size and put them in geometrically symmetrical positions. If it is necessary, use tubes filled with water.

- ◆ In a centrifuge it is very important the concept of weight so it is advisable to have a balance to compensate the weight of the different tubes. This is absolutely necessary because an upset of the rotor may produce vibrations that debilitate the centrifuge and headstocks increasing the rate of broken tubes and producing a worse sedimentation.

- ◆ Turn the centrifuge on by pressing the general power switch (1.4).

- ◆ Speed adjustment

Nahita centrifuge Digimax 2705 presents a speed range from 100 to 5000 r.p.m. with steps of 100 r.p.m. To select the desired working speed press ? ? buttons (2.2) and speed will increase or decrease on display until reaching the desired value.

- ◆ Time adjustment

Nahita centrifuge Digimax 2705 presents a time range from 1 to 99 minutes in steps of 1 minute. To select the desired time press ? ? buttons (2.4) and time will increase or decrease on display until reaching the desired value.

- ◆ Close de lid.

NOTE: Never turn the centrifuge on if the lid is open.

- ◆ Press RUN (2.5), the centrifuge begins to operate.
- ◆ On SPEED display (2.1) the speed value will increase until reaching the adjusted value.
- ◆ When the rotor reaches the selected speed, the timer starts the count down; the remaining operation time will be displayed on TIME display (2.3)

Note: When the centrifuge starts operating, the lid is blocked by an electromagnetic device. Never force the lid to open if the rotor is still in movement.

- ◆ Stop: when the selected time has passed the rotor speed begins to decrease.
- ◆ Do not take the samples out until the rotor is completely stopped.

- 36 Plastic tubes of 10 mL
- Net wire
- Instructions manual
- Warranty certificate

We will only accept any equipment return within 15 days after delivery and provided it comes in its original wrapping.

Installation

- ◆ Before using this instrument, it is convenient for you to familiarize with its components and basic essentials, as well as with its control functions.
- ◆ Please put the centrifuge on top of a horizontal, plane and stable table making a free space at least at 30 cm per side. During operation, people and dangerous material such as flammable or infectious liquids must be out of this area. Do not put the equipment near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc.
- ◆ The centrifuge is supplied with a Schuko standard wire 220-240 V, 50-60 Hz.
- ◆ Please insert the wire that feeds the AC electric current in the base of current provided with earth wire and to the other end to the centrifuge connector (1.5).

Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to persons suffered by the improper use of the electric connection. The tension should be 220 V, 50 Hz \pm 10%.

- ◆ If you are not using the centrifuge for a long period of time please make sure it is disconnected from the net and protected from dust (this way you will avoid accidents and will extend its working-life).

PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT WITH THE AIM OF ACHIEVING THE HIGHEST FEATURES AND THE MAXIMUM DURABILITY OF THE INSTRUMENT.

Placing the rotor

- ◆ The rotor is already placed; check that it is well fixed, for this, the grooves of the rotor must be well fitted in the axis of the rotor.
- ◆ Put the washer and the nut, and fix the rotor with a spanner.

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4. THEORETICAL FOUNDATIONS

Centrifugation is a separation technique based on the movement of the particles that are displaced to the distal end of the rotation axis depending on their shape and mass.

The Relative Centrifugal Force (R.C.F.) is the required force to separate particles. It is given as some value times that of gravity (g).

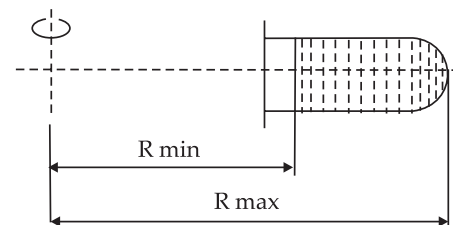
The R.C:F. is calculated with the following formula:

$$F.C.R. \ 1.118 \cdot 10^{-5} \cdot r \cdot n^2$$

1,118·10⁻⁵ is a constant.

r= radius, the horizontal distance in centimetres from the rotation axis to the bottom of the sample tube.

n= rotation speed expressed in revolutions per minute (r.p.m.)


5. INSTALLATION / SETTING UP

Preliminary inspection

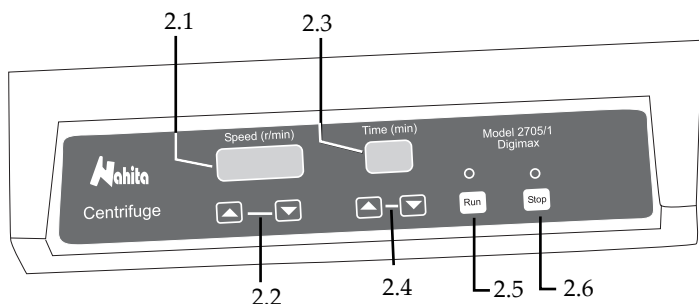- ◆ Unwrap the centrifuge; remove the upper cover by unscrewing the screws that hold it and then remove the four lateral covers that form the box by unscrewing the lateral screws.
- ◆ Take off the involving plastic and take off the polyspan protection in which it comes fitted.
- ◆ Check manually that the rotor spins freely without any obstacle and make sure the equipment does not present any damage because of the shipment. In case it presents any damage, tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.
- ◆ Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.
- ◆ Please check that all the accessories are enclosed with the equipment:
 - 4 Adapters for tubes of 250 mL
 - 4 Adapters for tubes of 50 mL
 - 4 Adapters for tubes of 10 mL
 - 4 Plastic tubes of 250 mL (approx.)
 - 8 Plastic tubes of 50 mL

- 2.1 SPEED display
 2.2 Speed increase/decrease button ▲ ▼
 2.3 TIME display
 2.4 Time increase/decrease button ▲ ▼
 2.5 RUN
 2.6 STOP

3. TECHNICAL SPECIFICATIONS

| | |
|------------------|--|
| Capacity | 4 adapters x 1 tube of 250 mL (67x105 mm) 4 adapters x 2 tubes of 50 mL (28x100 mm) 4 adapters x 9 tubes of 10 mL (16x80 mm) |
| Speed | 100-5000 r.p.m. (adjustable) |
| Max. R.C.F. | Tubes of 250 mL: 4420 g Tubes of 50 mL: 4700 g Tubes of 10 mL: 5000 g |
| Timer | 1-99.99 min |
| Rotor | Swing out |
| Speed adjustment | Digital |
| Time adjustment | Digital |
| Motor | Induction |
| Weight | 54 kg |
| Dimensions | 44x54x45 cm |
| Fuse | 8A |
| Power supply | 220-240 V ± 50-60 Hz |

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! ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

INDICE DE IDIOMAS

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INDEX OF CONTENTS

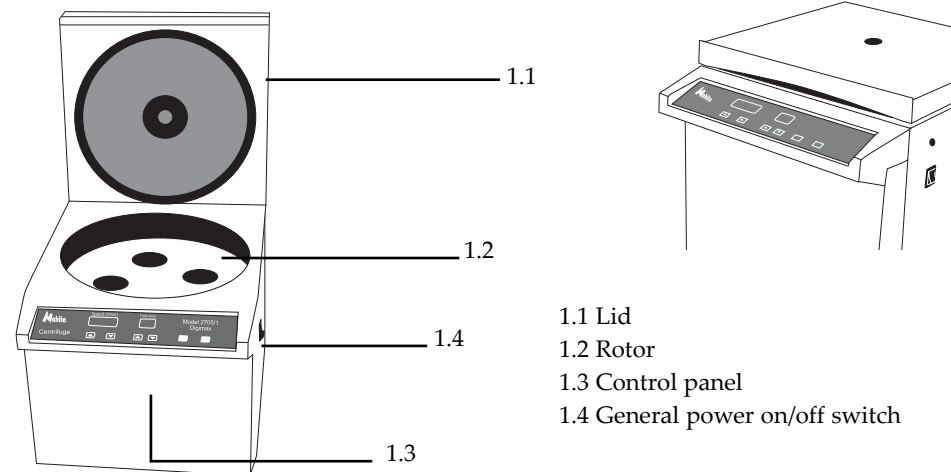
1. USES OF THE INSTRUMENT 13
 2. DESCRIPTION 13
 3. TECHNICAL SPECIFICATIONS 14
 4. THEORETICAL FOUNDATIONS 15
 5. INSTALLATION / SETTING UP 15
 6. MAINTENANCE AND CLEANING 18
 7. ANNEX I: CE CERTIFICATE 20

1. USES OF THE INSTRUMENT

Nahita centrifuge Digimax 2705 is a bench-top, silent and sturdy model without vibrations. It is equipped with three swing out adapters with capacity for 4 tubes of 250 mL, 8 tubes of 50 mL and 36 tubes of 10 mL and can reach a maximum speed of 5000 r.p.m.

Due to its simple conception and the quality of the materials, this centrifuge is specially indicated for any kind of analysis in Biology, Chemistry, Biochemistry, Medicine, Genetics, etc., since it can remain operating for long periods of time or keep a high daily rate of cycles without any difficulty

2. DESCRIPTION

- 1.1 Lid
- 1.2 Rotor
- 1.3 Control panel
- 1.4 General power on/off switch

Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita centrifuge Digimax 2705. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

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Please bear in mind the following:

- ◆ This manual is inseparable from the Nahita centrifuge Digimax 2705, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the centrifuge avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the centrifuge to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the equipment.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to asistencia@auxilab.es or by telephone: +34 807 117 040 (0.30 Euros/min).
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the equipment warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the centrifuge warranty.
- ◆ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Please do not forget filling the warranty certificate and send it before 15 days after the date of purchase.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

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