DECLARACIÓN DE CONFORMIDAD CE  
CE DECLARATION OF CONFORMITY  
DÉCLARATION DE CONFORMITÉ CE

El fabricante | The manufacturer | Le fournisseur:

AUXILAB, S.L.

Declara que el equipo | Declare that the equipment | Declare que l'appareil:

CENTRÍFUGA MINIVITE | CENTRIFUGE MINIVITE | CENTRIFUGEUSES MINIVITE

Código | Code | Code: 52716000

Modelo | Model | Modèle: 2716

Cumple las siguientes directivas | Meet the following directives | Accomplit les directives suivantes:

73/23/CE | Directiva de seguridad eléctrica  
Directive for electrical safety  
Directive de sécurité électrique

89/336/CE | Directiva de Compatibilidad electromagnética (CEM)  
Directive for electromagnetic compatibility (EMC)  
Directive de compatibilité électromagnétique (CEM)

Cumple las siguientes Normas: | Meet the following Standards | Accomplit les normes suivantes:

EN 61326 | Material eléctrico para medida control y uso en laboratorio. Requisitos de compatibilidad electromagnética (CEM).  
Electrical equipment for measurement, control and laboratory use. EMC requirements.  
Matériel électriques de mesure, de commande et de laboratoire. Prescriptions relatives à la CEM.

EN 61010-1 | Requisitos de seguridad de equipos eléctricos de medida, control y uso en laboratorio. Parte 1: Requisitos generales.  
Safety requirements for electrical equipments for measurement, control and laboratory use. Part 1: General requirements.  
Règles de sécurité pour appareils électriques de mesure, de régulation et de laboratoire. Partie 1: Prescriptions générales.

EN 61010-2-020 | Requisitos particulares para centrifugadoras de laboratorio.  
Particular requirements for laboratory centrifuges.  
Prescriptions particulières pour centrifugeuses de laboratoire.

Fdo: Alfonso Ainciburu Sanz  
DIRECTOR | GERENTE

BERIAIN a 28 de OCTUBRE de 2009

CENTRÍFUGAS MINIVITE  
MINIVITE CENTRIFUGES  
CENTRIFUGEUR MINIVITE

Nahita

Referencia / Code / Modèle 52716000

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For the cleaning of the different parts of the equipment, we recommend the use of neutral products with a pH value not higher than 6. These parts are also autoclavable.

- ◆ Pay special attention on aluminium alloy components; periodic treatment with antioxidant agents will prolong its life-span.
- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the Centrifuge and produce an early ageing of the equipment.
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

**ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

#### INSTRUCTIONS ON ENVIRONMENT PROTECTION

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

Thank you for choosing this equipment. We sincerely wish that you enjoy your centrifuge. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

**TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.**

Please bear in mind the following:

- ◆ This manual is inseparable from the Nahita Minivite Centrifuge, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the Centrifuge avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the Centrifuge to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the Centrifuge.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to [asistencia@auxilab.es](mailto:asistencia@auxilab.es).
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the Centrifuge warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the Centrifuge warranty.
- ◆ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

**ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

## INDEX OF LANGUAGES

Spanish .....	3-10
English .....	11-18
French .....	19-26

## INDEX OF CONTENTS

1. USES OF THE INSTRUMENT .....	12
2. DESCRIPTION .....	13
3. TECHNICAL SPECIFICATIONS .....	13
4. THEORETICAL FOUNDATIONS.....	14
5. INSTALLATION / SETTING UP .....	14
6. SECURITY .....	16
7. MAINTENANCE AND CLEANING .....	17
ANEXE I: CE CERTIFICATE .....	28

## 1. USES OF THE INSTRUMENT

The Nahita Minivite 2716 centrifuge is a bench-top, silent and sturdy model without vibrations. It is equipped with two angle headstocks with a capacity of 12 eppendorf-like tubes of 1.5 mL and 10 (53x13 mm) tubes of 5 mL, and presents an adjustable speed from 1000 to 16000 rpm.

Because of its simple conception and the quality of the materials this centrifuge is specially indicated for any kind of analysis in Biology, Chemistry, Biochemistry, Medicine, Genetics... It can remain operating for long periods of time or keep a high daily rate of cycles without any difficulty.

Examples of different activities in which the centrifuge acts as a separation technique:

- Separation of blood serum or plasma.
- Concentration of cells.
- Precipitation of nucleic acids.

## 7. MAINTENANCE AND CLEANING

To get the best results and a higher duration of this equipment it is essential to follow the processes of use.

*Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.*

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ Use the appropriate tubes resistant to the R.C.F. used.
- ◆ Balance the centrifuge by using tubes with the same weight, shape and size and put them in geometrically symmetrical positions. If it is necessary, use tubes filled with water. In a centrifuge it is very important the concept of weight so it is advisable to have a balance to compensate the weight of the different tubes. This is absolutely necessary because an upset of the rotor may produce vibrations that debilitate the centrifuge and rotors increasing the rate of broken tubes and producing a worse sedimentation.
- ◆ Never force the stoppage of the centrifuge, it is a risk of accident and debilitation and it may produce as well a mixing of the separated parts.
- ◆ Change of fuse: the fuse is situated near the socket.
- ◆ The change of brushes must be done by qualified staff when their length is less than 6 mm (after 500 h of operation).
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- ◆ The Centrifuge is supplied with a Schuko standard wire. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.
- ◆ Never try to repair the Centrifuge by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the Centrifuge (burns, hurts...) or damages in nearby equipments.
- ◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.

### Cleaning

When working at high speed there is a risk of rupture of the tubes. If this happens, the anodized aluminium rotors must be cleaned and disinfected with a disinfectant agent and the broken parts must be replaced; otherwise different components of the centrifuge will suffer important damages and due to the lack of cleaning mistakes in subsequent analysis will take place.

On speed display (2.1), the value of speed will increase until reaching the selected value.  
On time display (2.3), once the rotor spins at the selected r.p.m., time will decrease (the digit blinks) until reaching 00.

If you have selected the continuous mode, time display will show 00 (the digit blinks).

**Note:** When the centrifuge starts working, the lid is blocked by an electromagnetic lock(1.6). Never force its opening with the rotor running.

◆ Stop: when the programmed time finishes the rotor will decrease its speed until it stops.

**Note:** When working on continuous mode or if you need to stop the centrifuge before the programmed time, press Stop button (2.6). Never try to open the centrifuge before the rotor stops, the lid is blocked.

◆ Do not take the tubes until the rotor totally stops.

◆ Switch the equipment off by pressing the power on/off (1.4) switch and disconnect it from the main supply.

## 6. SECURITY

◆ The Centrifuge must be used by previously qualified staff that knows how the equipment works thanks to the user manual.

◆ Nahita centrifuges are not suitable for treatment of flammable liquids, explosive samples or samples capable of forming explosive mixtures.

◆ When working with toxic or pathologic substances, close hermetically the tubes and strictly observe the safety regulations corresponding to such substances.

◆ When you are not using the Centrifuge for a long period of time please make sure it is unplugged in order to avoid possible accidents.


◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).

◆ Never try to repair the Centrifuge by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).

◆ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the Centrifuge (see Maintenance).

◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

## 2. DESCRIPCIÓN

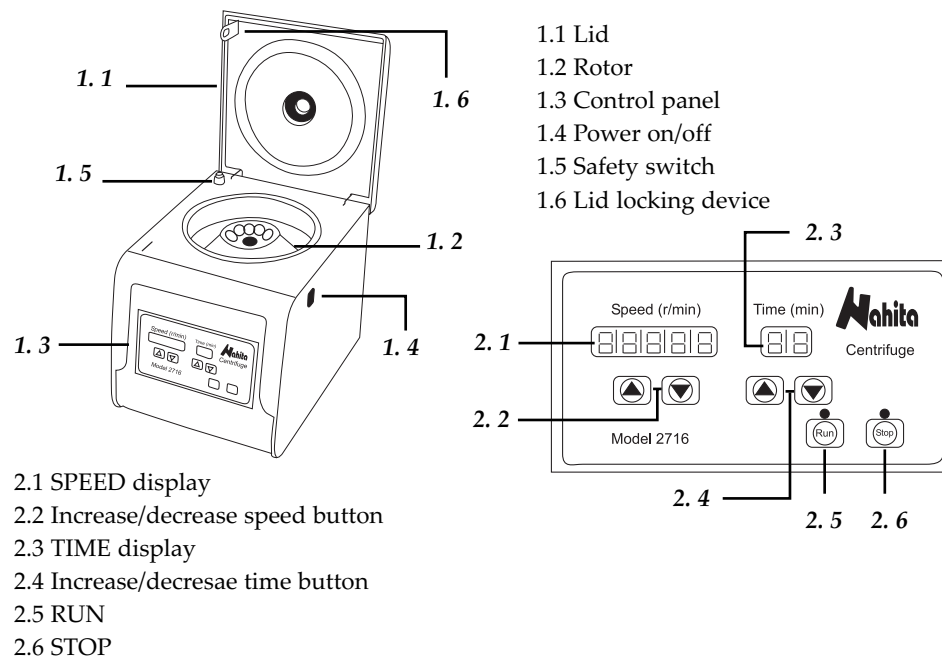## 3. TECHNICAL SPECIFICATIONS

Capacity	12 tubes of 1.5 mL. Eppendorf-like 10 tubes of 5mL.
Speed	2000-16000 r.p.m. (adjustable)
Max. F.C.R	17000 g
Timer	Continuous mode 1-99 min.
Rotors	Fixed angle: 12 tubes of 1.5 mL 10 tubes of 5 mL
Speed adjustment	Digital
Time adjustment	Digital
Motor	Elastic suspension with brushes
Weight	10.70 kg
Dimensions	32 X 27 X 24 cm
Fuse	5A
Power	220-240V ±50-60 HZ
Tubes size (length x diameter)	
1.5 mL tubes	41 x 10 mm
5 mL tubes	53 x 13 mm

#### 4. THEORETICAL FOUNDATIONS

Centrifugation is a separation technique based on the movement of the particles that are displaced to the distal end of the rotation axis depending on their shape and mass.

The Relative Centrifugal Force (R.C.F.) is the required force to separate particles. It is given as some value times that of gravity (g).


The R.C.F. is calculated with the following formula:

$$R.C.F.= 1.118 \cdot 10^{-5} \cdot r \cdot n^2$$

$1,118 * 10^{-5}$  is a constant.

$r$ = radius, the horizontal distance in centimetres from the rotation axis to the bottom of the sample tube.

$n$ = rotation speed expressed in revolutions per minute (r.p.m.).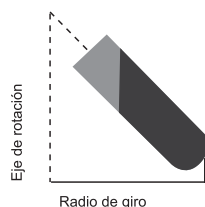

#### 5. INSTALLATION/SETTING UP

##### Preliminary inspection

Unwrap the centrifuge, take off the involving plastic and the take off the polispan protection in which it comes fitted.

Open the lid and check manually that the rotor spins freely, without any obstacle. Make sure that the equipment does not present any damage because of the shipment. In case it present any damage, tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.

Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.

Please check that all the accessories are enclosed with the equipment:

- Rotor for 12 eppendorf-like tubes of 1.5 mL (it comes placed inside the centrifuge).
- Rotor for 10 tubes of 5 mL (aprox.).
- Spanner.
- Schuko wire.
- Instructions manual.
- Warranty certificate.

**We will only accept any equipment return within 15 days after delivery and provided it comes in its original wrapping.**

##### Installation

Before using this instrument, it is convenient for you to familiarize with its components and basic essentials, as well as with its control functions. Please put the Centrifuge on top of a horizontal, plane and stable table making a free space at least at 30 cm per side. Do not put the Centrifuge near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc. During operation, people and dangerous material such as flammable or infectious liquids must be out of the working area.

The centrifuge is supplied with a standar Shuko wire.

Please insert the wire that feeds the AC electric current in the base of current 220V 50Hz  $\pm 10\%$  provided with earth wire and to the other end to the Centrifuge connector.

**Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to persons suffered by the improper use of the electric connection. The tension should be 220V 50Hz  $\pm 10\%$ .**

**PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT WITH THE AIM OF ACHIEVING THE HIGHEST FEATURES AND THE MAXIMUM DURABILITY OF THE INSTRUMENT.**

If you are not using the centrifuge for a long period of time please make sure it is disconnected from the net and protected from dust (this way you will avoid accidents and will extend its working-life).

##### Setting up

- ◆ Open the lid (1.1).
- ◆ Change the rotor if it is necessary (1.2).

**NOTE: The Nahita Minivite 2716 centrifuge is supplied with the headstock of 12 eppendorf-like tubes of 1.5 ml inside and ready to operate.**

If you wish to work with the 5 mL tubes rotor proceed as the following. With the spanner loosen the screw that holds the rotor to the rotor arm. Place the 10 tubes rotor by fitting the rotor grooves with the bolt of the rotor arm. Place the washer and tighten the screw.

- ◆ Put the tubes into their compartments. Make sure the tubes are well balanced and placed (See Maintenance and Cleaning).
- ◆ Switch the centrifuge on by pressing the power on/off (1.4).

##### Speed adjustment

The Nahita Minivite 2716 centrifuge presents an adjustable speed range from 2000-16000 r.p.m, in steps of 1000 r.p.m. To select the desired speed proceed as the following:

- ◆ Press ▲ ▼ (2.2) buttons, on display the speed value will increase or decrease in steps of 1000 r.p.m.

##### Time adjustment

The Nahita Minivite 2716 centrifuge presents an adjustable time range from 1-99 minutes, in steps of 1 minute. To select the desired time proceed as the following:

- ◆ Press ▲ ▼ (2.4) buttons, on display the time value will increase or decrease in steps of 1 minute.

**NOTE: To work on continuous mode select a 00 time value.**

- ◆ Close the lid (1.1).
- NOTE: Do not start the centrifuge without closing the lid.**
- ◆ Press RUN (2.5) , the centrifuge starts working.