AUXILAB S.L.

Material de laboratorio
Laboratory supplies

DECLARACIÓN DE CONFORMIDAD CE
CE DECLARATION OF CONFORMITY
DÉCLARATION DE CONFORMITÉ CE

El fabricante | The manufacturer | Le fournisseur:

AUXILAB, S.L.

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PROBLEMA	CAUSA	SOLUCIÓN
La centrifuga hace demasiado ruido	The surface on which the equipment is placed is unstable	Put the equipment on a plane and totally stable surface
	Mechanical components loose	Contact the Nahita Technical Assistance Department
	Motor is damaged	Contact the Nahita Technical Assistance Department
Unstable spinning speed	Failure of the control circuit	Contact the Nahita Technical Assistance Department
Unbalance protection	Samples are not well balanced	Put again the tubes into the rotor so as they are perfectly balanced
	Dirtiness on the rotor surface or into the holes for the tubes	Clean the rotor properly
	Failure of the motor	Contact the Nahita Technical Assistance Department

Instructions on environment protection

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location

Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita high speed centrifuge 2718. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

Please bear in mind the following:

- ◆ This manual is inseparable from the Nahita high speed centrifuge 2718, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the centrifuge avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the centrifuge to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the equipment.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to asistencia@auxilab.es or by telephone: +34 807 117 040 (0.30 Euros/min).
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the equipment warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the centrifuge warranty.
- ◆ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Please do not forget filling the warranty certificate and send it before 15 days after the date of purchase.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

INDEX OF LANGUAGES

Inglés	15-26
Spanish	4-14
French	27-38

INDEX OF CONTENTS

1. USES OF THE INSTRUMENT	16
2. DESCRIPTION	17
3. TECHNICAL SPECIFICATIONS	18
4. THEORETICAL FOUNDATIONS	18
5. INSTALLATION / SETTING UP	19
6. MAINTENANCE AND CLEANING	24
7. TROUBLESHOOTING	25
8. ANNEX I: CE CERTIFICATE	40

1. USES OF THE INSTRUMENT

Nahita high speed centrifuge model 2718 has been especially designed to be used in the field of Biochemistry, Molecular and Cellular Biology, Medicine and Research to carry out several routine techniques. With a solid design, quiet and easy-to-use, it presents a wide versatility of use thanks to the different rotors with which it can be used.

Among the different techniques that can be performed with this centrifuge it is worth mentioning the following:

- Separation of serum or plasma.
- Concentration of cells, bacteria or cellular organelles.
- Purification of nucleic acids, proteins, lipids.
- Protocols of Molecular biology, Genetics engineering, etc.

- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.
- ◆ The Nahita centrifuge 2718 can reach 18000 rpm. If any tube breaks, the interior of the centrifuge must be cleaned and disinfected with a disinfectant agent and the broken parts must be removed; otherwise different components of the centrifuge will suffer important damages and due to the lack of cleaning mistakes in subsequent analysis will take place.

ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

7. TROUBLESHOOTING

PROBLEM	CAUSE	SOLUTION
No display or disorderly display	Fuse blown	Change the fuse
	Problems with the cable that connects to display board	Contact the Nahita Technical Assistance Department
	Problems with the chip of the equipment	Contact the Nahita Technical Assistance Department
Display present but centrifugation can not be carried out	Failed to press Enter key for confirmation after setting parameters	Press Enter and try again
	Centrifuge lid is not well closed	Close the lid properly
	Speed fails to rise due to insufficient power supply	Change the power supply
	Failed to open the lid after last centrifugation finishes	Open the lid, close it properly and try again
	When pressing Start, the button makes no well contact	Contact the Nahita Technical Assistance Department
High noise	The equipment is in a tilting state	Put the equipment completely horizontal and levelled

In case you have any doubts do please immediately unplug the centrifuge (see Maintenance).

◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

6. MAINTENANCE AND CLEANING

To get the best results and a higher duration of this equipment it is essential to follow the processes of use.

NOTE: *All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.*


- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- ◆ The centrifuge is supplied with a Schuko standard wire. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.
- ◆ Never try to repair the centrifuge by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the centrifuge (burns, hurts...) or damages in nearby equipments.
- ◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.
- ◆ Use the appropriate tubes resistant to the R.C.F. used.
- ◆ Balance the centrifuge by using tubes with the same weight, shape and size and put them in geometrically symmetrical positions. If it is necessary, use tubes filled with water. In a centrifuge it is very important the concept of weight so it is advisable to have a balance to compensate the weight of the different tubes.
- ◆ This is absolutely necessary because an upset of the rotor may produce vibrations that debilitate the centrifuge and rotor increasing the rate of broken tubes and producing a worse sedimentation.
- ◆ Never force the stoppage of the centrifuge, it is a risk of accident and debilitation and it may produce as well a mixing of the separated parts.
- ◆ Change of fuse: unplugged the centrifuge. The fuse compartment is located at the base of the socket. With the help of a sharp object remove the fuse compartment and change the fuse.

Cleaning


- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the centrifuge and produce an early ageing of the equipment.

2. DESCRIPTION

- 1.1 Lid.
- 1.2 Rotor.
- 1.3 Control panel.
- 1.4 Lid opening button.
- 1.5 General switch.
- 1.6 Socket.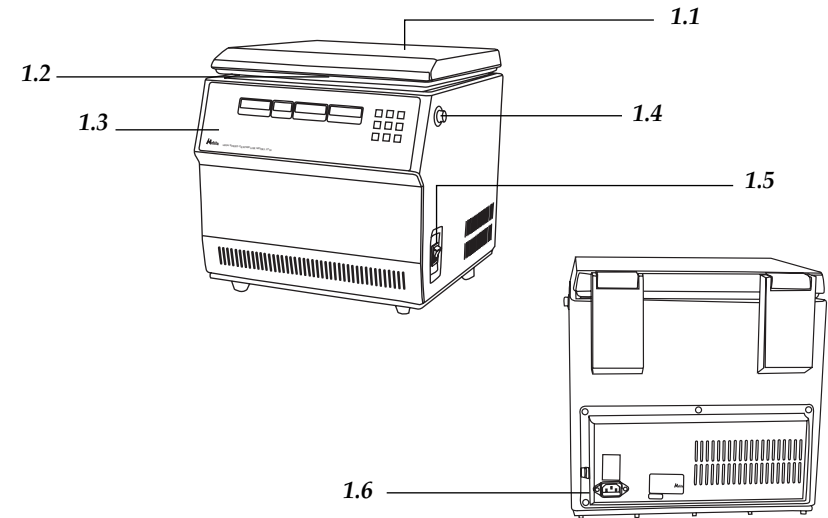

- 2.1 Rotor / program display.
- 2.2 Time display.
- 2.3 Speed / acceleration/braking rates display.
- 2.4 RCF display.
- 2.5 Time set.
- 2.6 Speed set.
- 2.7 RCF set.
- 2.8 Program set.
- 2.9 Parameter increase (▲) / decrease (▼) button.
- 2.10 Enter.
- 2.11 Start.
- 2.12 Stop.

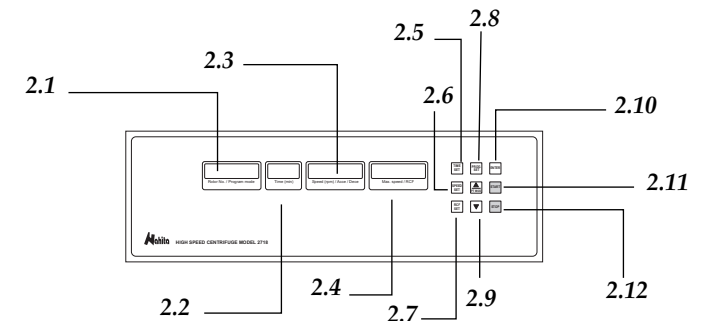3. TECHNICAL SPECIFICATIONS

<i>Model</i>	2718
<i>Rotor</i>	Not included
<i>Automatic detection of rotor</i>	Yes
<i>Maximum speed</i>	18.000 rpm (see rotor chart)
<i>Maximum RCF</i>	23545 g (see rotor chart)
<i>Timer</i>	1-999 min; continuous mode
<i>Memory</i>	12 programs
<i>Acceleration rate</i>	10 levels
<i>Braking rate</i>	10 levels
<i>Motor</i>	Induction, maintenance-free
<i>Noise</i>	< 68 dB
<i>Power</i>	220-230 V, 50 Hz
<i>Dimensions</i>	38x45x39 cm
<i>Weight</i>	42 Kg

ROTOR CHART

<i>Code</i>	<i>Name</i>	<i>Capacity</i>	<i>Max. speed.</i> <i>Max RCF</i>
92718001	H 240.2	24 x 0.2 mL, Eppendorf-like	18000 rpm / 18474 g
92718002	H 240.5	24 x 0.5 mL, Eppendorf-like	18000 rpm / 23907 g
92718003	H 2402	24 x 1.5/2 mL, Eppendorf-like	14000 rpm / 17465 g
92718004	H 1204	12 x 5 mL (12x75 mm)	14000 rpm / 15339 g
92718005	H 1211	12 x 10 mL (16x100 mm)	12000 rpm / 14489 g
92718006	P 361.5	24 microhaematocrit capillaries	12000 rpm / 16421 g

4. THEORETICAL FOUNDATIONS

Centrifugation is a separation technique based on the movement of the particles that are displaced to the distal end of the rotation axis depending on their shape and mass.

The Relative Centrifugal Force (R.C.F.) is the required force to separate particles. It is given as some value times that of gravity (g).

- With ▲ ▼ (2.9) choose the desired number of program (ctr 01, 02, 03, etc.) on program display (2.1).

- Press Enter (2.10) to select it.

Start centrifugation

Once time and speed values or a certain program have been selected, proceed as follows:

- ◆ Put the tubes into the rotor so as they are perfectly balanced and put the lid of the rotor.
- ◆ Close the lid of the centrifuge.
- ◆ Press Start (2.11) to start centrifugation.
- ◆ The centrifuge will stop after the selected time has passed and an acoustic signal will sound. When working on continuous mode or to stop the centrifuge before the operation time has passed, press Stop (2.12).

NOTE: if the parameter increase button (▲) is pressed, the centrifuge will be locked. To unlock it turn the equipment off and then on again by pressing the general switch (1.5).

Security

- ◆ The centrifuge must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ◆ Put the centrifuge on top of a horizontal, plane and stable table making a free space at least at 30 cm per side.
- ◆ Do not put the centrifuge near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc. Avoid vibrations, dust and dry environments.
- ◆ During operation, people and dangerous material such as flammable liquids or pathological material should be out of the working area.
- ◆ Nahita centrifuges are not suitable for treatment of flammable liquids, explosive samples or samples capable of forming explosive mixtures.
- ◆ As a safety mechanism the centrifuge is provided with a double mechanical lock with double switch that avoids operation if the lid is open and that automatically stops centrifugation in case of opening the lid during operation. Never try to open the lid while centrifuge is operating.
- ◆ When you are not using the centrifuge for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ◆ Never try to repair the centrifuge by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Try not to spill any liquid on the control panel, though it is properly insulated.

- On speed display (2.3) the value of speed (rpm) will change at the same time as RCF is selected.

- Press Enter (2.10) to confirm the RCF selected.

Programs

The centrifuge 2718 presents a memory to save up to 12 different centrifugation programs. On each program time, speed or RCF, and acceleration and braking rates can be adjusted.

◆ To save or modify a program proceed as follows:

- Press Prog. Set (2.8).

- With ▲ ▼ (2.9) select the desired number of program (ctr 01, 02, 03, etc.) shown on program display (2.1).

- Press again Prog. Set (2.8).

- Time display (2.2) will start flashing. With ▲ ▼ (2.9) select the desired time and press Enter (2.10).

- Speed display (2.3) will start flashing. If you want to select speed, enter the desired value with ▲ ▼ (2.9) and press Enter (2.10).

- RCF display (2.4) will start flashing. If you want to select RCF, enter the desired value with ▲ ▼ (2.9) and press Enter (2.10).

- The level of the acceleration rate will be shown on speed display (2.3). This parameter is related with the time that the rotor needs to reach the pre-selected speed. The centrifuge 2718 is provided with 10 acceleration rate levels, being the level 1 the fastest and the level 10 the slowest. With ▲ ▼ (2.9) select the desired level and press Enter (2.10).

- The level of the braking rate will be shown on speed display (2.3). This parameter is related with the time that the rotor needs to stop when spinning at the selected speed. The centrifuge 2718 is provided with 10 braking rate levels, being the level 1 the fastest and the level 10 the slowest. With ▲ ▼ (2.9) select the desired level and press Enter (2.10).

NOTE: When using a rotor for the first time, the centrifuge will use as default the first program saved; if there are no programs saved then the acceleration and braking rates will be set at level 5 as default.

- All parameters have been selected and the program has been saved.

NOTE: If during the process of setting a program Enter (2.10) is not pressed for a long period of time, the centrifuge will automatically save the data of the program.

◆ To select a program previously saved on the memory of the centrifuge, proceed as follows:

- Press Prog Set (2.8).


The R.C:F. is calculated with the following formula:

$$F.C.R. = 1.118 \cdot 10^{-5} \cdot r \cdot n^2$$

- $1.118 \cdot 10^{-5}$ is a constant..

- r = radius, the horizontal distance in centimetres from the rotation axis to the bottom of the sample tube.

- n = rotation speed expressed in revolutions per minute (r.p.m.).

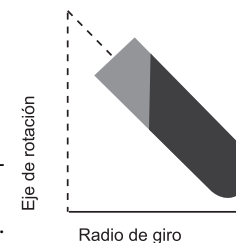5. INSTALLATION / SETTING UP

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Unwrap the centrifuge, take off the involving plastic and take off the polyspan protection in which it comes fitted. Take off all protections and, without connecting the centrifuge to the net, check manually that the rotor spins freely, without any obstacle. Make sure that the equipment does not present any damage because of the shipment. In case it presents any damage, tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.

Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.

Please check that all the accessories are enclosed with the equipment:

- ◆ Schuko wire.
- ◆ User's manual.
- ◆ Warranty.

We will only accept any equipment return within 15 days after delivery and provided it comes in its original wrapping.

Installation

Before using this instrument, it is convenient for you to familiarize with its components and basic essentials, as well as with its control functions.

PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT WITH THE AIM OF ACHIEVING THE HIGHEST FEATURES AND THE MAXIMUM DURABILITY OF THE INSTRUMENT.

◆ Put the centrifuge on top of a horizontal, plane and stable table making a free space at least at 30 cm per side.

◆ Do not put the equipment near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc. Avoid vibrations on the working area.

◆ During operation, people and dangerous material such as flammable or infectious liquids must be out of the working area.

◆ Please insert the Schuko wire that feeds the AC electric current in the base of current 220 V, 50 Hz \pm 10% provided with earth wire and to the other end to the centrifuge socket (1.6).

Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to persons suffered by the improper use of the electric connection. The tension should be 220 V, 50 Hz \pm 10%.

◆ **Placing the rotor (Fig. 3)**


Carefully insert the rotor in the rotor arm so as the two pins (3.1) located in the rotor arm fit with the two holes (3.2) of the base of the rotor.

◆ Put the supplied nut (3.3) and tighten it firmly with a spanner.

Finally, put the lid of the rotor (3.4).

NOTE: When closing the lid of the centrifuge and turning the equipment on, the rotor will be automatically detected and its name and maximum speed and RCF of use will be displayed.

If you are not using the centrifuge for a long period of time please make sure it is disconnected from the net and protected from dust (this way you will avoid accidents and will extend its working-life).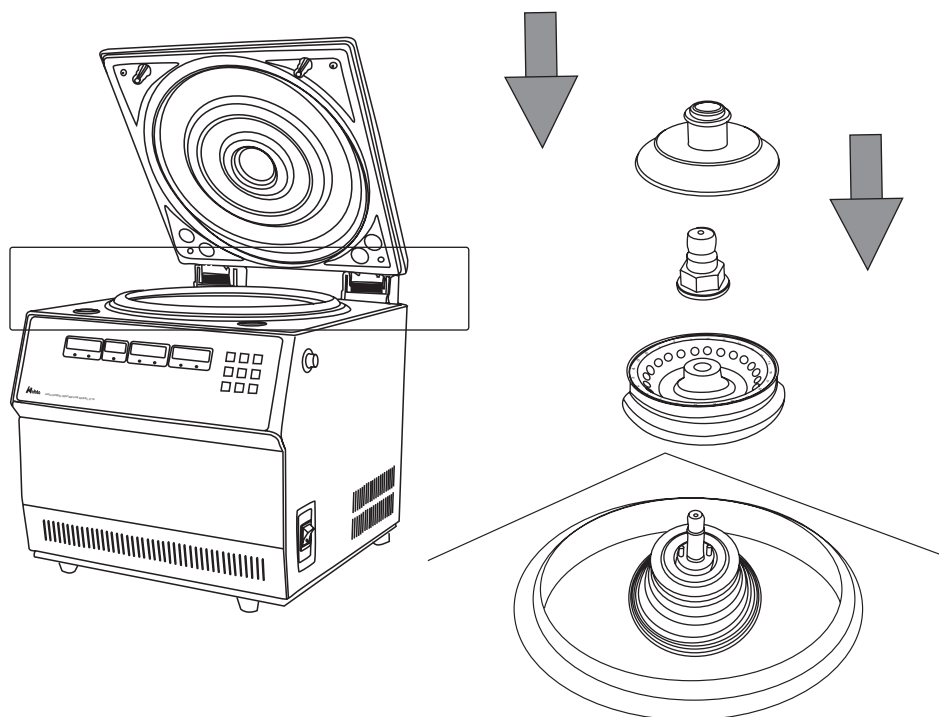

Operation mode

Press the general switch (1.5) to turn the equipment on. Then, proceed with the selection of the different parameters of the centrifugation protocol.

Time adjustment

The centrifuge 2718 can operate on continuous mode or during adjustable periods of time within a range of 1-999 min.

◆ To work on continuous mode, proceed as follows:

- If time display shows a number, press Time set (2.5) twice; time display (2.2) will show "-----" flashing.

- Press Enter (2.10); time display (2.2) will show "-----".

The centrifuge is ready to work on continuous mode; it will keep on working until pressing Stop (2.12).

◆ To work during concrete periods of time, proceed as follows:

- If time display shows a number, press Time set (2.5) once; time display (2.2) will show the previously selected time flashing.

NOTE: If time display (2.2) shows "-----" (continuous mode) press twice Time set (2.5) to adjust the working time to a certain period of time.

- Select the working time with the buttons \blacktriangle \blacktriangledown (2.9) within a range of 1-999 min in steps of 1 min.

- Press Enter (2.10) to confirm the selected time.

NOTE: when programming the centrifuge take into account that countdown begins when centrifugation starts not when selected speed is reached

Speed or RCF adjustment

The centrifuge 2718 allows adjustment of speed (rpm) or RCF (g), which depend on the type of rotor used (see Rotor chart). The centrifuge automatically detects the type of rotor and sets the upper limit of the speed and RCF of use.

◆ For speed adjustment proceed as follows:

- Press Speed set (2.6).

- With \blacktriangle \blacktriangledown (2.9) select the value of the centrifugation speed (in steps of 100 rpm) shown on speed display (2.3).

- On RCF display (2.4) the value of RCF will change at the same time as speed is selected.

- Press Enter (2.10) to confirm the selected speed.

◆ For RCF adjustment proceed as follows:

- Press FCR set (2.7).

- With \blacktriangle \blacktriangledown (2.9) select the value of the RCF (in steps of 1 g) shown on RCF display (2.4).