**MANTAS CALEFACTORAS  
HEATING MANTLES  
CHAUFFE BALLONS**

**Nahita**

**Serie / Series / Série 655**

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Thank you for choosing this equipment. We sincerely wish that you enjoy your heating mantles Nahita series 655. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user. The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of *Maintenance*, where we expose the security notes.

**TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.**

Please bear in mind the following:

- ◆ This manual is inseparable from the heating mantles, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the heating mantles avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the heating mantles to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the heating mantles .
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to [asistencia@auxilab.es](mailto:asistencia@auxilab.es) or by telephone: +34 807 117 040 (0,30 Euros/min).

◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).

◆ Overhaul is not covered by the heating mantles warranty.

Operations made by non-qualified staff will automatically produce a loss of the heating mantles warranty.

Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.

◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.

◆ Please do not forget filling the warranty certificate and send it before 15 days after the date of purchase.

Manufacturer reserves the right to modify or improve the manual or equipment.

**ATTENTION!**

**IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

**INDEX OF LANGUAGES**

Spanish .....	2-7
English .....	8-13
French .....	14-19

**ANNEX I: CE CERTIFICATE**

**AUXILAB S.L.**

**CE DECLARATION OF CONFORMITY  
HEATING MANTLES of AUXILAB, S.L. for the Directive of Machines (89/392/CEE modified) and the regulations adopted for their transposition**

NAME OF THE MANUFACTURER / IMPORTER:

**AUXILAB, S.L.**

ADDRESS:

**Polígono Morea Norte, 8  
31191 Beriáin (Navarra)**

WE STATE THAT:

**NAHITA HEATING MANTLES SERIES 655  
(MODELS 50655220, 50655250, 50655310, 50655320, 50655330 and 50655350)**

Are designed and manufactured according to:

- ◆ Directive 89/392/CEE, including the modifications and the national regulations that transpose them.
  - ◆ Directive 73/23/CEE modified over the electric security.
  - ◆ Directive 89/336/CEE modified over the electromagnetic compatibility.
- And that the following harmonized rules have been applied (or part of them):

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BERIAIN 21st June 2004

Signed by: ALFONSO AINCIBURU SANZ  
DIRECTOR/MANAGER

Polígono Morea Norte, 8 31191 Beriain (Navarra) - Spain. Tel. 948 310 513 Fax 948 312 071  
Internet: [www.auxilab.es](http://www.auxilab.es) · Email: [correo@auxilab.es](mailto:correo@auxilab.es)

**5. MAINTENANCE AND CLEANING**

To get the best results and a higher duration of this equipment it is essential to follow the processes of use.

*Note:* All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- ◆ The heating mantles is supplied with a Schuko standard wire and it should be connected to a current wire provided with an earth wire, it should be handy to be disconnected in case of emergency.
- ◆ Never try to repair the heating mantles by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the heating mantles (burns, hurts...) or damages in nearby equipments.
- ◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.

Cleaning

- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the heating mantles and produce an early ageing of the equipment.
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

 <p style="margin: 0;"><b>ATTENTION!</b> IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.</p>
---


**INDEX OF CONTENTS**

1. USES OF THE INSTRUMENT .....	9
2. DESCRIPTION .....	9
3. TECHNICAL SPECIFICATIONS .....	9
4. INSTALLATION / SETTING UP .....	10
5. MAINTENANCE AND CLEANING .....	12
ANNEX I: CE CERTIFICATE .....	13

**1. USES OF THE INSTRUMENT**

Nahita heating mantles with electronic temperature regulation are basic instruments in every laboratory to be used in distillation, digestion, evaporation, extraction and boiling processes.

**2. DESCRIPTION**

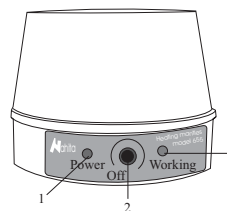1. *Power:* Current intake indicator, when we plug the instrument, the red pilot lightens.
2. *Off:* Temperature adjustment knob with on/ off functions.
3. *Working:* Green pilot which shows temperature's intensity (the more bright is the light, the more temperature reaches the sample).

**3. TECHNICAL SPECIFICATIONS**

Nahita heating mantles with electronic temperature adjustment are basic instruments in every laboratory. Chrome- nickel heater elements are homogeneously distributed around all the glass wool fabric to make a better and wider contact with the flask thanks to its semi-spherical shape. This way you can obtain a quicker and more homogeneous heating and lower energy consumption.

The external furniture is made of furnace-painted steel and the inner chamber is thermally sealed with mineral fiber wool. An electronic adjustment knob with switch enables the mantle's energy variation, therefore the temperature applied on the flask. The equipment is provided with connection pilot light, On pilot light and earth connection.

The protective fuse varies depending on the model supplied.

Resistance's temperature range goes from room temperature up to 360-380° C and the electric supply is of 220V, 50Hz ±10%.

The following models are available:

Code	Capacity (ml)	Consumption (w)
50655210	100	100
50655220	250	150
50655250	500	250
50655310	1000	350
50655320	2000	450
50655330	3000	600
50655350	5000	800

#### 4. INSTALLATION / SETTING UP

##### Preliminary inspection

- ◆ Unwrap the heating mantle, take off the involving plastic and take off the polispan protection in which it comes fitted. Take off all the protective items and, without connecting the heating mantle to the net, make sure that it does not present any damage because of the shipment. In case the equipment presents any damage tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.
- ◆ Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.
- ◆ Please check that all the accessories are enclosed with the equipment:
  - Rod
  - Instruction manual
  - Warranty certificate

*We will not accept any equipment in return period unless it comes in its original wrapping.*

##### Installation

- ◆ Before using this instrument, it is convenient for you to familiarize with its components and basic essentials.

**PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT.**

- ◆ Please put the heating mantle on top of a horizontal, plane and stable table making a free space at least at 30 cm per side. Do not put it near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc.
- ◆ Avoid inflammable or toxic substances in the working area.
- ◆ The heating mantle is supplied with a Schuko standard wire
- ◆ Please insert the wire that feeds the AC electric current in the base of current 220V 50Hz  $\pm 10\%$  provided with earth wire and to the other end to the heating mantle connector.

*Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to persons suffered by the improper use of the electric connection. The tension should be 220V 50Hz  $\pm 10\%$ .*

- ◆ If you are not using the heating mantle for a long period of time please make sure it is disconnected from the net and protected from dust (this way you will avoid accidents and will extend its working-life).

##### Setting up

- ◆ Before starting the mantle, make sure there is nothing inside that can melt.
- ◆ Plug the equipment to the net; red pilot light (1) will lighten showing energy reaches the equipment and will stay lighten as the cable is connected to the net.
- ◆ Move the temperature knob (2) to the right until you hear a "click"; this shows the heating mantle is switched on. To adjust the temperature, keep moving the knob to the right.
- ◆ Green pilot light (3) -Working- shows the temperature variation. Thus, the more intense is the light, the more temperature the sample reaches.
- ◆ Once you have finished working with the equipment he highly recommend you to unplug it, as its adapter electric components make a little energy consumption. This consumption can cause the early ageing of the heating mantle, as well as provoke accidents due to overheating or an unnecessary waste of energy.

##### Security

- ◆ The heating mantles must be used by previously qualified staff that knows how it works thanks to the user manual.
- ◆ You should put the heating mantles in a horizontal plane stable table, having a safety area of at least 30 cm per side.
- ◆ Do not place the heating mantles near any warm supply (burners, blowlamps, etc), nor expose it directly to the sun. Avoid vibrations, dust and dry environments.
- ◆ During its functioning dangerous materials such as flammable or pathological substances must be out of the safety area.
- ◆ When you are not using the heating mantles for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ◆ Never try to repair the heating mantles by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the heating mantles (see *Maintenance*).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.