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This manual should be available for all users of these equipments. To get the best results and a higher duration of this equipment it is advisable to read carefully this manual and follow the processes of use.

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AUXILAB S.L.

CE DECLARATION OF CONFORMITY

Multifunction electronic balances serie 5021 of Auxilab,S.L for the Directive of machines (89/392/CEE modified) and the regulations adopted for its transposition

NAME OF THE MANUFACTURER / IMPORTER:

AUXILAB, S.L.

ADDRESS:

**Polígono Morea Norte, 8
31191 Beriain (Navarra)**

WE STATE THAT:

**MULTIFUNCTION ELECTRONIC BALANCES SERIES 5021
Models 55021220, 55021230, 55021250, 55021260, 55021310**

Are designed and manufactured according to:

- ◆ Directive 89/392/CEE, including the modifications and the national regulations that transpose them.
- ◆ Directive 73/23/CEE, modified over the electric security.
- ◆ Directive 89/336/CEE, modified over the electromagnetic compatibility.
- ◆ And that the following harmonized rules have applied (or part of them):
UNE 292-1/-2/-2/A1, UNE-EN 55022, UNE-EN 55024, UNE-EN 61000, UNE-EN 61010, UNE-EN 61326.

BERIAIN 31th May 2004

Signed by: ALFONSO AINCIBURU SANZ
DIRECTOR/MANAGER

Polígono Morea Norte, 8 31191 Beriain (Navarra) - Spain. Tel. 948 310 513 Fax 948 312 071
Internet: www.auxilab.es · Email: correo@auxilab.es

Note: According to the in force legislation regarding "Non-automatic weighting instruments" in which balances are included, by means of writ dating from 22nd October, 1994 (BOE 1/3rd/95), these balances must not be used for:

- Comercial transactions
- Calculating taxes, tariffs, rates, indemnities and other similar canons
- Judicial surveys
- Pharmaceutical medicine preparations, as well as analysis made in medical or pharmaceutical laboratories
- Determining the price or total amount in sale price or in pre-packaged preparations

INSTRUCTIONS ON ENVIRONMENT PROTECTION

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

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**8. TROUBLESHOOTING**

TROUBLE	CAUSE	SOLUCIÓN
Digits are not displayed	The AC adapter is not connected The battery is run down	Plug the adapter in the electrical current Change the battery
The value on display is inconstantly shown	Low battery	Change the battery
The displayed value is wrong	The balance is not calibrated The balance has not been tared before weighing	Calibrate the balance Tare the balance before each weighing
The value on display does not stabilizes	The adapter is damaged	Calibrate the balance Tare the balance before each weighing
The balance does not calibrate	The weight used for calibration is not the correct one Zero is not adjusted	Use the correct weight See Tare function
Err1	Reading error	Turn the balance on again
Err2	The calibration weight is not the correct one	Use the correct weight
Err4	The load is too large to 125% the capacity	Decrease the load
Full	The load is too large to 125% the capacity Hardware error due to a load too high	Decrease the load Adjust the zero. If error remains, contact to Nahita Technical Assistance Department
FAIL	Hardware error due to a load too low	Adjust the zero. If error remains, contact to Nahita Technical Assistance department

7. MAINTENANCE AND CLEANING

To get the best results and a higher duration of the balance it is essential to follow the processes of use.

Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ Prevent the balance from sudden movements and knocks, as well as from direct sunlight or air flows. The balance is a precision instrument, you must handle it carefully.
- ◆ Place the balance in a location with adequate ventilation to prevent heat accumulation. Never use the balance in a wedged location as for example a shelf.
- ◆ The balance is supplied with an adapter. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.
- ◆ Never unplug the adapter by pulling the wire, do it from the base.
- ◆ If you use batteries, remove them when you are not using the balance for a long period of time.
- ◆ Never use sharp objects as pens, etc... to press the buttons of the control panel; only use your fingers.
- ◆ Never place on the pan an object heavier than the maximum capacity of the balance, the sensor could be damaged.
- ◆ Neither submerge the balance nor spill liquids on it.
- ◆ If any liquid comes into contact with the electric parts of the balance please do immediately disconnect it from the net and send it to the technical service, as soon as possible, for its checking and adjustment.
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.

Cleaning

- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the balance and produce an early ageing of the equipment.
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

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T (tare) key (1.4) again and 0.0 g (or the corresponding zero) will be displayed.

NOTE: The balance presents a tare range the same as its maximum capacity. This maximum must be taken into account since if the maximum capacity is tared nothing more could be weighted. (For example: if 450 g are tared in a balance with a maximum capacity of 500 g, only 50 g more will be weighted).

Piece counting-function

- ◆ Make sure that there is nothing on the pan and check that 0.0 g is displayed.
- ◆ Put 10 pieces on the pan since it is the minimum quantity needed to use this function.
- ◆ Press the N (counting) key (1.8), at first "COUNT ON" will be displayed and then "10 pcs". Remove the pieces from the pan and the balance is ready to count pieces.
- ◆ Put the pieces to be count on the pan and automatically the quantity will be displayed.
- ◆ Once the operation is finished, press the N (counting) key (1.8) to return to weighing mode; "COUNT OFF" will be displayed and the balance will be ready to weight.
- ◆ When you are not using the balance for a long period of time, make sure it is disconnected from the net or in case of using the batteries remove them from the corresponding compartment.

Security

- ◆ The balance must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ◆ Put the balance on top of a horizontal, plane and stable table making a free space of at least 30 cm per side.
- ◆ During operation, dangerous products such as flammable liquids or pathological material must be out of the working area.
- ◆ Do not put the balance near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc. Avoid vibrations, dust and dry environments.
- ◆ When you are not using the balance for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ◆ Never try to repair the balance by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the equipment (see Maintenance).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

Important!! Do not move the balance during calibration

- ◆ Make sure that there is nothing on the pan (1.3) and check that 0.0 g is displayed. Press C (calibration) key (1.7) and "C--- 0" will be displayed.
- ◆ Press the C (calibration) key again (1.7). "Hold" will be displayed and after a few seconds, the weight of the calibration weight needed to calibrate the balance will be shown on display.

TABLE 2: "Calibration weight"

Code	Calibration weight (NOT INCLUDED)
55021220	200 g
55021230	200 g
55021250	500 g
55021260	500 g
55021310	1000 g

- ◆ Put the corresponding calibration weight on the pan and press C key again; "Hold" will be displayed while the balance is calibrating.
- ◆ Wait until the weight of the calibration weight is displayed, then, the balance automatically enters in weighing mode. Remove the weight from the pan and 0.0 g will be displayed.
- ◆ After calibration the balance is ready to be used.

6. FUNCTIONS OF THE BALANCE**Tare function**

Tare function is used when a substance contained in a recipient is needed to be weight; for example liquids or solid reagents contained in a beaker to prepare a solution. For this function, proceed as follows:

- ◆ Make sure that there is nothing on the pan (1.3) and check that 0.0 g is displayed (when using a weighing unit different from grams, its corresponding zero will be displayed).
- ◆ Put the recipient on the pan (1.3) and press T (tare) key (1.4). The display will shown 0.0 g (when using a weighing unit different from grams, its corresponding zero will be displayed), wait until a circle is displayed on the lower left side and then, the balance is ready to weight.
- ◆ Once the weighing operation is finished, remove the recipient from the pan and its weight in negative will be displayed. To continue with the normal use of the balance, press

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tions or mistakes in weighing.

- ◆ Make sure that there is nothing on the pan and press the power switch (1.1) and then the ON/OFF key (1.5). The display will light up (1.9) and the balance will make an auto-test that finishes when 0.0 g is visualized on display (when using a weighing unit different from grams, its corresponding zero will be displayed).
- ◆ By pressing the unit configuration key (1.6), select the desired weighing unit: grams (g), pounds (lb), ounces (oz), troy ounces (ozt), grains (gn), carats (ct), pennyweights (dwt), tael (tl). The following chart shows the unit conversion.

TABLE 1: "Unit conversion"

Abbreviation	Name	Conversion
g	Gram	1 g
Kg	Kilogram	1000 g
oz	Ounce	28.3495231 g
Ozt	Troy ounce	31.10347668 g
dwt	Pennyweight	1.55517384 g
GN	Grain (UK)	0.06479789 g
lb	Pound	453.7205 g
ct	Carat	0.2 g
tl	Tael (HK)	37.7994 g

- ◆ The balance will be ready to be used when a black circle is displayed on the lower left side of the screen (it indicates that the balance is stabilized).

NOTE: To prevent possible oscillations in weighing, it is important to avoid vibrations on the table in which the balance is situated.

- ◆ Switch the balance off once operation is finished. Press the ON/OFF key (1.5) and then the power switch (1.1).

NOTE: The balance automatically switches off after five minutes of no use (approx.).

5. CALIBRATION

Before starting using the balance and whenever observing important variations in weighing or not correct weighings, balance should be calibrated. Some factors such as variation of gravity depending on latitude, change of working place as well as an inappropriate use of the equipment lead to the necessity of recalibrate the balance.

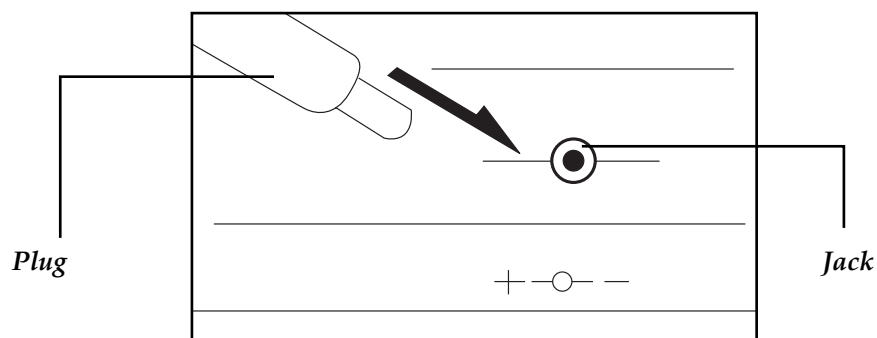
- ◆ Please put the balance on top of a horizontal, plane and stable table making a free space of at least 30 cm per side. Do not put the balance near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc.
- ◆ Put the stainless steel pan on the weighing sensor (Se puede hacer un dibujo explicando)
- ◆ Nahita electronic balances series 5021 can operate with batteries or by connecting them to the electric current by an AC adapter.

Battery installation

- ◆ Before installing the battery, remove the pan to prevent damage to the balance
- ◆ Put the balance upside down and remove the lid of the battery compartment by pressing up the tab.
- ◆ Place the 9 V battery inside the compartment bearing in mind to put it correctly according to its polarity.
- ◆ Replace the lid of the compartment

AC adapter installation

- ◆ Connect the adapter +5 V/+9 V into the jack on the back side of the balance. Then, plug the adapter into the electric current provided with earth.

- ◆ Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to people suffered by the improper use of the electric connection. The tension should be 220 V, 50 Hz \pm 10%.

Setting up

NOTE: Before using the balance, it must be calibrated according to "Calibration" point of this manual. Recalibration must be performed whenever observing important varia

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4. INSTALLATION / SETTING UP**Preliminary inspection**

Unwrap the balance, take off the involving plastic and take off the polispán protection in which it comes fitted. Take off all the protective items and, without connecting the balance to the net, make sure that it does not present any damage because of the shipment. In case the balance presents any damage tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.

Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.

Please check that all the accessories are enclosed with the equipment:

- Pan
- Adapter 220/240 V,50-60 HZ, 9 V AC
- Battery 9 V
- Instructions manual
- Warranty

We will only accept any equipment return within 15 days after delivery and provided it comes in its original wrapping.

Installation

◆ Before using the balance, it is convenient for you to familiarize with its components and basic essentials, as well as with its control functions.

PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT WITH THE AIM OF ACHIEVING THE HIGHEST FEATURES AND THE MAXIMUM DURABILITY OF THE INSTRUMENT.

ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

INDEX OF LANGUAGES

Spanish 2-12
 English..... 13-23

INDEX OF CONTENTS

1. USES OF THE INSTRUMENT 14
 2. DESCRIPTION 14
 3. TECHNICAL SPECIFICATIONS 15
 4. INSTALLATION / SETTING UP 15
 5. CALIBRATION 17
 6. FUNCTIONS OF THE BALANCE 18
 7. MAINTENANCE AND CLEANING 20
 8. TROUBLESHOOTING 21
 ANNEX I: CE CERTIFICATE 23

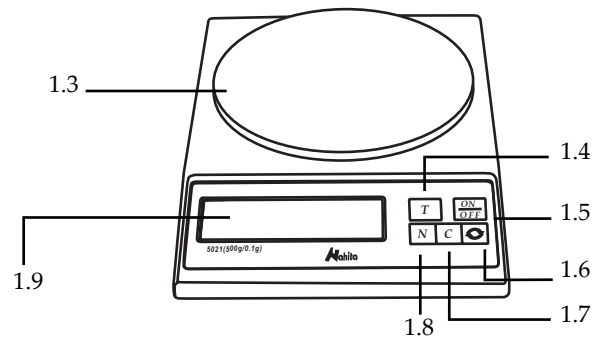
1. USES OF THE INSTRUMENT

Nahita multifunction electronic balances are essential instruments at any laboratory when weighing tasks are needed to be carried out. Developed with the most advanced technology they are reliable and durable instruments, with a solid, clean and ergonomic design suitable for industrial applications.

Nahita balances series 5021 are equipped with piece-counting function and weighing in grams, pounds and ounces (among other units) which together with an easy installation will allow to work in a wide range of fields (from educational and research fields to industry).

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Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita electronic balance, series 5021. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

Please bear in mind the following:

- ◆ This manual is inseparable from the Nahita electronic balance series 5021, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the balance avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the balance to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the balance.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to asistencia@auxilab.es or by telephone: +34 807 117 040 (0.30 Euros/min).
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the equipment warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the warranty.
- ◆ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Please do not forget filling the warranty certificate and send it before 15 days after the date of purchase.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.