DECLARACIÓN DE CONFORMIDAD CE
CE DECLARATION OF CONFORMITY
DÉCLARATION DE CONFORMITÉ CE

El fabricante | The manufacturer | Le fournisseur:

AUXILAB, S.L.

Declara que el equipo | Declare that the equipment | Declare que l'appareil:

BALANZA DIGITAL | DIGITAL BALANCES | BALACES DIGITAL

Código | Code | Code: 55041220, 55041250, 55041320, 55041350

Modelo | Model | Modèle: 5041

Cumple las siguientes directivas | Meet the following directives | Accomplit les directives suivantes:

73/23/CE Directiva de seguridad eléctrica	89/336/CE Directiva de Compatibilidad electromagnética (CEM)
Directive for electrical safety	Directive for electromagnetic compatibility y (EMC)
Directive de sécurité électrique	Directive de compatibilité électromagnétique (CEM)

Cumple las siguientes Normas: | Meet the following Standars | Accomplit les normes suivantes:

EN 61326 | Material eléctrico para medida control y uso en laboratorio
Requisitos de compatibilidad electromagnética (CEM.)
Electrical equipment for measurement, control and laboratory use
EMC requirements.
Matériel électriques de mesure, de commande et de laboratoire
Prescriptions relatives à la CEM.

EN 61010-1 | Requisitos de seguridad de equipos eléctricos de medida, control y uso en laboratorio
Parte 1: Requisitos generales
Safety requirements for electrical equipments for measurement, control and laboratory use
Part 1: General requierements.
Règles de sécurité pour appareils électriques de mesurage, de régulation et de laboratoire
Partie 1: Prescriptions générales.

Fdo: Alfonso Ainciburu Sanz
DIRECTOR | GERENTE

BERIAIN a 05 de MARZO de 2009

Nahita

Serie / Series / Série 5041

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7. TROUBLESHOOTING

TROUBLE	CAUSE	SOLUTION
Digits are not displayed	The AC adapter is not connected	Plug the adapter in the electrical current
	Batteries are run down	Change the batteries
The value on display is inconsistently shown	Low batteries	Change the batteries
The value on display does not stabilize	The adapter is damaged	Use batteries; if the balance operates correctly, change the adapter. If the problem remains, contact to the Nahita Technical Assistance Department
The displayed value is wrong	The balance is not calibrated	Calibrate the balance
	The balance has not been tared before weighing	Tare the balance before each weighing
The balance does not calibrate	The weight used for calibration is not the correct one	Use the correct weight
	Zero is not adjusted	See Tare function

Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita digital balance, series 5041. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

Please bear in mind the following:

- ◆ This manual is inseparable from the Nahita digital balance series 5041, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the balance avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the balance to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the balance.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to assistance@auxilab.es or by telephone: +34 807 117 040 (0.35 Euros/min).
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the equipment warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the warranty.
- ◆ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Please do not forget filling the warranty certificate and send it before 15 days after the date of purchase.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

 ¡ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

INDEX OF LANGUAGES

English 11-18
 Spanish 2-10
 French 19-27

INDEX OF CONTENTS


1. USES OF THE INSTRUMENT 12
 2. DESCRIPTION 12
 3. TECHNICAL SPECIFICATIONS 13
 4. INSTALLATION / SETTING UP 13
 5. FUNCTIONS OF THE BALANCE 15
 6. MAINTENANCE AND CLEANING 16
 7. TROUBLESHOOTING 18
 ANNEX I: CE CERTIFICATE 28

1. USES OF THE INSTRUMENT

Nahita digital balances are essential instruments at any laboratory when weighing tasks are needed to be carried out. Developed with the most advanced technology they are reliable and durable instruments, with a solid, clean and ergonomic design suitable for industrial applications.

2. DESCRIPTION

1. PAN
2. DIGITAL DISPLAY
3. ON/OFF
4. TARE
5. MODE: to select the weighing unit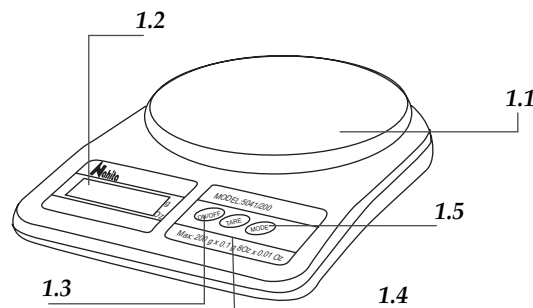

- ◆ Neither submerge the balance nor spill liquids on it.
- ◆ If any liquid comes into contact with the electric parts of the balance please do immediately disconnect it from the net and send it to the Technical Assistance Department, as soon as possible, for its checking and adjustment.
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.

Cleaning

- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the balance and produce an early ageing of the equipment.
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives

 ¡ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

NOTE: According to the in force legislation regarding “Non-automatic weighting instruments” in which balances are included, by means of writ dating from 22nd October, 1994 (BOE 1/3rd/95), these balances must not be used for:

- Comercial transactions.
- Calculating taxes, tariffs, rates, indemnities and other similar canons.
- Judicial surveys.
- Pharmaceutical medicine preparations, as well as analysis made in medical or pharmaceutical laboratories.
- Determining the price or total amount in sale price or in pre-packaged preparations.

INSTRUCTIONS ON ENVIRONMENT PROTECTION

Nahita digital balances series 5041 can operate with batteries; once they are run down, please always deposit them in special containers used for this purpose or in a collection point. Batteries contain heavy metals (zinc, lithium, mercury, etc.) that could contaminate water, soils, etc.

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

After calibration, the balance is ready to be used.

Security

- ◆ The balance must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ◆ Put the balance on top of a horizontal, plane and stable table making a free space of at least 30 cm per side.
- ◆ Do not put the balance near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc. Avoid vibrations, dust and dry environments.
- ◆ When you are not using the balance for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ◆ Never try to repair the balance by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the equipment (see Maintenance).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

6. MAINTENANCE AND CLEANING

To get the best results and a higher duration of the balance it is essential to follow the processes of use.

Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ Prevent the balance from sudden movements and knocks, as well as from direct sunlight or air flows. The balance is a precision instrument, you must handle it carefully.
- ◆ Place the balance in a location with adequate ventilation to prevent heat accumulation. Never use the balance in a wedged location as for example a shelf.
- ◆ Optionally, the balance is supplied with an adapter. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.
- ◆ Never unplug the adapter by pulling the wire, do it from the base.
- ◆ If you use batteries, remove them when you are not using the balance for a long period of time.
- ◆ Never use sharp objects as pens, etc... to press the buttons of the control panel; only use your fingers.
- ◆ Never place on the pan an object heavier than the maximum capacity of the balance, the sensor could be damaged.

3. TECHNICAL SPECIFICATIONS

Among the main features of Nahita digital balances series 5041 we point out the following:

- ◆ Rapid starting, weighing and stabilization
- ◆ Easy-to-use and easy-to-read LCD display
- ◆ Tare function
- ◆ Weighing in grams and ounces
- ◆ Hermetic keyboard to avoid liquid entrance
- ◆ Electrical current connection by adapter (optional) or operation with batteries (not included)
- ◆ Automatic disconnection after two minutes of no use

MODEL	55041220	55041250	55041320	55041350
Capacity	200 g	500g	2000 g	5000 g
Readability	0.1 g	0.1 g	1 g	2 g
Pan diameter	145 mm			
Power source	Batteries (4 x 1.5 V AA) or adapter 220 V (optional)			

4. INSTALLATION / SETTING UP

Preliminary inspection

Unwrap the balance, take off the involving plastic and take off the polispán protection in which it comes fitted. Take off all the protective items and, without connecting the balance to the net, make sure that it does not present any damage because of the shipment. In case the balance presents any damage tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.

Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.

Please check that all the accessories are enclosed with the equipment:

- Instructions manual
- Warranty

We will only accept any equipment return within 15 days after delivery and provided it comes in its original wrapping.

Installation

Before using the balance, it is convenient for you to familiarize with its components and basic essentials, as well as with its control functions.

PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT WITH THE AIM OF ACHIEVING THE HIGHEST FEATURES AND THE MAXIMUM DURABILITY OF THE INSTRUMENT.

Please put the balance on top of a horizontal, plane and stable table making a free space of at least 30 cm per side. Do not put the balance near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc.

Nahita digital balances series 5041 can operate with batteries or by connecting them to the electrical current by an AC adapter.

AC adapter installation

Connect the adapter into the jack on the back side of the balance. Then, plug the adapter into the electrical current provided with earth.

Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to people suffered by the improper use of the electrical connection. The tension should be 220 V, 50 Hz \pm 10%.

BATTERY INSTALLATION

- ◆ Put the balance upside down and remove the lid of the battery compartment by pressing up the tab.
- ◆ Place four 1.5 V batteries type AA inside the compartment bearing in mind to put them correctly according to their polarity.
- ◆ Replace the lid of the compartment

SETTING UP

1. Press the ON/OFF button (1.3) to switch the balance on.
2. When "0.0" is displayed, the balance is ready to weight.
3. By pressing the ON/OFF button (1.3) the balance will switch off. After two minutes (approx.) of no use, the balance will automatically switch off.

By pressing the MODE button (1.5) select the desired weighing unit (grams or ounces).

Abbreviation	Name	Conversion
g	gram	1 g
Oz	ounce	28.3495231 g

5. FUNCTIONS OF THE BALANCE

Tare function

Tare function is used when a substance contained in a recipient is needed to be weight; for example liquids or solid reagents contained in a beaker to prepare a solution.

Switch the balance on and put the empty recipient on the pan. Then press the TARE button (1.4) and "0.0" will be displayed. The symbol "▲" will appear at the lower left side of the display indicating that the balance is tared. The balance is ready to weight.

NOTE: The balance presents a tare range the same as its maximum capacity. This maximum must be taken into account since if the maximum capacity is tared nothing more could be weighted. (For example: if 450 g are tared in a balance with a maximum capacity of 500 g, only 50 g more will be weighted).

- ◆ To extend its life-span, the balance automatically switches off after two minutes of no use.
- ◆ When you are not using the balance for a long period of time, make sure it is disconnected from the net or in case of using the batteries remove them from the corresponding compartment.

Calibration

- ◆ Before starting using the balance and whenever observing important variations in weighing or not correct weighings, balance should be calibrated. Some factors such as variation of gravity depending on latitude, change of working place as well as an inappropriate use of the equipment lead to the necessity of recalibrate the balance.

!Important!! Do not move the balance during calibration

- a. Make sure that there is nothing on the pan (1.1) and check that the balance is switched off.
- b. Press the ON/OFF (1.3) + MODE (1.5) buttons simultaneously for a few seconds until four digits with a minus sign are displayed (e.g. -3029).
- c. Press the ON/OFF button (1.3)
- d. On display, "Zero" will be visualized (for a few seconds) and then the weight that should be used to perform calibration.
- e. Put the indicated calibration weight on the pan.
- f. Press the ON/OFF button (1.3). "Cal" will be displayed and then "PASS"; after a few seconds, and depending on the model, "PASS" or the weight used for calibration will be displayed. Then, remove the weight and switch the balance off by pressing the ON/OFF button (1.3).
- g. The balance is calibrated.

Code	Calibration weight (Not included)
55041220	100 g
55041250	200 g
55041320	2000 g
55041350	5000 g