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Thanks for choosing a NAHITA Precision Electronic Balance. Its high mechanical quality will allow you to make use of this equipment for a long time. Nevertheless, to get the best results and a longer life out of this instrument, it is highly advisable to read this manual and follow the enclosed instructions.

BEFORE CONNECTING OR OPERATING WITH THIS INSTRUMENT, READ THESE INSTRUCTIONS THOROUGHLY.

In order to use this equipment, bear in mind the following points:

- ♦ This manual is an integral part of the balance and must be available to all users of this equipment.
- ♦ As it is a precision instrument, it should be handled carefully avoiding sudden movements and knocks.
- ♦ Never dismantle the different pieces of the balance to repair them yourself. In addition to losing the warranty it could lead to a faulty performance of the equipment.
- ♦ In order to prevent fire or electrical unloads, avoid atmospheres excessive humidity.
- ♦ Any doubts can be clarified by contacting the NAHITA Technical Service, by e-mail asistencia@auxilab.es, by telephone to the 948310513 or by fax 948310500 - 948312071.

ATTENTION !!

IF AN EQUIPMENT IS NOT PROPERLY CLEAN AND DISINFECTED, IT WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

INDEX OF LANGUAGES

Spanish	2-7
English	8-13
French	14-19

INDEX OF CONTENTS

1. USES OF THE INSTRUMENT	8
2. DESCRIPTION	9
3. TECHNICAL SPECIFICATIONS	9
4. OPERATING INSTRUCTIONS	10
5. FUNCTIONS OF THE BALANCE	11
6. MAINTENANCE	12
7. TROUBLESHOOTING	13

1. USES OF THE INSTRUMENT

The Nahita Electronic Balance is an essential instrument in laboratory work when weighing tasks need to be carried out. It is a reliable and long-lasting instrument, which has been developed with the latest technology. Its simple, sturdy, ergonomic design makes this balance suitable for any industrial use, and its smooth surface makes the cleaning easier.

Thanks to its piece-counting function, it can be used in a variety of fields ranging from education or laboratories, to pharmacy, chemistry as well as many others.

7. TROUBLESHOOTING

PROBLEM	CAUSE	SOLUTION
- Digits do not appear on the display.	- Power adapter not connected. - The fuse does not work.	- Connect power adapter . - Replace the fuse.
- Incorrect weight reading.	- Balance was not rezeroed before weighing. - Balance has not been properly calibrated	- Press T with no weight on the pan, then weigh the item. - Recalibrate correctly.
- Calibration procedure does not work.	- Incorrect weights being used - Zero adjustment may be incorrect.	- Use correct weights. - See taring.

Note: According to the in force legislation regarding *Non-automatic weighing instruments* in which balances are included, by means of writ dating from 22nd October, 1994 (BOE 3/1/95), these balances must not be used for:

- Commercial Transactions.
- Calculating taxes, tariffs, rates, indemnities and other similar canons.
- Judicial surveys.
- Pharmaceutical medicine preparations, as well as analysis made in medical or pharmaceutical laboratories.
- Determining the price or total amount in sale price or in prepackaged preparations.

Counting function


- ◆ This function allows the balance to calculate and display the total number of pieces, as well as the total weight of the pieces. In piece-counting, the balance assumes a reasonably uniform weight among the parts.
- ◆ To exit the piece-counting function, press *N* again. The display will now show *Count Off*. Then, wait until the desired weighing unit is displayed. The balance will come back to its normal weighing mode.
- ◆ Make sure there is nothing on the pan, press button "*T*" and the display shows *0.00 g*.
- ◆ Place 10 items on the pan.
- ◆ Press Counting (*N*), the display will show "*c 1.0*". Remove the pieces. The balance is now ready to start weighing.
- ◆ Place the whole amount of pieces that need to be counted. The display will show the total amount .
- ◆ To exit the piece-counting function, press "*N*" again.
- ◆ The balance is now ready to start weighing .

6. MAINTENANCE

- ◆ Preserve the balance from heavy shocks. Look for a dry place, away from dust, direct sun beams or air flows. Since it is a precision instrument, it should be handled carefully.
- ◆ Never remove any of its mechanical components. In addition to losing the warranty, it could also lead to a faulty performance of the balance.
- ◆ Save all packing material until you are sure that the instrument works correctly. All defective items must be returned in the original packing together with the supplied accessories.
- ◆ It is prohibited to use strong chemical solvents to clean the surface of the balance. Don't dip the balance into water, or leave it out in the rain. In case it needs cleaning, don't wash it with water but rather use a cloth damped with a mild detergent.
- ◆ Always leave space around the balance for its correct ventilation.
- ◆ You have this manual always by hand, so that any person can consult it.
- ◆ Follow the instructions and warnings relative to this manual.
- ◆ Do not press the display or any key with a pointed object such as a ballpoint pen or a screwdriver, as this may scratch or damage the key.
- ◆ It does not cause vibrations when this weighing.
- ◆ It is prohibited to use strong chemical solvents to clean the surface of the balance. Don't dip the balance into water, or leave it out in the rain. In case it needs cleaning, don't wash it with water but rather use a cloth damped with a mild detergent.
- ◆ Never remove any of its mechanical components. In addition to losing the warranty, it could also lead to a faulty performance of the balance.
- ◆ Do not throw away the packing material. It might be useful in case you need to move it, to send it back for repairing or for long-term storage.

2. DESCRIPTION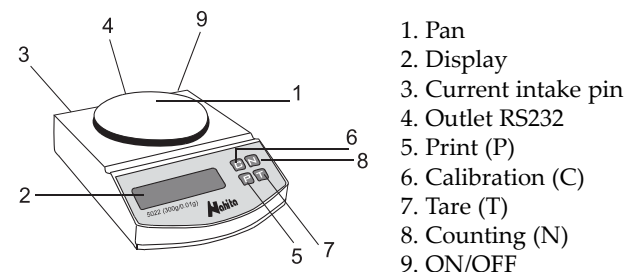

3. TECHNICAL SPECIFICATIONS

Among the main features, we point out the following:

- ◆ External metallic structure, stainless steel weighing pan and hermetic keyboard to avoid liquids entrance.
- ◆ Stable and rapid weighing.
- ◆ Clear reading through the LCD display, with lightweight.
- ◆ Subtractive tare.
- ◆ Automatic calibration.
- ◆ Alternative weighing modes: grams, ounces, carat.
- ◆ Parts counting function.
- ◆ It Includes calibration weight.
- ◆ It includes weighing pan of 130 mm Ø.
- ◆ External AC 220V power supply adapter included.
- ◆ Automatic disconnection after five minutes without using it.
- ◆ RS 232 interface.

Code	55022220	55022230	55022250
Capacity	200 g	300 g	500 g
Readability	0.01 g	0.01 g	0.01 g
Tare range	200 g	300 g	500 g
Pan size	130 mm Ø		
Stabilization time	3 seconds		
Working temperature	5° C - 40° C		
Dimensions	215 x 184 x 50 mm		
Weight	1 Kg		
Power source	220-240V, 50-60Hz (AC)		

4. OPERATING INSTRUCTIONS

Preliminary examination

- ◆ Remove the balance from the packing material and examine it carefully to make sure no damage has occurred during shipping. If there is any damage, notify to your dealer/forwarding agency.
- ◆ Keep all packing material until you are sure that the instrument works correctly. All defective items must be returned in the original packing together with any other supplied accessories.
- ◆ Verify you have received all the accessories that come with the balance:
 - Pan
 - Weight for calibration
 - Cable
 - 2 Fuse 0.5 A
 - Instruction manual
 - Warranty

Installation

BEFORE CONNECTING OR OPERATING WITH THIS INSTRUMENT, READ THESE INSTRUCTIONS THOROUGHLY.

- ◆ The instrument must be placed on a flat, steady, horizontal desk. Keep it away from humidity, heat, vibrations or airflows. The natural light must not be too bright. Keep it away from any magnetic force or electric field or gadget which can produce a high frequency field.
- ◆ The balance is provided with a plug. This plug must be connected into an outlet which is connected to ground. The outlet must be at hand so as to disconnect it in case of an emergency.

Neither the manufacturer nor the supplier will assume any responsibility for damages caused to the balance, or for any other damages suffered by any person due to the neglect of the correct electrical connection process.

Starting the balance

- ◆ With no load on the pan, turn the power switch on. Press the ON/OFF key (9). When the balance is first switched on, all digits on the display will light up. After a few seconds, it will show 88888, E- 1/ E- 9 and finally, the display will show 0.00 g.
- ◆ Stop to stabilize the balance during 10 minutes.
- ◆ Select the desired weighing unit by pressing "T". Available units: grams (g), ounce (oz), carat (quilate):
 - Press the button counting "N" and it will show "-----".
 - Press "T" and choose the unit. grams ("-----"), carate ("C") ounces("---0---").
 - Press "N" and show display 0.00 g (grams), 0.00 C(carat) and 0.0000 o (ounces).

UNITS CONVERSION

Abbreviation	Name	Conversion
g	Gramme	1 g
oz	Ounze	28.3495231 g
ct	Carat	0.2 g

5. FUNCTIONS OF THE BALANCE

Taring

- ◆ Taring enables the balance to automatically subtract the weight of the container, when the material to be weighed must be held in a container (such is the case of liquids). Proceed as follows:
 - ◆ Make sure there is nothing on the pan.
 - ◆ Put the recipient on the balance pan, press button "T", in display show 0.00 g. If you select another weighing unit, 0 will be relative to the selected unit. The balance is now ready to start weighing.
 - ◆ Place the material to be weighed into the container. When the material is added, its net weight will be displayed.
 - ◆ When the container and its contents are removed from the pan, the tared weight (the container's weight) will be displayed as a negative number.
 - ◆ To recover the balance's usual mode, press "T" again.
 - ◆ The combined weight of the container and the material must not exceed the capacity of the balance.

Calibration

- ◆ The balance has been calibrated before shipment. However, it should be checked before use and recalibrated if necessary. Calibration could be influenced by factors such as:
 - Variations in the earth gravitational field at different latitudes of the world.
 - Rough handling.
 - Changes in work location.

IMPORTANT!
DO NOT MOVE THE BALANCE DURING THE PROCESS OF CALIBRATION

- ◆ From time to time it is advisable to calibrate the balance according to the performance it is going to give:
 - ◆ Switch the balance on and once the pan is empty, press on button "T". You will read value 0.00 g from the display, then, press key "C" and letter "C" will appear. If "CE" is read, switch it off and repeat the calibration process once 0.00 g has appeared.
 - ◆ Place the calibration weight taking into account the balance's model, it should be 200 g for balances of a 200 and 300 g capacity, and 500 g for 500 g, balances. Wait and check the screen will show the calibration's weight and listen to the pib indicating the balance is already calibrated.
 - ◆ Remove the calibration weight and the balance will automatically recover the usual weighing mode. The display will turn into 0.00 g.
 - ◆ Once the calibration is finished, the balance is already prepared to start working.