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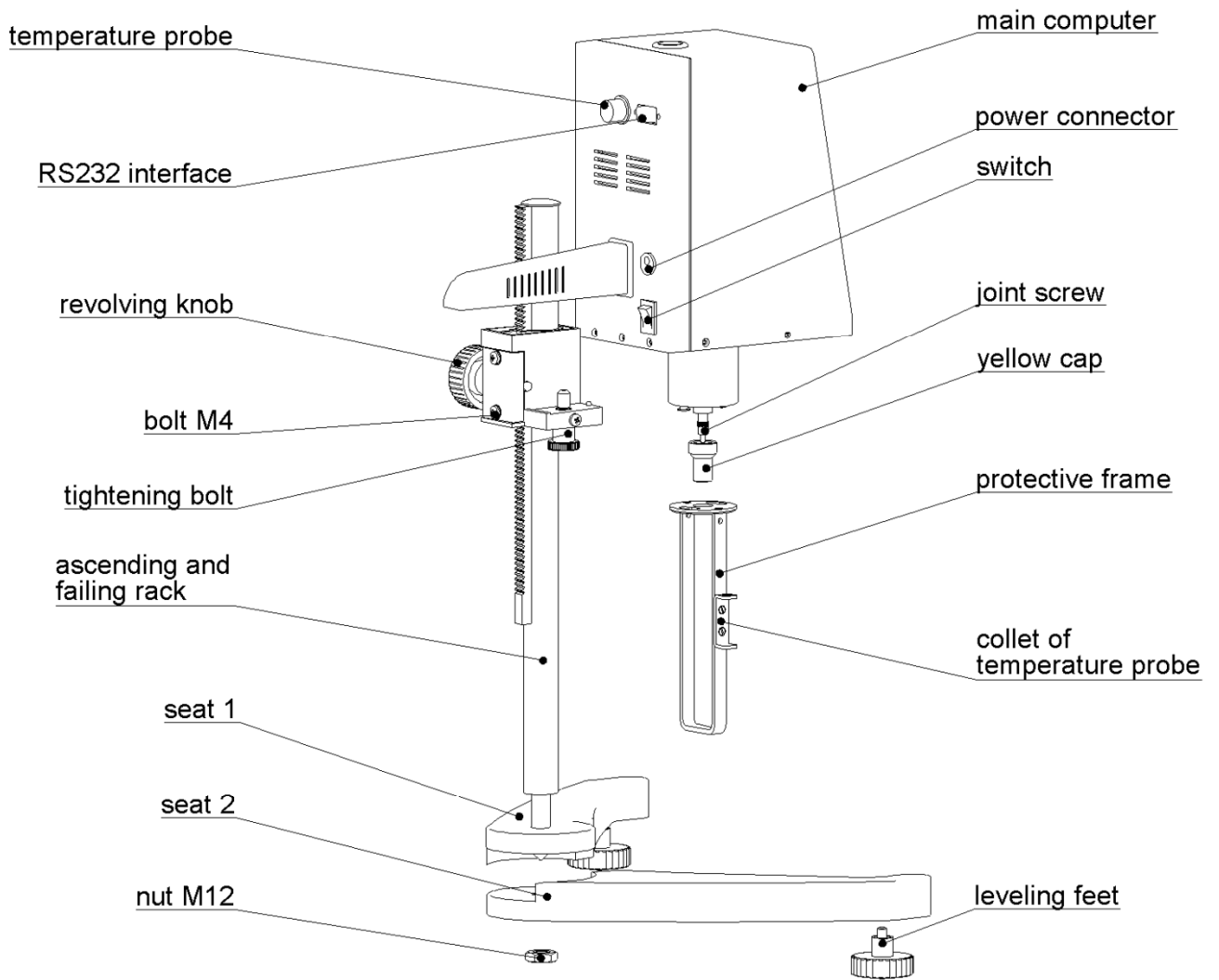
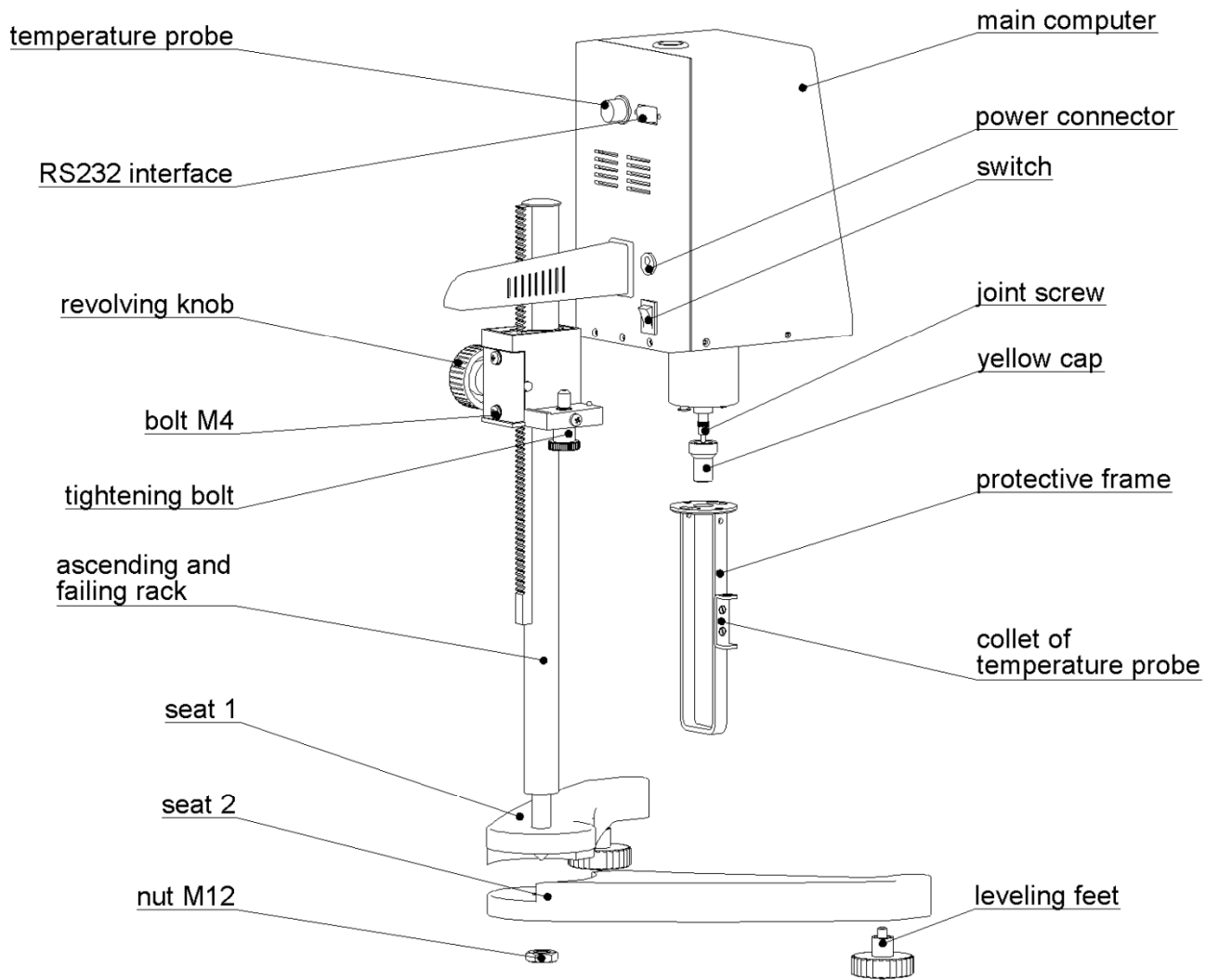
**USER'S MANUAL  
DIGITAL VISCOSIMETERS  
SERIES 802/807**


Nahita digital viscosimeters are equipped with a high subsection driven step motor, a 16 bit micro-computer controlled processor and a back-lit LCD display. The viscosimeter is stable and accurate in motion, definitude in key demonstration, programmable in design, easy for operation. The display directly demonstrates the viscosity, rotating speed, rotor number and the maximum viscosity measured for the selected rotor and the current rotating speed. The main controlling board and subsection driven board are all manufactured by adopting the Surface Mount Technology (SMT). The circuit is adopted by the micro processor that is most advanced at present, with compact and reasonable structure. Provided with R232, the space for printing can be set up by the user. The full range and linearity at points are adjusted by PC interface. Its performance and functions have reached to the advanced level compared with other similar products.

Nahita digital viscosimeters are used in checking the viscous resistance and dynamical viscosity of liquid. They are widely used in measuring the viscosity of various liquid such as grease, painting, foodstuff, dope, paper making, cosmetics, chemical industry, capsule stickiness agent and medicines.

# Mail Installing Chart

## Rotor Kits

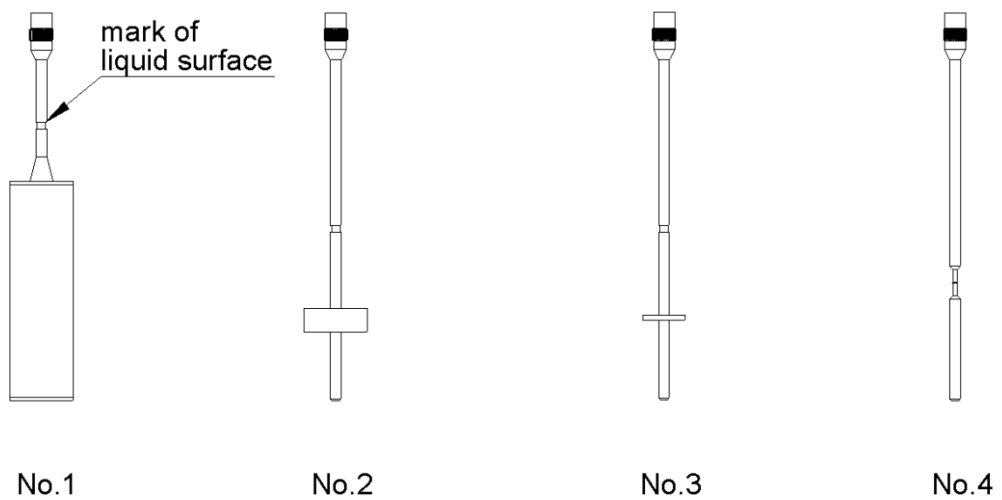
## Technical specifications

Model	802	807/1	807/2
Measuring range ( mPa·s )	1-100000	1—600000	1—6000000
Rotating speed (RPM)	6, 12, 30 o 60	1—60	0.1—99.9
Rotor number	Rotor N <sup>o</sup> .1,2,3,4 (supplied with equipment)		
Accuracy	±2.0%		
Return capacity	0.5%		
Power	Adapter (input 220 V; 50±1 Hz; output 15 V, 1.2 A)		

## 2. Structure and principle of operation

As shown in the picture, the pointer of the load cell driven by the high subsection driven step-motor is driving the rotor to go round and round through the hairspring and rotating axis. If the rotor is not resisted by the liquid, the pointer of the hairspring load cell is in the same position with the pointer of the load cell for the step-motor. On contrary, if the rotor is being resisted by viscous liquid, the hairspring produces the wresting square and repel and balance with the viscous resistance for reaching the balance at large. At this moment, by transmitting the output signal by the photoelectrical load cell to 16 bit micro processor for data processing, the viscous value (mPa·s) will be displayed on back-lit LCD screen.
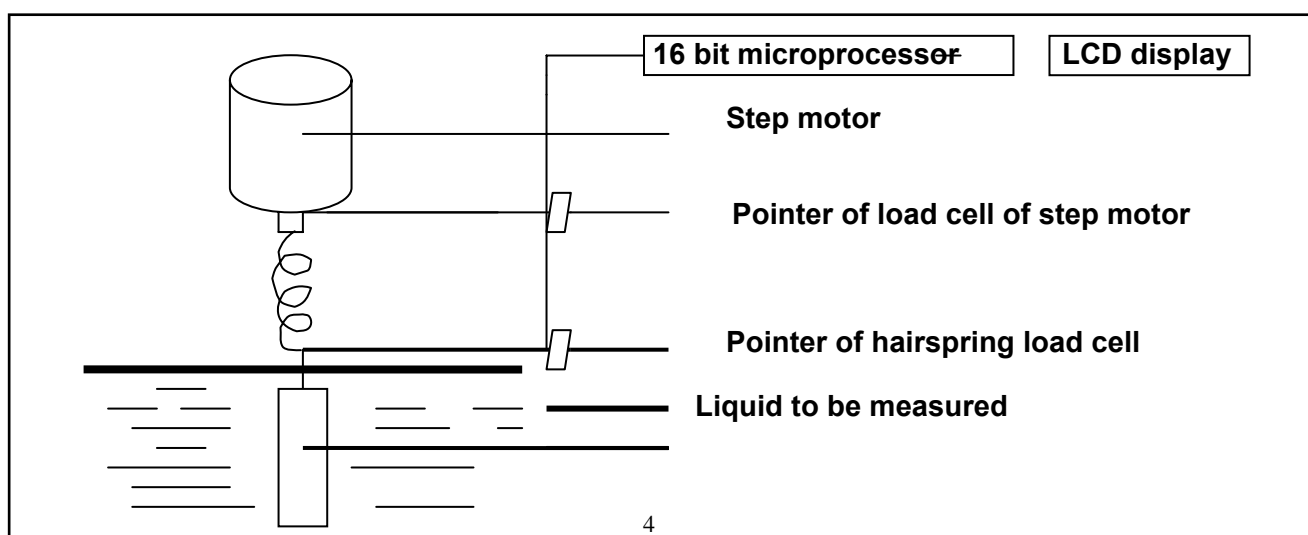
## **Rotor**

### **3. Supplied items**

- (1) Digital viscosimeter
- (2) Ascending and falling rack
- (3) Base
- (4) Rotors ( No.1,2,3,4)
- (5) Rotor protective fram
- (6) Temperature probe (Series 807)
- (7) Power adapter
- (8) Spanner
- (9) Screw driver

### **4. Installation (Please refer to the installing chart)**

- (1) Take the base, the ascending and falling rack, the main body of viscosimeter, the rotors and the protective frame out from the case.
- (2) Put the ascending and falling rack into the hole of both parts of the base so as the opening direction of the base and the upper rack face user; the tighten firmly the M12 nut with the supplied spanner.
- (3) By turning the knob of the ascending and falling rack check that the clamp moves up and down easily; if movement is too loosen or tighten, then adjust the M4 screw at the back by using the screw driver supplied. It is recommended that the screw is firmly tighten so as avoiding the viscosimeter to fall down during operation
- (4) Install the viscosimeter body on the clamp of the ascending and falling rack and hold it firmly with the tightening bolt. Check that the rotor joint screw is well protected with the yellow cap (always put this cap in case of transporting the equipment or if it is not going to be used for a long time).
- (5) Adjust the two leveling feet so as the equipment is well leveled (the level bubble at the upper part of equipment must be well centered).
- (6) Connect the temperatura probe.
- (7) Make sure that the switch is in off position and connect the equipment to the main supply through the adapter.

## 5. Function of keys

**Reset:** Stop measuring / Reset the applied program

**Print:** Printing function

**Run:** Start measuring

**Spindle:** Rotor selection (maximum viscosity value measured depending on the selected speed)

**Speed:** selection of rotating speed

*Series 807*

↑: Increase speed value

↓: Decrease speed value

→: Go to next digit (during speed selection)

←: Back to previous digit (during speed selection)

## 6. Description of displayed data

*Series 807*

<b>20.0 °C</b>	<b>60.0 RPM</b>	<b>45.7%</b>
		<b>SP 4</b>
	<b>888888.8 mPa·s</b>	

**20.0 °C:** Temperature measured by the temperature probe in °C

**SP:** Number of selected rotor (e.g. SP 4 means selection of rotor nº 4)

**60.0 RPM:** Rotating speed of motor in RPM

**888888.8 mPa·s:** Maximum viscosity measured at the selected rotating speed and with the selected rotor in mPa·s.

**45.7 %:** Percentage of measurement or wetting square. Its value must be between 20% and 90%; if it is out of this range, it will be necessary to change the rotating speed and the rotor since the selected ones are not adequate for the viscosity of the liquid to be analyzed.

## 7. Operation mode

- (1) Install the equipment according to *Installation chapter* instructions and put the viscosimeter at an proper place.
- (2) Put the rotor protective frame (turn it rightwards to put and leftwards to remove)
- (3) Install the rotor in the joint screw (turn it leftwards to put and rightwards to remove); previously, the yellow cap must be removed.
- (4) Turn the equipment on by pressing the on/off switch; the motor will start running.
- (5) Select the rotor number by pressing the rotor selection key (Spindle); every time the key is pressed number will change following this sequence 1→2→3→4→0→. When the desired rotor number is displayed release the selection key.
- (6) Select the rotating speed by pressing the corresponding key (Speed); use keys with arrows to increase or decrease the speed value. Press the rotation speed selection button (Speed) to save the input value.
- (7) By turning the knob of the ascending and falling rack, the viscosimeter will go up or down slowly to let the rotor immerse in the liquid to be measured; the surface of the liquid must be at the level of the mark in the rotor. Check the equipment is well leveled.
- (8) Press the button to start measurement (Run); rotor will start rotating and the viscosity measured at those conditions and the percentage of measurement will be displayed at the same time.
- (9) If rotor is needed to be changed during operation, press directly the stop key (Reset); motor will stop and equipment will be kept on. Once the rotor is changed, you can continue with measurement as indicated in steps 6 and 8 mention above.
- (10) Data printing
  - a. First of all, select the time interval for data printing. For this, after pressing printing key (*Print*) the interval S : 00 : 05 will be displayed. The time interval for printing can be modified by pressing printing key (*Print*) successively, so as the following sequence will be displayed 00 : 05→00 : 10→00 : 20→00 : 30→01 :

00→01 : 30→02 : 00→05 : 00→. Being, S : 00 : 05 data printing every 5 seconds,

00 : 10 data printing every 10 seconds and so on.

b. After connecting the viscosimeter with printer, press printing key (*Print*) once more. The command “on” will be printed on paper and then, data will start to be printed. Press printing key (*Print*) again to stop printing data; the command “Off” will be printed on paper.

## **8. Measurement of a liquid with unknown viscosity**

- (1) General principle of measurement: for samples with high viscosity, chose a small rotor (n° 3 or 4) and a low rotation speed. For samples with high viscosity, choose a big rotor (n° 1 or 2) and a high rotation speed. If during measurement, the percentage of measurement is in a normal value between 20% and 90%, then the measured viscosity of the sample is correct.
- (2) First of all the viscosity range of the sample to be measured must be estimated and then, choose the smallest rotor and low speed for samples with high viscosity or choose the biggest rotor and high speed for samples with low viscosity. In most cases, the rotor should be chosen at first and then the rotating speed. For example, when choosing rotor n° 1 and rotating speed is 60 rpm, the maximum viscosity that can be measured will be 100 mPa·s (shown on display); if rotating speed is changed into 6 rpm, then the maximum viscosity that can be measured will be 1000 mPa·s.
- (3) If the viscosity of the sample cannot be estimated, then the rotor and rotating speed should be chosen in base of the percentage of measurement (wresting square). A percentage in the range between 20% and 90% is a normal value. If percentage is out of this range, an alarm will sound to remind that the speed and rotor must be changed. Never forget to input the correct rotor n° once it has been changed.

## 9. Cautions

- (1) Be careful when installing or uninstalling the rotor; rise softly the joint screw and never force it. Never put the rotor to receive the force horizontally to prevent it from being bended.
- (2) Never lie down or tilt the viscosimeter once it is mounted and installed.
- (3) Always keep clean the joint screw and the nut of the rotor; otherwise, measurements and correct operation could be affected.
- (4) Hold the ascending and falling rack with the hand every time the viscosimeter is moved up or down to avoid it from falling due to its weight.
- (5) After changing rotor, input the correct rotor number. Always clean the rotor before installing; never clean the rotor once it is installed in the equipment.
- (6) When changing the sample, clean and dry the rotor and the protective frame to avoid mistakes in measurement due to the mixture of samples.
- (7) Each viscosimeter has been calibrated and adjusted with its own set of rotors, never interchange rotors between different viscosimeters.
- (8) Never dismantle neither change any spare part of the equipment.
- (9) When moving or transporting the viscosimeter, do not forget to put the yellow cap to protect the joint screw; use the case to transport all the pieces properly protected.
- (10) Once the rotor is installed, do not make it turn for a long time without liquid to avoid it from being damaged.
- (11) In non Newtonian liquids, values of viscosity vary with speed and time so as the results obtained with different rotors and speed can vary. This is normal in this kind of liquids; it is not a problem of the equipment. Usually the rotor, rotating speed and measuring time must be adjusted to each non Newtonian liquid.
- (12) To obtain a high accuracy of results, please, bear in mind the following points:
  - a . More precisely control of the temperature of the liquid to be measured
  - b . Immerse the rotor in the liquid to be measured for an enough long time and keep it always at the same temperature.

- c . Keep the uniformity and homogeneity of the liquid to be measured
- d . To perform measurements, the rotor should be placed at the centre of the recipient and the protective frame should properly placed
- e . Make sure the rotor is perfectly clean
- f . Turn the equipment off when changing from high to low speed or keep it for a longer time at low speed to avoid the error caused by the liquid inertia.
- g . When working with low viscosity, use rotor n° 1 and when working at high viscosity use rotor n° 4.
- h . The measure time for the viscosity measured at low speed is relatively longer.
- i . Raise the viscosimeter by turning the knob of the ascending and falling rack every time the rotor or sample is going to be changed.

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