AUXILAB S.L.

DECLARATION OF CONFORMITY
NAHITA CENTRIFUGE 2924 of Auxilab, S.L for the Directive
of Machines (89/392/CEE modified) and the regulations
adopted for their transposition

NAME OF THE MANUFACTURER / IMPORTER:
AUXILAB, S.L.

ADDRESS: **Polígono Morea Norte, 8**
31191 Beriáin (Navarra)

WE STATE THAT:

NAHITA ROTARY EVAPORATOR MODEL 503
CODE 59000503

Are designed and manufactured according to:

- ◆ Directive 89/392/CEE, including the modifications and national regulations that transpose them.
- ◆ Directive 89/336/CEE modified over the electromagnetic compatibility.
- ◆ Directive 73/23/CEE modified over the electric security.

And that the following harmonized rules have applied (or part of them):

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BERIAIN 11th November 2006

Signed by: **ALFONSO AINCIBURU SANZ**
DIRECTOR/MANAGER

Polígono Morea Norte, 8 31191 Beriain (Navarra) - Spain. Tel. 948 310 513 Fax 948 312 071
Internet: www.auxilab.es · Email: correo@auxilab.es

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This manual should be available for all users of these equipments. To get the best results and a higher duration of this equipment it is advisable to read carefully this manual and follow the processes of use.

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- ◆ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the equipment (see Maintenance).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

5. MAINTENANCE AND CLEANING

To get the best results and a higher duration of the rotary evaporator it is essential to follow the processes of use.

Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

- ◆ Please follow the processes of use of this manual.
- ◆ This manual should be available for all users of this equipment.
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- ◆ The rotary evaporator is supplied with an adapter. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.
- ◆ Never try to repair the rotary evaporator by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the rotary evaporator (burns, hurts...) or damages in nearby equipments.
- ◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.

Cleaning

- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the rotary evaporator and produce an early ageing of the equipment.
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

INSTRUCTIONS ON ENVIRONMENT PROTECTION

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

Remember: grease or spread with vaseline the ground joints of all the glass pieces to avoid breakage when disassembling.

Fill the water bath.

Note: never connect the bath without water inside, the resistance will break down.

The Nahita rotary evaporator presents the possibility of turning the glass assembly into an angle of 0-45° by a screw that acts as a turning device. Put the glass assembly into the desired angle and fix such position with the screw.

Once the evaporating system is ready, connect the equipment to the power supply.

The rotary evaporator is supplied together with a Schuko standard wire. Please insert the wire that feeds the AC electric current in the base of current 220V 50Hz ±10% provided with earth wire and to the other end to the rotary evaporator connector.

Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to people suffered by the improper use of the electric connection. The tension should be 220V 50Hz ± 10%.

7. Select the desired temperature of the water bath by the analogical adjustment knob.

8. Press the power ON.

9. Adjust the turning speed of the evaporating flask by the analogical adjustment knob.

NOTE: When turning the equipment on, put the bath cover to avoid spilling when working at high temperatures.

10. Connect the condenser to the water flux. The condenser presents two connections that must be respectively connected to the water inlet and outlet hoses.

11. The vacuum connection of the system is situated at the upper part of the condenser; it must be connected to a water jet pump or a vacuum pump.

NOTE: In case you have to send the equipment for its repair, DO NOT send the glass components since due to their fragility they can be damaged. We do not take responsibility for the broken pieces that arrive to our Technical Assistance Department.

Security

- ◆ The rotary evaporator must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ◆ Put the rotary evaporator on top of a horizontal, plane and stable table making a free space at least at 30 cm per side.
- ◆ Do not put the rotary evaporator near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc. Avoid vibrations, dust and dry environments.
- ◆ Durante su funcionamiento el material peligroso como líquidos inflamables o material patológico, deben estar fuera del área de trabajo.
- ◆ When you are not using the rotary evaporator for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ◆ Never try to repair the rotary evaporator by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).

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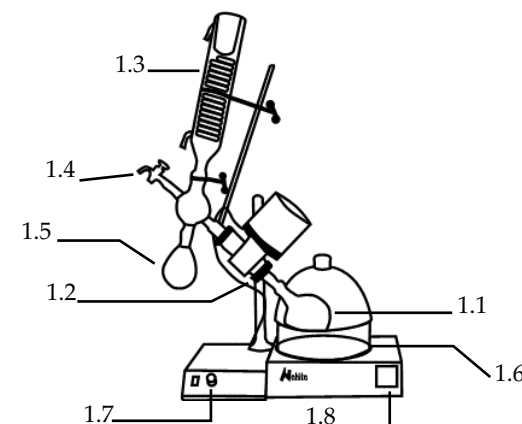
PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT WITH THE AIM OF ACHIEVING THE HIGHEST FEATURES AND THE MAXIMUM DURABILITY OF THE INSTRUMENT.

Please put the rotary evaporator on top of a horizontal, plane and stable table making a free space at least at 30 cm per side. Do not put the equipment near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc.

During operation, dangerous material such as flammable or infectious liquids must be out of the working area.

When you are not using the rotary evaporator for a long period of time please make sure it is unplugged and protected from dust, in order to avoid possible accidents and extend its life-span.

1. Assemble the evaporating system (see drawing)

Previously, put the rod on the motor support with the four supplied screws. Place the motor on its corresponding bosshead; you have received a fixing screw, insert it bearing in mind that it presents two slots that must fit with the rod.

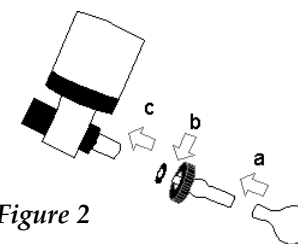
Figure 2

2. Put the connector (white Teflon piece) into the fixed evaporating system by introducing the narrower part of the connector inside the system. Screw the metal ring and the black piece of the piece to fix the evaporating flask, in the ground joint adapter neck. Finally, join this set with the white Teflon piece previously inserted into the system and use the black piece to fix all the components (see figure 2).

3. To assemble the left part of the rotary evaporator, place the rod with the four supplied screws (figure 3a). Arrange the bossheads (figure 3b) and fixing rings (figure 3c) that will hold the condenser, placing the narrower ring at a lower position and the wider ring at an upper position. Put the joining glass piece in the position where the rest of the glass components (condenser, receiving flask and inlet tube) converge (figure 4).

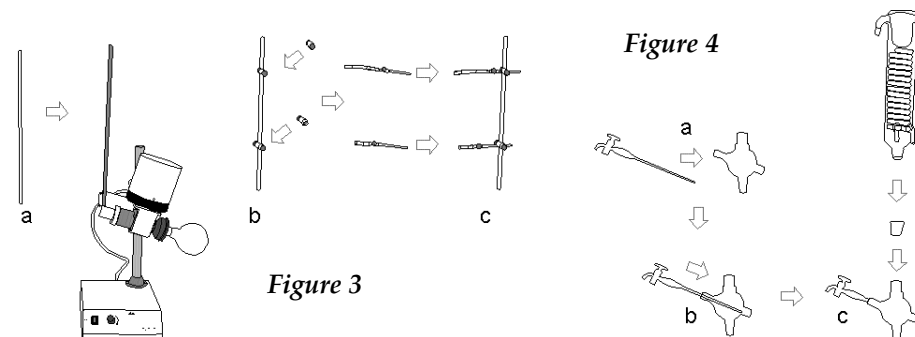
Figure 3

Figure 4

3. TECHNICAL SPECIFICATIONS

Manual raising mechanism	Up to 150 mm
Turning angle of the glass assembly	45°
Speed range	0-90 r.p.m
Speed adjustment	Analogical
Temperature range of the water bath	RT - 100 °C
Speed adjustment	Analogical
Precision	+/- 1 °C
Bath power	1100 W
Rotor power	40 W
Power supply	220-240 V, 50-60 Hz +/- 10%

4. INSTALLATION / SETTING UP

Preliminary inspection

Unwrap the rotary evaporator, take off the involving plastic and take off the polispán protection in which it comes fitted. Take off all protections and without connecting the rotary evaporator to the power supply, make sure that the equipment does not present any damage because of the shipment. In case it presents any damage, tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.

Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.

Please check that all the accessories are enclosed with the equipment:

- Rotary evaporator
- Thermostatic water bath
- 1 Glass condenser (ground joint 29/32)
- 1 Glass joining piece with three necks
- 1 Inlet tube
- 1 Ground joint adapter
- 1 Round bottom flask, 500 mL
- 1 Pear-shaped flask, 250 mL
- 2 Rings, 2 Bossheads
- 2 Plastic clips for ground joints 29/32
- 1 Metal clamp
- 1 Metal piece to fix the evaporating flask
- 1 Rod to hold the glass assembly
- 1 Rod to hold the motor
- Warranty
- Instructions manual

We will only accept any equipment return within 15 days after delivery and provided it comes in its original wrapping.

Installation

Before using the rotary evaporator, it is convenient for you to familiarize with its components and basic essentials, as well as with its control functions.

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Spanish 2-9
 English..... 10-16

INDEX OF CONTENTS

1. USES OF THE INSTRUMENT 11
 2. DESCRIPTION 11
 3. TECHNICAL SPECIFICATIONS 12
 4. INSTALLATION / SETTING UP 12
 5. MAINTENANCE AND CLEANING..... 15
 6. ANNEX I: CE CERTIFICATE 16

1. USES OF THE INSTRUMENT

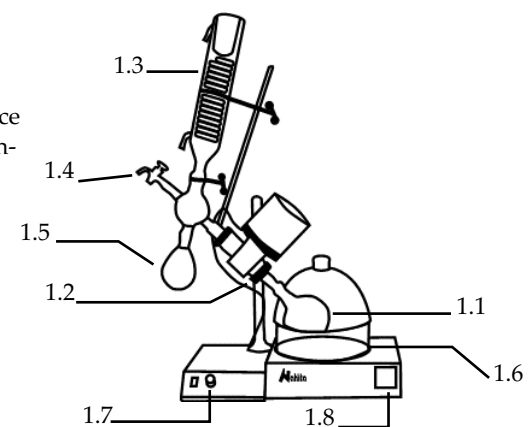
Nahita rotary evaporators are essential instruments in Biochemistry, Chemistry and Pharmacy laboratories to carry out those operations in which rapid distillation or sample concentration is required.

Rotary evaporator is one of the most important laboratory instruments due to its capacity of increase evaporation speed twice as much as that got with traditional methods. This is obtained by two different mechanisms:

- Increase in evaporating surface by continuously heating and turning the flask into the water bath.
- Application of vacuum inside the system.

2. DESCRIPTION

- 1.1 Evaporating flask
- 1.2 Ground joint adapter with a piece to fix the evaporating flask and the connector.
- 1.3 Condenser
- 1.4 Inlet tube
- 1.5 Receiving flask
- 1.6 Thermostatic water bath
- 1.7 Speed adjustment knob
- 1.8 Temperature adjustment knob

Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita rotary evaporator 503. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

- ◆ Please bear in mind the following:
- ◆ This manual is inseparable from the Nahita rotary evaporator 503, so it should be available for all the users of this equipment.
- ◆ You should carefully handle the equipment avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the equipment to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the rotary evaporator.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to asistencia@auxilab.es.
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the equipment warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the warranty.
- ◆ Neither fuses nor accessories (including their loss), are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

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