DUCHAS Y LAVAOJOS DE EMERGENCIA EMERGENCY SHOWER AND EYE WASHER

Nahita

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This manual should be available for all users of these equipments. To get the best results and a higher duration of this equipment it is advisable to read carefully this manual and follow the processes of use.

4. INSTRUCTIONS OF USE

Note: Showers and eye washers must be use only in case of emergency, not for a daily use. Activate the shower and eye washer at least weekly to assure that it will operate properly when needed.

- ◆ To activate the sprays push the lateral lever clockwise; to activate the shower pull the handle downwards.
- ◆ In case of using the eye washer, keep eyes opened with the help of hands so as the damaged zone is abundantly washed with water.
- ◆ To stop water flow, press the lateral lever counter-clockwise /eye washer) or push the handle upwards (shower).

Note: The time of wash must be adequate to the type of damage produced. The emergency showers and eye washers can reduce the degree of damage produced but they are never substitutive of a medical treatment. Apply the appropriate medical treatment.

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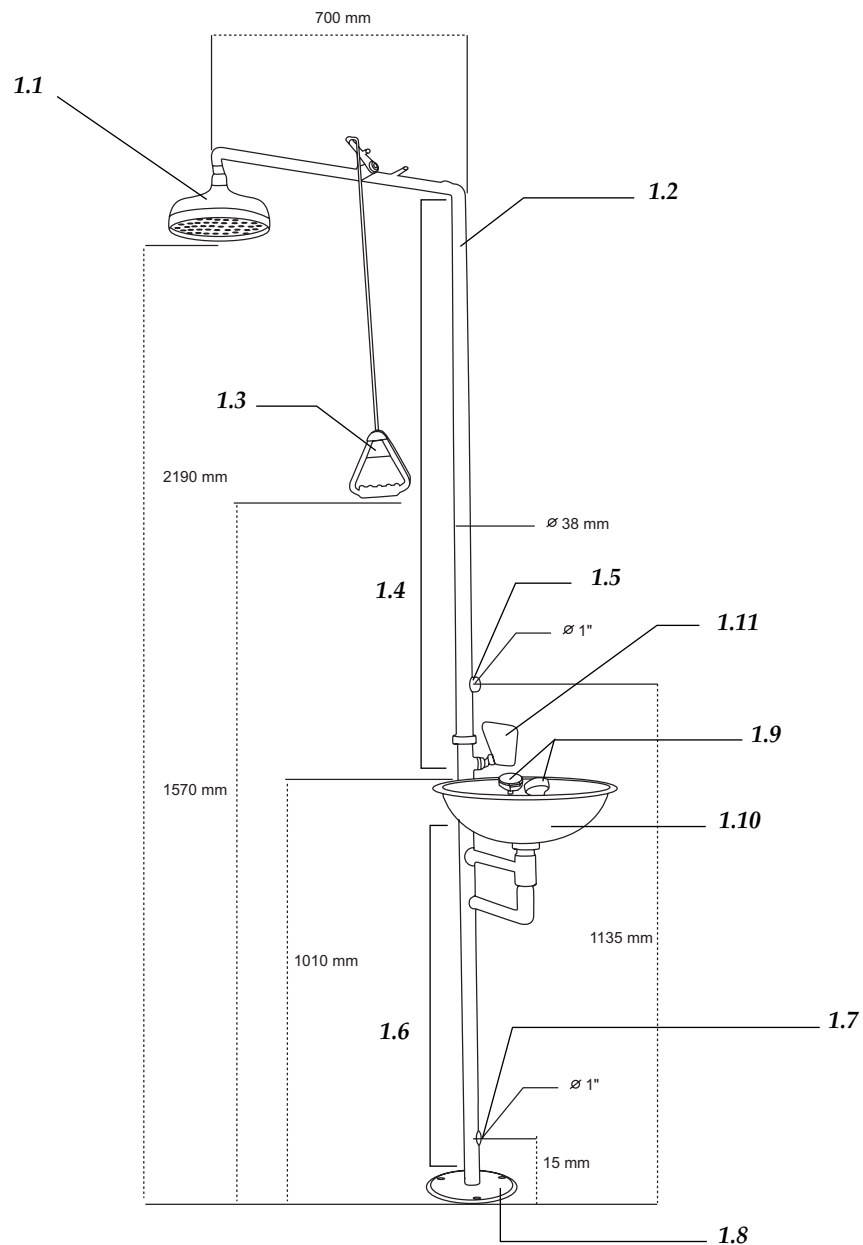
- Tank
- Upper half part with water inlet
- Lower half part with water outlet
- Shower pipe
- Base
- Eye washer starter lever
- Shower starter handle
- User's manual

We will only accept the product return within 15 days after delivery provided it comes in its original wrapping.

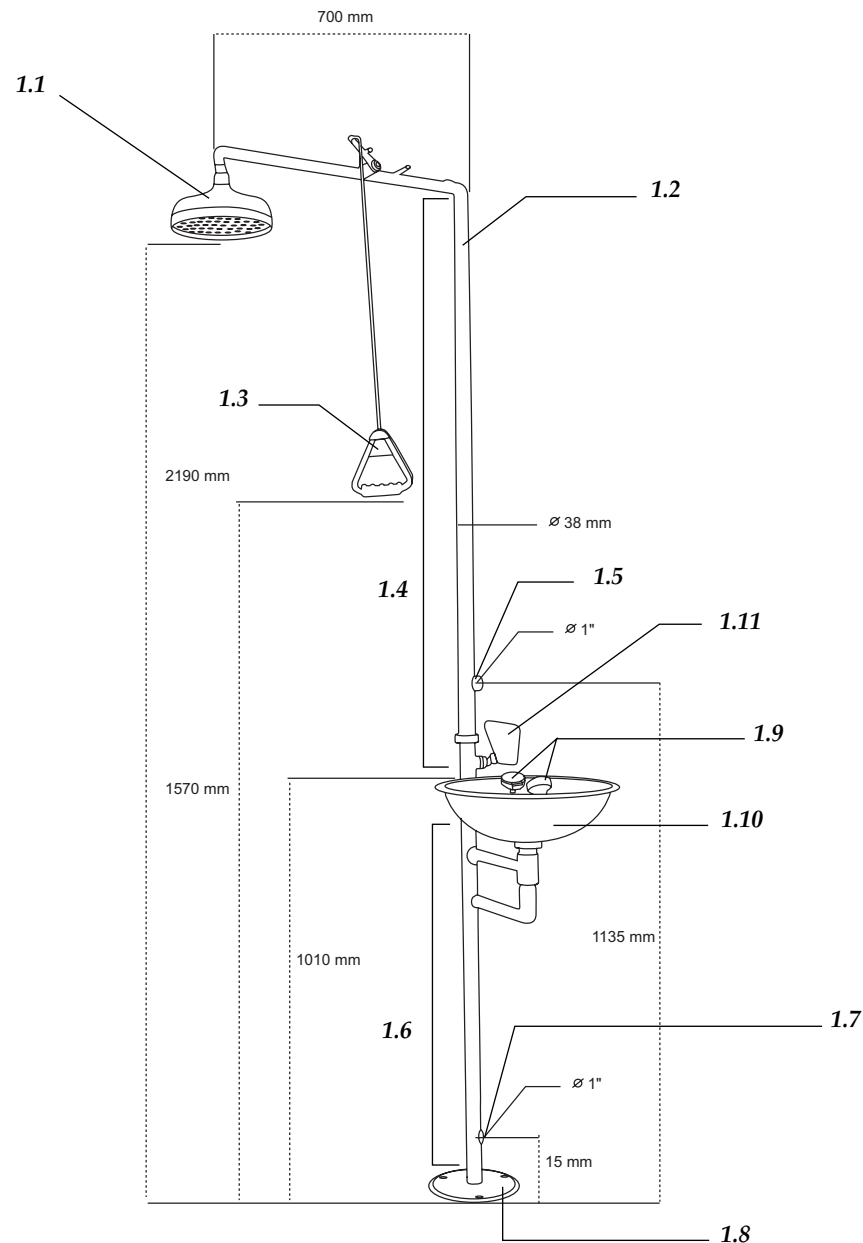
Installation

IMPORTANT: *The place chosen to install the shower and eye washer must be adequate, so as it is perfectly visible and accessible from the region where the risk of accident exists. The distance must not be higher than 15 m or 10 s walking from the zone of risk.*

Wrap all threads with teflon tape or similar to assure all joints are perfectly sealed and avoid water leaks. Then, assemble the different pieces of the shower and eye washer according to the diagram shown:

- ◆ Screw the lower half part of the pipe in the base supplied so as water outlet is downwards.
- ◆ Assemble the upper half part with the shower pipe.
- ◆ Assemble the upper and lower half parts of the pipe.
- ◆ Install the eye washer tank and then the head with two sprays. Put the starter lever so as it opens the valve when being pushed clockwise.
- ◆ Screw the shower head and install the starter handle.
- ◆ Connect the shower and eye washer to the water inlet and outlet.

- 1.1 Shower head
- 1.2 Shower pipe
- 1.3 Shower starter handle
- 1.4 Upper half part
- 1.5 Water inlet
- 1.6 Lower half part
- 1.7 Water outlet
- 1.8 Base
- 1.9 Sprays
- 1.10 Tank
- 1.11 Eye washer starter lever

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1. USES

Emergency showers and eye washers are a necessary security measure with high utility in most of companies and laboratories. They can reduce the damage caused by the splashing in the eyes, face and the body with acids, bases, organic compounds and other toxic or corrosive components, as well as in case of fire. It is the easiest and cheapest way of decontamination in case of accident.

2. TECHNICAL SPECIFICATIONS

		Pressure		Flujo (L/min)		Temperatura
Nominal pressure	Seal pressure	Working pressure	Ducha	Lavaojos	Agua tibia	
0.4	0.45	0.20-0.40	≥ 75.5	≥ 11		

- ◆ Made of stainless steel resistant to corrosion.
- ◆ The base can be fixed to the floor to assure a high stability.
- ◆ Wide diameter shower head that provides a high and circular water flow.
- ◆ High flow sprays of low pressure water to avoid damages. Protected from dust and other contaminants with caps that are removable without the intervention of user.
- ◆ The eye washer unit is designed with enough space to keep eyes opened with the help of hands while water flows towards the eyes.
- ◆ The valves are automatically opened when pressing the lateral lever (eye washer) or pulling the handle (shower); they allow a permanent water flow until they are closed manually by pressing the lever to the opposite direction or by pushing the handle upwards.

3. INSTALLATION

Preliminary inspection

- ◆ Unwrap the shower and eye washer, take off all the protective items and make sure that it does not present any damage because of the shipment. In case it presents any damage tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.
- ◆ Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.
- ◆ Please check that all the parts are enclosed:
 - Shower head
 - Eye washer head with two sprays