**BAÑOS POR ULTRASONIDOS
ULTRASONIC BATHS
BAIN À ULTRASONS**

Serie / Series / Séries / 610-620-626 (Digital)

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Thank you for choosing this equipment. We sincerely wish that you enjoy your Nahita ultrasonic baths. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products according to the CE marking regulations as well as emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

Please bear in mind the following:

- ◆ This manual is inseparable from the Nahita ultrasonic bath series 610, 620, so it should be available for all the users of this equipment.
- ◆ Follow all required applicable laws about safety and health (electrical security devices, inflammable, toxic and/or pathological products, etc.).
- ◆ You should carefully handle the equipment avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the equipment to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the ultrasonic bath.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to asistencia@auxilab.es.
- ◆ This equipment is protected under the Warranties and consumer goods regulation (10/2003).
- ◆ Overhaul is not covered by the equipment warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the warranty.
- ◆ Neither fuses nor accessories, including their loss, are covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

 **ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

INDEX OF LANGUAGES

Spanish 2-8
 English..... 9-15
 French 16-22

INDEX OF CONTENTS

1. GENERAL INFORMATION 10
 2. DESCRIPTION 11
 3. TECHNICAL SPECIFICATIONS 12
 4. INSTALLATION / SETTING UP 12
 5. MAINTENANCE AND CLEANING..... 14
 ANNEX I: CE CERTIFICATE 23

1. GENERAL INFORMATION

Nahita ultrasonic baths allow cleaning any kind of instruments, even those complicated pieces, without the need to dismantle them; cavitation will penetrate everywhere where the cleaning solution is in contact with the piece's surface.

With Nahita ultrasonic baths, only a few minutes are necessary to get a perfect cleaning of holes or inaccessible cavities of different material; this involves a time saving compared with the traditional cleaning methods. Nahita ultrasonic baths can be used to obtain an absolute cleaning in:

- ◆ **Laboratories:** scalpels, ground glass, filters, syringes, sieves, chromatography columns, optics, etc.
- ◆ **Industry:** semiconductors, precision mechanical pieces, etc.
- ◆ **Others:** instrumental used in odontology, medicine, optics, clockmaking, etc.

Ultrasonic cleaning is based on the use of high frequency waves (40 / 50 Hz) that produce a cavitation effect inside the tank that results in the generation of millions of low pressure microscopic bubbles that exert a molecular cleaning; thus removing impurities, contaminants and dirt from the pieces or material that must be cleaned.

Nahita ultrasonic baths are made of metal body with stainless steel tank. Those models with a capacity more than 6 L present a device to empty the tank.

- ◆ The bath is supplied with a Schuko standard wire; it should be connected to a current wire provided with an earth wire, and should be handy to be disconnected in case of emergency.
- ◆ Never try to repair the bath by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the bath (burns, hurts...) or damages in nearby equipments.
- ◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.

Cleaning

- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the bath and produce an early ageing of the equipment.
- ◆ Use a fluff-free cloth dampened with soaped water that does not contain abrasives.

 **ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.**

INSTRUCTIONS ON ENVIRONMENT PROTECTION

 At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

Depending on the series of ultrasonic baths:

Series 620: it is provided with two different intensities of ultrasonic frequencies. To select one or the other press ON/OFF (1.2).

Series 610: it is provided with heating function. To switch this function on press ON/OFF (2).

Series 626: it is also possible to select both working temperature (2.2) and power (2.3) as well as DEGAS (2.4) function if necessary. SV/PV button shows programmed T and real T alternately.

Security

- ◆ The bath must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ◆ Put the bath on top of a horizontal, plane and stable table making a free space of at least 30 cm per side.
- ◆ Do not put the bath near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc. Avoid vibrations, dust and dry environments.
- ◆ During operation, dangerous material such as flammable or infectious liquids must be out of the working area.
- ◆ When you are not using the bath for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece (e.g. replacement of a fuse).
- ◆ Never try to repair the bath by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Avoid the entry of liquid into the control panel, though it is properly insulated from the bath's tank. In case you have any suspect that any liquid has entered, do please immediately unplug the equipment (see Maintenance).
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

5. MAINTENANCE AND CLEANING

To get the best results and a higher duration of the ultrasonic bath it is essential to follow the processes of use.

Note: All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

Please follow the processes of use of this manual


- ◆ This manual should be available for all users of this equipment.
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.

2. DESCRIPTION

610 and 620 series. Analogical

1.1 Timer

1.2 ON/OFF: ultrasonic frequency (series 620) or heating (series 610)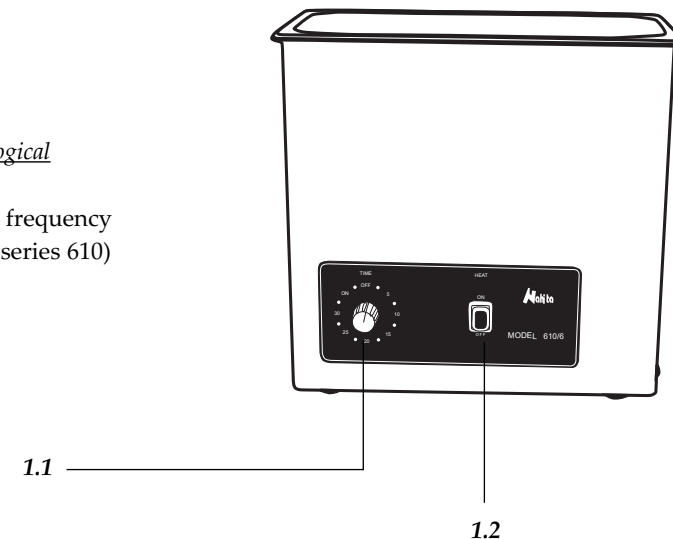



626 Series. Digital

2.1 Time display and selector / Start / Stop Button

2.2 Temperature display and selector / SV / PV Button

2.3 Power selector 0-100%

2.4 DEGAS Selector

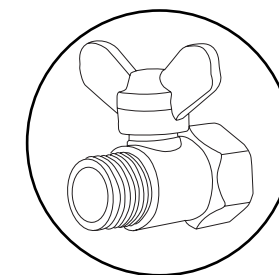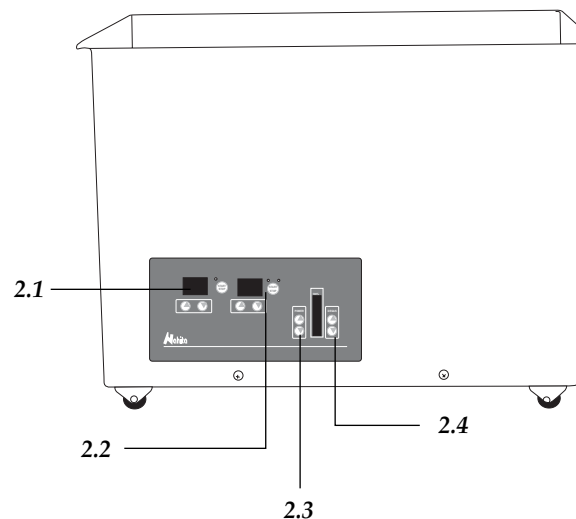Example of the drainage valve (models with more of 6 L).

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Code	Model	Capacity	Useful dimensions (mm)	Timer	Heating	Power	Drainage valve
Series 626 Digital							
50626003	626/3	3 L	235x135x100	1-99min	RT-95°C	120 (0-100%)	-
50626006	626/6	6 L	300x150x150	1-99min	RT-95°C	150 (0-100%)	-
50626010	626/10	10 L	300x250x150	1-99min	RT-95°C	260 (0-100%)	Yes
50626020	626/20	20 L	500x300x150	1-99min	RT-95°C	500 (0-100%)	Yes
Series 620 Analogical							
50620002	620/2	2 L	150x140x100	1-30 min	--	50 W	-
50620003	620/3	3 L	240x140x100	1-30 min	--	100 W	-
50620006	620/6	6 L	300x150x150	1-30 min	--	400 W	-
50620010	620/10	10 L	300x250x150	1-30 min	--	250 W	Yes
50620022	620/20	22 L	500x300x150	1-30 min	--	600 W	Yes
Series 610 Analogical							
50610002	610/2	2 L	150x140x100	1-30 min	~70°C (fixed)	50 W	-
50610003	610/3	3 L	240x140x100	1-30 min	~70°C (fixed)	100 W	-
50610006	610/6	6 L	300x150x150	1-30 min	~70°C (fixed)	400 W	-
50610010	610/10	10 L	300x250x150	1-30 min	~70°C (fixed)	250 W	Yes
50610022	610/20	22 L	500x300x150	1-30 min	~70°C (fixed)	600 W	Yes

4. INSTALLATION / SETTING UP

Preliminary inspection

- ◆ Unwrap the bath and take off the polyspan protection in which it comes fitted. Take off the involving plastic and without connecting the bath to the power supply, make sure that the equipment does not present any damage because of the shipment. In case it presents any damage, tell it immediately to your transport agent or dealer so that they can make the claims in the correct time limit.
- ◆ Please keep the original wrapping; you will always need it for returns enclosed with all the accessories supplied.
- ◆ Please check that all the accessories are enclosed with the equipment:
 - Schuko wire
 - Lid and basket
 - User's manual
 - Warranty certificate

We will only accept any equipment return within 15 days after delivery and provided it comes in its original wrapping.

Installation

Before using the bath, it is convenient for you to familiarize with its components and basic essentials, as well as with its control functions.

PLEASE READ THOROUGHLY THE INSTRUCTIONS BEFORE CONNECTING AND OPERATING WITH THIS EQUIPMENT WITH THE AIM OF ACHIEVING THE HIGHEST FEATURES AND THE MAXIMUM DURABILITY OF THE INSTRUMENT.

- ◆ Please put the bath on top of a horizontal, plane and stable table making a free space of at least 30 cm per side. Do not put the bath near any warm supply (burners, blowlamps...), nor expose it directly to the sun, etc. During operation, dangerous material such as flammable or infectious liquids must be out of the working area.
- ◆ The bath is supplied together with a Schuko standard wire. Please insert the wire that feeds the AC electric current in the base of current 220 V, 50 Hz \pm 10% provided with earth wire and to the other end to the bath connector.

Neither the manufacturer nor the distributor will assume any responsibility for the damages produced to the equipment during its installation or damages to people suffered by the improper use of the electric connection. The tension should be 220 V, 50 Hz \pm 10%.

Setting up

- ◆ Important! Before connecting the equipment to the electrical current, make sure that the tank is filled with cleaning solution (water, solvent, etc.) up to 2/3 of its total capacity.

NOTE: Take into account that the water level will rise when the material to be cleaned is introduced.

- ◆ Nahita recommends the use of the Nahita cleaning solution for ultrasonic baths (code 90600900). Do not use abrasive solutions, the detergent must be phosphate free; the use of inappropriate solvents will damage the stainless steel tank and produce a loss of the product warranty.
- ◆ Connect the ultrasonic bath to the electrical current (the tension should be 220-240 V, 50-60 Hz).
- ◆ To start the operation select the time with the timer (1.1 or 2.1 depending on which models). The operation time will depend on the dirtiness of the objects to be cleaned; once this time has passed the equipment will be automatically turned off.