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6. TROUBLESHOOTING

FAILURE	PROBABLE CAUSE	RESOLUTION
The centrifuge won't run	1. The power cord is unplugged 2. Motor carbon brushes lost conductivity because of a long term inventory 3. Foreign objects wind around the rotor shaft 4. The motor is near the end of its expected life	1. Plug the power cord in 2. Turn the rotor five or six times gently by hand 3. Remove the foreign objects after unplugging the centrifuge 4. Call Nahita technical assistance department
Abnormal noise The centrifuge vibrates abnormally	1. Imbalanced loads 2. Foreign objects are sticking on the rotor or chamber 3. Defective motor	1. Balance the load properly 2. Remove the foreign articles after unplugging the centrifuge 3. Call Nahita technical assistance department
The rotor won't be installed properly	1. The rotor is seated upside down 2. Damaged or deformed rotor or motor shaft boss	1. Seat the rotor properly 2. Call Nahita technical assistance department

INSTRUCTIONS ON ENVIRONMENT PROTECTION

At the end of its life cycle, please, does not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

Thank you for choosing this equipment. We sincerely wish that you enjoy your mini centrifuge Nahita 2507. We highly recommend looking after this equipment according to what is stated in this manual.

Nahita develops its products emphasizing the ergonomics and security for its user.

The correct using of the equipment and its good quality will permit you to enjoy this equipment for years.

The improper use of the equipment can cause accidents and electric discharges, circuit breakers, fires, damages, etc. Please read the point of Maintenance, where we expose the security notes

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

Please bear in mind the following:

- ◆ This manual is inseparable from the centrifuge Nahita series 2507, so it should be available for all the users of this equipment.
- ◆ Follow all required applicable laws about safety and health (electrical security devices, inflammable, toxic and/or pathological products, etc.).
- ◆ You should carefully handle the centrifuge avoiding sudden movements, knocks, free fall of heavy / sharp objects on it. Avoid spilling liquids inside the equipment.
- ◆ Never dismantle the different pieces of the centrifuge to repair it yourself, since it could produce a defective use of the whole equipment and a loss of the product warranty, as well as injuries on people that handle the centrifuge.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ If you have any doubt about setting up, installation or functioning do not hesitate in contacting your wholesaler. You can also tell us any doubts or suggestions you have by contacting Nahita Technical Assistance Department by email to asistencia@auxilab.es.
- ◆ This equipment is protected under the Warranties and consumer goods regulation .
- ◆ Overhaul is not covered by the centrifuge warranty.
- ◆ Operations made by non-qualified staff will automatically produce a loss of the centrifuge warranty.
- ◆ Accessories (including their loss), are not covered by the product's warranty. The warranty neither covers piece's deterioration due to the course of time.
- ◆ Please make sure you keep the invoice, either for having the right to claim or asking for warranty coverage. In case you have to send the equipment to Nahita Technical Assistance Department you should enclose the original invoice or a copy as guarantee.
- ◆ Manufacturer reserves the right to modify or improve the manual or equipment.

ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

INDEX OF LANGUAGES

Spanish.....	2-10
English.....	11-18
French	19-26

INDEX OF CONTENTS

1. USES OF THE INSTRUMENT.....	12
2. DESCRIPTION	13
3. TECHNICAL SPECIFICATIONS	13
4. INSTALLATION / SETTING UP.....	14
5. MAINTENANCE AND CLEANING.....	16
6. TROUBLESHOOTING	18

1. USES OF THE INSTRUMENT

Mini centrifuges series 2507 are compact table top centrifuges, ultra quiet, very stable and easy to use. Great versatility of use due to the different rotors and adapters for use with Eppendorf tubes 0.2, 0.5 and 1.5 mL making it ideal for techniques such as PCR, HPLC or microfiltration.

rotor is damaged, do not use the centrifuge and call Nahita technical assistance department.

- ◆ Inspect the lid is able to open and close smoothly, and inspect the lid is locked when pressing down.
- ◆ Always use original components and supplies. Other devices can be similar but they can damage the equipment.
- ◆ Never try to repair the centrifuge by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the centrifuge (burns, hurts...) or damages in nearby equipments.
- ◆ Inspect the power plug and outlet for a loose connection and dust.
- ◆ In the event of breakdown please contact your distributor to overhaul through Nahita Technical Assistance Department.

Cleaning

- ◆ Check and clean your centrifuge at regular intervals or immediately after spills occur to prevent premature wear and safety hazards
- ◆ Never use scourers or substances that can grate for cleaning metallic parts such as stainless steel, aluminium, coatings, etc. as they damage the centrifuge and produce an early ageing of the equipment.
- ◆ Wipe the rotor with a soft cloth moistened with a neutral detergent with extreme care and remove the detergent with a clean clamp cloth.

ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.

Security

- ◆ The centrifuge must be used by previously qualified staff that knows how the equipment works thanks to the user manual.
- ◆ With wet hands do not touch the plug nor the power switch.
- ◆ Follow all required applicable laws about safety and health (electrical security devices, inflammable, toxic and/or pathological products, etc.).
- ◆ Do not place the centrifuge near any warm supply (burners, blowlamps, etc), nor expose it directly to the sun. Avoid vibrations, dust and dry environments.
- ◆ To prevent fire or electric discharges avoid dry or dusty environments. In case it may happen unplug the equipment immediately.
- ◆ Keep the centrifuge out of direct sunlight, dust and poorly ventilated areas.
- ◆ The environmental conditions required are: room temperature: 5 to 35 °C and relative humidity: 30 to 85 %.
- ◆ During its functioning dangerous materials such as flammable or pathological substances must be out of the safety area.
- ◆ When you are not using the centrifuge for a long period of time please make sure it is unplugged in order to avoid possible accidents.
- ◆ It is essential to have the equipment switched off and unplugged from the net before cleaning, checking components or replacing any piece.
- ◆ Never try to repair the centrifuge by yourself, since you will lose the warranty and may provoke damages to the general operating system or the electrical installation, as well as injuries to the people that usually handle the equipment (burns, hurts...).
- ◆ Try not to spill any liquid on the control panel, though it is properly insulated. In case you have any doubts do please immediately unplug the centrifuge (see Maintenance).
- ◆ The centrifuge is supplied with a Schuko standard wire. It has to be plugged to an earth connection and the socket should be handy and ready to unplug the equipment in case of emergency.
- ◆ Made under the European regulations for electrical security, electromagnetic compatibility and security on machines.

5. MAINTENANCE AND CLEANING


All the processes of use mentioned below will not have any value unless you keep a continued and careful maintenance.

To get the best results and a higher duration of this equipment it is essential to follow the processes of use:


Please follow the processes of use of this manual.

- ◆ This manual should be available for all users of this equipment.
- Visually inspect the rotor for damage such as dents or chips from being dropped. If the

2. DESCRIPTION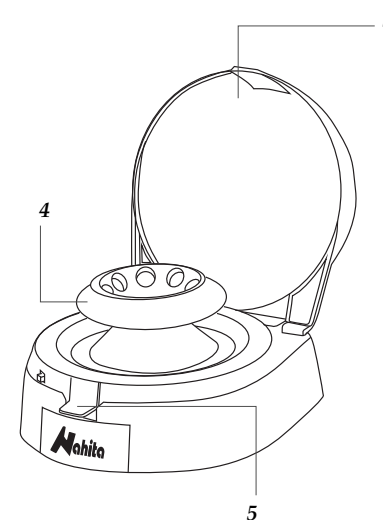

- 3. Lid
- 4 Rotor
- 5. Opening lid

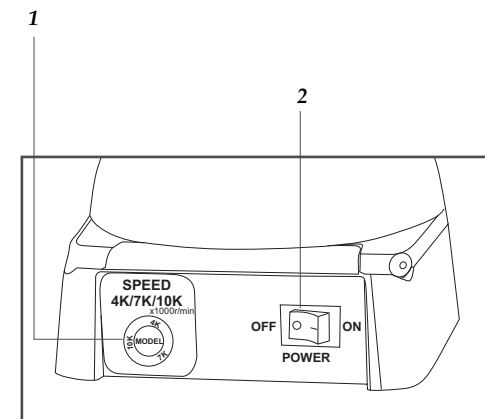- 1. Speed button (only code 52507015)
- 2. ON / OFF

3. TECHNICAL SPECIFICATIONS

Code	52507007	52507011	52507015
Model	2507/7	2507/10	2507/15
Capacity			
Angle rotor	8x1,5 mL 8x0,5 mL* 8x0,2 mL*		
Strip rotor	2 tiras de 8 x 0.2 mL		
Speed	7200 rpm	10000 rpm	4000/7000/10000 rpm
FCR	3260 g	5610 g	990/ 2910/ 5610 g
Dimensions	155 x 175 x 125 mm		
Weight	±0,8 Kg		
Power supply	AC 85-240V, 50/60Hz		

**Using the included adapters