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INDEX OF CONTENTS

PREFACE AND WARRANTY.....	5
1. SAFETY INSTRUCTIONS.....	5
2. PRODUCT DESCRIPTION	6
3. OPERATING INSTRUCTIONS.....	7

PREFACE AND WARRANTY

Thank you for purchasing a Nahita Blue mini magnetic stirrer. Users should read this manual carefully, follow the instructions and procedures, and beware of all the cautions when using this instrument.

If it is necessary, contact your dealer to get the necessary help and advises for a proper use of the equipment. For this, user must give information about the serial number of the equipment, the problems found, the procedures followed to try to solve the problem and contact information.

Nahita Blue mini magnetic stirrer is warranted to be free from defects in materials and workmanship under normal use and service, for a period of 24 months from the date of invoice. It shall not apply to any product or parts which have been damaged on account of improper installation, improper connections, misuse, accident or abnormal conditions of operation.

For claim under the warranty please contact your distributor in order to send the equipment to Nahita Technical Service Department.

1. SAFETY INSTRUCTIONS

Read the user's manual before operating with the equipment. This equipment must be used only by qualified staff.

This equipment is forbidden to use with inflammable, explosive, poisonous and strong corrosive solutions.

Connect the device to an earthed power supply to ensure safety of machine and experiment; the voltage indicated on the equipment must correspond to that of the power supply used.

Non-authorized personnel are not allowed to disassemble and repair this machine. If the equipment presents a problem, contact to your distributor

Make sure the equipment is placed and installed in a plain, horizontal and stable surface.


Pay attention to overloading of material in relation to the capacity of the equipment. An overloading could produce overheating and damages in the equipment.

2. PRODUCT DESCRIPTION

<i>Speed range</i>	100-2000rpm
<i>Maximum volume of stirring</i>	1L
<i>Plate dimensions</i>	110mm
<i>Power</i>	AC adapter 220V/12V
<i>Dimensions (LxWxH)</i>	120x120x50mm
<i>Weight</i>	0.4Kg


FRONT

- 1.1. Off/on switch
- 1.2. Speed knob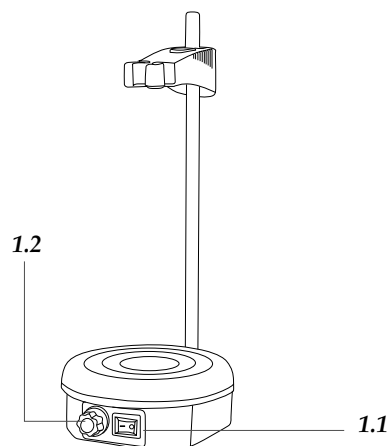

BACK

- 2.1. Rod holder
- 2.2. Adaptor plug

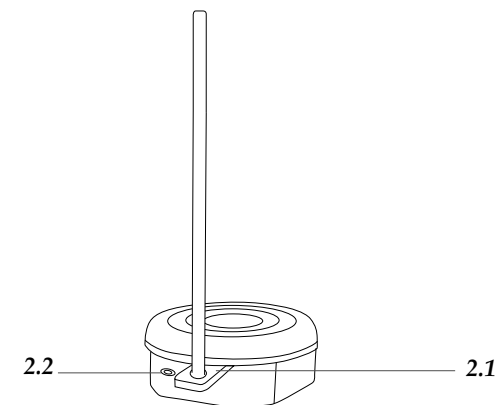Accessories:

- Power adapter (output 12V; input 220V)
- Rod and support
- Magnetic bar

3. OPERATING INSTRUCTIONS

- ◆ Screw the rod with the support in the rod holder of the stirrer (2.1).
- ◆ Connect the power adapter to the stirrers adaptor plug (2.2) and then to the main.
- ◆ Introduce the magnetic bar into the vessel where the liquid that is going to be stirred is placed. When you fill the vessel, do not fill it totally and leave enough space to prevent spillage during stirring.
- ◆ Place the vessel on top of the stirrer.
 - Step 1: Turn on the equipment using the off/on switch(1.1).
 - Step 2: Adjust the speed using the knob (1.2), slowly until the required stirring speed is achieved.
- ◆ Maintain it under stirring the necessary time. Once the process has finished, decrease the speed to zero (1.2.) and turn on the equipment (1.1).

INSTRUCTIONS ON ENVIRONMENT PROTECTION

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.