DECLARACIÓN DE CONFORMIDAD CE
CE DECLARATION OF CONFORMITY
DÉCLARATION DE CONFORMITÉ CE

El fabricante | The manufacturer | Le fournisseur:
AUXILAB, S.L.

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Thank you for purchasing a Nahita Blue multi-function vortex mixer. Users should read this manual carefully, follow the instructions and procedures, and beware of all the cautions when using this instrument.

If it is necessary, contact your dealer to get the necessary help and advises for a proper use of the equipments. For this, user must give information about the serial number of the equipment, the problems found, the procedures followed to try to solve the problem and contact information.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

Nahita Blue stirrers are warranted to be free from defects in materials and workmanship under normal use and service, for a period of 24 months from the date of invoice. It shall not apply to any product or parts which have been damaged on account of improper installation, improper connections, misuse, accident or abnormal conditions of operation.

For claim under the warranty please contact your distributor in order to send the equipment to Nahita Technical Service Department.

ATTENTION!! IF EQUIPMENTS ARE NOT PROPERLY CLEAN AND DISINFECTED THEY WOULD NOT BE ALLOWED TO REPAIR BY OUR TECHNICAL SERVICE.


INDEX OF LANGUAGES

Spanish	2-8
English	9-15
French	16-22

INDEX OF CONTENTS

1. DESCRIPTION	10
2. SAFETY INSTRUCTIONS	10
3. INSPECTION.....	11
4. OPERATING MODE.....	13
5. CHANGE OF ATTACHMENT	13
6. TECHNICAL SPECIFICATIONS.....	14
7. MAINTENANCE AND CLEANING	15
ANNEX I: CE CERTIFICATE.....	24

1. DESCRIPTION

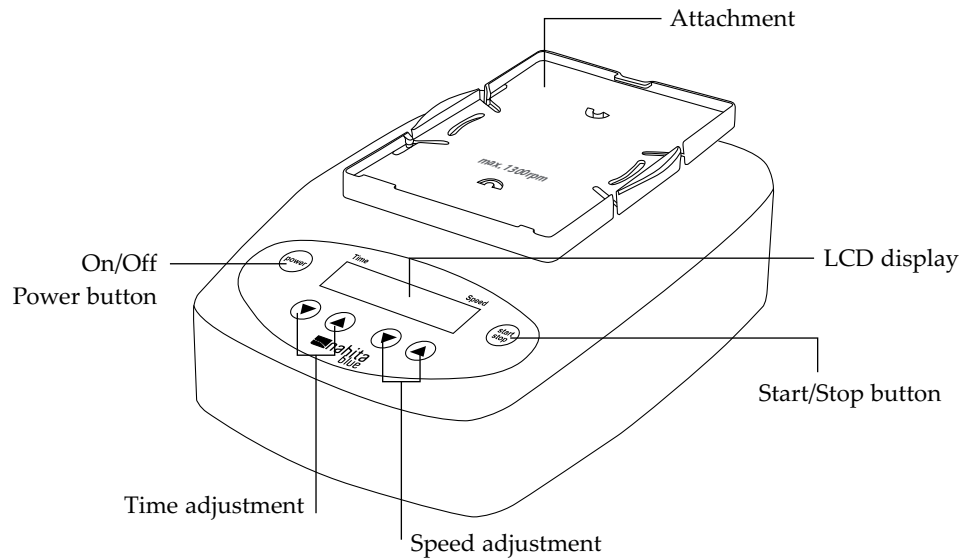2. SAFETY INSTRUCTIONS

- ◆ Read the user's manual before operating with the equipment. This equipment must be used only by qualified staff.
- ◆ This manual should be available for all users of this equipment.
- ◆ Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations indicated in this manual.
- ◆ The voltage stated on the label must correspond to the main power supply.
- ◆ The device must only be operated with the original plug-in power supply unit.
- ◆ The instrument can only be disconnected from the main power supply by pulling out the connector plug, not the cable.
- ◆ During operation wear the necessary personal protections according to the hazard category of the medium to be processed; otherwise there can be risk of:
 - Splashing liquids
 - Projected particles
 - Body parts, hairs, clothing or jewels to get caught.
- ◆ Make sure the equipment is placed and installed in a spacious area on an even, stable, clean, non-slip, dry and fire-proof surface.
- ◆ Respect the minimum safety distances between devices, between the device and the wall and above the assembly (min.100 mm).

7. MAINTENANCE AND CLEANING

The equipment is maintenance-free.

- ◆ Do not allow moisture to get into the equipment when cleaning.
- ◆ Unplug the power line when cleaning.
- ◆ Use only the following products:
 - For dyes: Isopropyl alcohol
 - For construction materials: Water containing tenside / Isopropyl alcohol
 - For cosmetics: Water containing tenside / Isopropyl alcohol
 - For food: Water containing tenside
 - For fuels: Water containing tenside
- ◆ Wear the proper protective gloves during cleaning of the instrument.
- ◆ Antes de utilizar algún otro método de limpieza o descontaminación, contacte con su distribuidor para evitar dañar el equipo.
- ◆ The instrument must be cleaned and put it into the initial packaging carton before sending to service for repair, avoiding the contamination of hazardous.
- ◆ Use the instrument in a dry clean room and stable temperature environment.

INSTRUCTIONS ON ENVIRONMENT PROTECTION

At the end of its life cycle, please, do not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non adequate disposal would damage the environment.

The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.

Please inquire at the community administration for the authorized disposal location.

6. TECHNICAL SPECIFICATIONS

Power supply

<i>Input</i>	100-240 V AC; 1000 mA; 50-60Hz
<i>Output</i>	24 V DC; 24 W.LPS; 50-60 Hz

Agitador

<i>Operating voltage</i>	24 V DC
<i>Power consumption, normal operation</i>	20 W
<i>Power consumption, standby operation</i>	2 W
<i>Motor-output power</i>	8 W
<i>Drive</i>	Brushless
<i>Speed</i>	300-3000 rpm
<i>Orbital diameter</i>	4.5 mm
<i>Shaking motion</i>	Orbital
<i>Timer</i>	
<i>Mode seconds</i>	1-59 s
<i>Mode minutes</i>	1-59:99 min
<i>Mode hours</i>	1-24 h
<i>Mode continuous</i>	Continuous operation
<i>Display</i>	LCD
<i>Environ. temperature of use</i>	5-40°C
<i>Environ. relative humidity</i>	80% RH
<i>Protection grade</i>	IP21 (according to DIN EN 60529)
<i>Contamination</i>	Level 2
<i>Overvoltage category</i>	II
<i>Operation at terrestrial altitude</i>	max. 2000 m
<i>Dimensions</i>	210x150x65 mm
<i>Weight</i>	2.5 Kg
<i>Max. Shaking weight</i>	0.5 Kg (with attachment)

- ◆ The feet of the equipment must be clean and not damaged.
- ◆ If the shaking movement of the equipment increases (resonance), reduce the speed or pass through the critical phase as quickly as possible.
- ◆ Reduce the speed if:
 - The liquid splashes out of the vessel because the speed is too high.
 - The equipment does not operate smoothly.
 - The equipment moves because of dynamic forces.
- ◆ Secure firmly the attachments and vessels in the corresponding place; otherwise the vessel could be damaged or projected out.
- ◆ Check the equipment and accessories before each use. Never use damaged components.
- ◆ Place one single vessel in the center or several vessels evenly spread out.
- ◆ Vessels with sharpened edges cause wear debris on the attachments.
- ◆ Beware the risk of:
 - Flammable materials
 - Glass breakage as a result of shaking mechanical forces.
- ◆ Only process material that will not react dangerously with the extra energy produced during shaking. This also applies to any extra energy produced by other methods, as for example light exposure.
- ◆ Do not use the equipment in explosive atmospheres, with hazardous substances or under water.
- ◆ Safe operation is only guaranteed with the accessories described on the "Supplied accessories" chapter.
- ◆ Always disconnect the equipment before fitting accessories.
- ◆ The appliance does not start up again automatically following a cut in the power supply.
- ◆ Protect the equipment and accessories from knocks and impacts.
- ◆ The equipment may heat up during use.
- ◆ Non-authorized personnel is not allowed to disassemble and repair this machine. If the equipment presents a problem, contact to your distributor.
- ◆ Keep away from high magnetic field.

3. INSPECTION


Unpack the equipment carefully and check for any damages which may have arisen during transport. If it happens, please contact your distributor.

Note: If there is any apparent damage on the equipment, do not plug it into the power line.



Supplied accessories

Nahita Blue multi-function vortex mixer is supplied with the following accessories:



<i>Accessory</i>	<i>Description</i>	<i>Continuous mode</i>	<i>Touch mode</i>	<i>Speed</i>
Multi-function mixer	Mixer	Yes	Yes	300-3000
Standard attachment	Test tubes and small vessels; $\varnothing$ up to 50 mm	Yes	Yes	300-3000
Round attachment	For the different rubber foam inserts	No	Yes	300-1300
Microplate attachment	For 1 microplate	No	Yes	300-1300
Microdish insert	For inserting in the round attachment	No	Yes	300-1300
Test tube insert 1	Holes can be made depending on user necessities	No	Yes	300-1300
Test tube insert 2	For 14 test tubes; $\varnothing$ 10mm	No	Yes	300-1300
Test tube insert 3	For 6 test tubes; $\varnothing$ 12 mm	No	Yes	300-1300
Test tube insert 4	For 4 test tubes; $\varnothing$ 16 mm	No	Yes	300-1300
Power supply	Input: 100-240V AC 50-60Hz Output: 24V DC; 1000 mA	—	—	—
Connector plug	Chinese standard	—	—	—
Connector plug	American standard	—	—	—
Connector plug	European standard	—	—	—
Connector plug	British standard	—	—	—

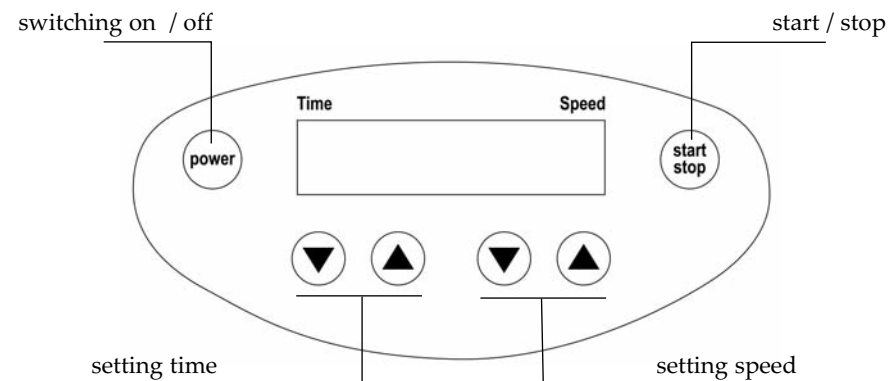4. OPERATING MODE**5. CHANGE OF ATTACHMENT**

Standard attachment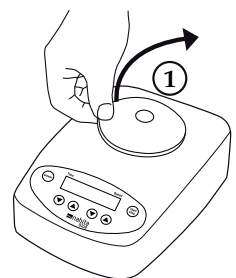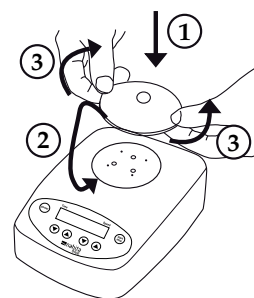

Round attachment

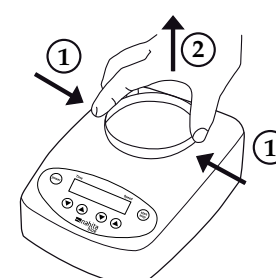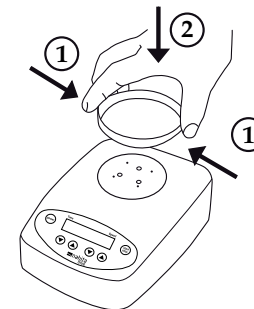Microplate attachment